

# Design of a Device to Cyclically Stretch Isolated Cells

Colleen Brinkmann

Anthony Corbo

Amy Morin

**Gregory Trumbull** 

Advisors: Professor Kristen Billiar

Professor Domhnull Granquist-Fraser

## **Table of Contents**

## **Contents**

Table of Contents	1
Table of Figures and Tables	3
Abstract	4
Authorship Page	5
1 – Introduction	6
2- Background	8
2.1- Mechanobiology	8
2.1.1 External Mechanical Forces	10
2.1.2 Internal Mechanical Forces	10
2.1.3 Mechanobiological Responses	11
2.1.4 Cellular Chemical Signaling (Mechanotransduction)	14
2.2 Cell Stretch Devices	14
2.2.1 Pressure to Achieve Stretch	14
2.2.2 Vacuum to Achieve Stretch	15
2.2.3 Motor Driven Stretch Devices	20
3 – Design Approach	22
3.1 – Design Alternatives	22
3.1.1 – Four Motors System	23
3.1.2 – Two Motors Systems	24
3.1.3 – Vacuum (Flexcell)	28
3.1.4 – One motor system	28
3.2 – Design Process	30
3.3- Final Design	32
3.4 - Device Description	33
3.4.1 – Motor Selection	34
3.4.2 – Gear and Pulley System	36
3.4.3 – Linear Guide System to Directly Stretch Strex Well	37
3.4 – Microscopy Methods Currently Used	42

Objectives	42
Constraints	43
Best-Fit Microscope	43
3.5 – Incubation Methods Currently Used	44
4 – Device Validation	46
4.1 Test Trials – High Density Mapping	46
4.2 Device Validation	49
5 – Conclusions and Recommendations	54
5.1 Conclusions	54
5.2 Recommendations	54
References	58
6.1 – Appendix A – Pairwise Comparison Chart	61
6.2 – Appendix B – Functions Means Tree	62
6.3 – Appendix C- Budget Breakdown	63
6.4 – Appendix D- Strain Parameter Examples	64
6.5 - Appendix E- General Use Manual	65
6.6 – Appendix F- Design Drawings	73
6.7 – Appendix G- Ordered Parts	83
6.8 - Appendix H- Custom Parts	104
6.9 - Appendix I- HDM Matlab Compile Function	122
6.9 - Appendix J - HDM Displacement in the X Direction Raw Data	127
6.11 - Appendix K- HDM Displacement in the Y Direction Raw Data	133
6.12 - Appendix L- HDM Strain in the X Direction Raw Data	136
6.13 - Appendix M- HDM Strain in the Y Direction Raw Data	139
6.14 – Appendix N- HDM Shear Strain Raw Data	142
6.15 – Appendix O- Instron Data of Strex Well Stretch	145
6.16 – Appendix P- Cells imaged on the stretch device	146

## **Table of Figures and Tables**

Figure 2.1: Schematic illustration of the "mechanical nature" of cellular mechanotransduction	on
mechanisms (Taylor and Francis, 2008).	13
Figure 2.6: Strex Cell Stretch Device (B-Bridge International, Inc. 2009).	21
Figure 3.1 3-D Model Drawing of Strex design (top) and view of Strex device (bottom)	25
Figure 3.2: Linear Bearings One Motor design (equibiaxial set up)	29
Table 3.1 Comparison of Strex and One motor design	30
Figure 3.2: Design in equibiaxial (top) and uniaxial (bottom)CAD	32
Figure 3.3: Our device on the Zeiss microscope (top) and close up with cells seeded on the	well
	33
Figure 3.4 –Exploded View of Crank Shaft	37
Figure 3.5 – Igus Linear Rail and Carriage System	39
Figure 3.6 – Four Stretch Arms	40
Figure 3.7 – Exploded View of X and Y Guided System	41
Figure 3.8 – Exploded View of Corner Assembly	42
Figure 4.1: Hand stretched Strex well to 20%	47
Figure 4.2: HDM results in x axis (top), y axis (middle) and shear strain (bottom)	48
Figure 4.3: Stretch Device on Zeiss Microscope	50
Figure 4.4: Rat Aorta Cells at 40X magnification with Hoescht and Phase Views	51
Table 4.1: Heat Test Results	52
Figure 4.4: Graphical Results of Heat Test Over Time	53
Figure 5.1: Heim Joints (Lunsford, 2010).	56
Figure 5.2: Igus plain linear bearing (left) and Igus enclosed linear bearing (right) (Igus, 202	10).
	57

#### **Abstract**

Cells in the musculoskeletal system are constantly subjected to numerous mechanical forces in vivo. Years of research in the field of mechanobiology has shown mechanical forces, including tension and compression, significantly impact various cellular functions such as gene expression, cell proliferation and differentiation, and secretion of matrix proteins. Mechanical signals are also converted into a cascade of cellular and molecular events, initialized by mechanotransduction mechanisms. Multiple commercially produced and custom-built devices are available to study cellular responses to substrate strain. However, real-time data analysis of dynamic cellular responses is not available or easily feasible. WPI has both commercially manufactured and custom built stretch devices available and these devices are currently being used by Professor Billiar. However, they unfortunately do not allow for real-time analysis of cellular response to mechanical stimuli. We have developed a novel stretch device optimized for live-cell imaging. The device assembles on a standard inverted Zeiss microscope and can apply constant cyclic stretch for extended periods of time on the silicone STREX 4cm<sup>2</sup> X-Y culture wells (B-Bridge International). The magnitude and rate of stretch are variable, with maximum stretch of 20%, and a stretch frequency of 1.9 Hz (114 rpm). Interchangeable arms, and a movable linear guide allow for equibiaxial and strip biaxial (pure uniaxial) stretch profiles. Strain analysis achieved by HDM validates and shows reproducible applications of stretch. This device will be used in the future to study molecular dynamics, and cell response to mechanotransduction. Post interpretations verify the stretch device is a novel tool and is compatible with live-cell microscopy for studying dynamic structural remodeling and cellular responses under mechanical strain.

## **Authorship Page**

Key: All-Colleen Brinkmann, Anthony Corbo, Amy Morin and Gregory Trumbull, CB-Colleen Brinkmann, AC-Anthony Corbo, AM- Amy Morin, GT- Gregory Trumbull

Sections	<b>Author/Editor</b>
Abstract	AC
Introduction	All
Background	All
Mechanobiology	AC
Cell Stretch Devices	All/AM
Microscopy	GT/AC
Incubation	GT/AC
Design Approach	AM, GT/AC
Design Options	AM, GT/AC
Design Process	GT
Final Design	AM, GT/AC
Results	AC, CB/AM, GT
Design Validation	AC, CB/AM, GT
Conclusions	AM/AC
Recommendations	AC, CB/AM
References	All
User Manual	AC/GT, AM

#### 1 - Introduction

The human body is a dynamic example of constant changes. On a macroscopic level, people grow, age, and mature. On a microscopic level, cells are growing, dying, proliferating, and constantly changing. These cells are responding to various mechanical forces, and the study of mechanobiology suggests that these forces are crucial to the creation of specific tissue types in the body.

Bone cells for instance need to undergo a compressive force in order to create cortical bone that can withstand the normal forces bones are exposed to on a regular basis. Muscle cells also need an external force constantly acting upon them in order to maintain mechanical strength. In order to ensure these tissues are able to function under the same conditions present in the body, the mechanical properties need to be tested. It is important to see how the individual cells respond to this external force because tissues are not isotropic so the force will not be distributed evenly amongst the tissue (Mizutani et. al, 2006).

Ideally, these different mechanical forces created in the body could be replicated in a laboratory setting; however, no machine can function exactly like the body. For quite some time, devices to stretch cells have been created that attempt to mimic the in vivo conditions that would be present in the body. Stresses can be applied to test the mechanical properties in tension, compression, and shear situations. Cells can be stretched in many different directions so that they can experience similar stresses and forces that would be present in the body. The device allows those properties to be tested, and see how much stress or strain in any loading situation the cells can handle. Most devices focus on one specific type of stretch such as uniaxial, equibiaxial or strip biaxial, and some contain viewing platforms to see the individual cell responses to this

stretch while others do not. By stretching the cells various ways, we will be able to view how the cells respond to that stretch and how they change in orientation and other pathways because of that stretch.

In order to keep cells alive during testing, tissue stretching devices have been equipped with various mechanisms to maintain cell homeostasis so a realistic cell response can be seen from the different stretches the cells are subject to. Each individual cell responds to the forces acted on it during a stretch, so it is important to see not only the collective movements of a tissue membrane, but also the movements of each individual cell. This method will help to see if there are any trends in the movement of cells or if some cells experience more stress or strain than others and causes them to respond to stimuli in a different manner than the collective tissue. In order to see these cell responses, the cells need to be examined on the microscopic level while it is experiencing the stretch.

The goal of this project is to design, build, and test a cell stretching device that can stretch cells biaxially from equibiaxial to strip biaxial strain. The device must allow observation of individual cell responses to the stretch applied to the membrane before, during (through the use of cameras), and after the stretch to see how they reorient over that period of time. The machine should stretch along two axes in order to assure an area with uniform strain to see the true movement of the cells to the stretch. This project will be accomplished this with our budget of \$1000 and our time constraint of completion of the project of April 2011.

## 2- Background

### 2.1- Mechanobiology

Mechanical stimuli have been found to play a crucial role in the development of many tissues, and these mechanical stimuli effect tissues at the cellular level (Wang D. et. al, 2010). The interactions at the cellular level can cause tissues to develop differently if they were somehow changed. Different microenvironments have the potential to initiate different signal pathways in cells, which can effect gene expression.

In vitro model systems are commonly used to explore the effects of mechanical stimulus on cells. Generally, deformable materials with smooth culture surfaces are used to apply cyclic mechanical stretching to cells (Wang, J. et. al, 2005). Cyclic stretching model systems have shown mechanical stimulus induces various responses, including cell reorientation, actin cytoskeletal remodeling, and altered cell proliferation, gene expression, and protein synthesis (Wang D. et. al, 2010).

These model systems have some limitations, the most significant limitation regarding the smooth deformable material. Previous cell stretch studies on smooth substrates have shown cells randomly orient initially, and then reorient in a direction with minimal deformation after stretching (Wang, H. et. al, 1995). This induces potential for heterogeneous strain on the cells, which changes depending on their orientation to the stretching direction. To avoid heterogeneous strain on the cells equibiaxial stretch systems have been developed. These systems offer isotropic strain gradients on smooth culture substrates, allowing the cells to experience the same strain despite their orientation (Wang, J. et. al, 2005). However, equibiaxial stretching does not physiologically mimic what many cells experience. For example, tendon and ligament fibroblasts are subjected to uniaxial stretching in vivo (Wang, J. et. al, 2005). Another limitation of these systems is the variation in organization and shape. Cells in vivo are well organized and have

defined shapes, whereas cells in vitro follow a random trend. Considering cell organization and cell shape influence cell function and phenotype, it is quite possible to recognize variation in data (Chen et. al, 1997). Currently to avoid these issues culture systems are being developed which promote/control cell organization, shape and mechanical conditions more precisely.

Cells in the musculoskeletal system are subjected to various mechanical forces in vivo. Years of research have shown that these mechanical forces, including tension and compression, greatly influence various cellular functions such as gene expression, cell proliferation and differentiation, and secretion of matrix proteins. Cells also use mechanotransduction mechanisms to convert mechanical signals into a cascade of cellular and molecular events (Wang, J. et. al, 2010).

Mechanical forces act on humans at different levels, from the body as a whole to individual organs, tissues, and cells. It is well known that appropriate mechanical loads are beneficial to bone and muscle by enhancing their mass and strength. On the other hand, excessive mechanical forces can also be detrimental; for example, excessive mechanical loading of tendons plays a major role in the development of tendinopathy (Fredberg et. al, 2008)(Wang, JH. et. al, 2006). Thus, mechanical forces have a profound effect on tissue homeostasis and pathophysiology. The central players in the human body's response to mechanical forces are various types of mechano-sensitive cells. Examples of such cells include tenocytes in tendons, fibroblasts in ligaments and skin, osteocytes in bone, chondrocytes in articular cartilage, and endothelial cells in blood vessels. Mechanical forces induce a wide range of cellular events, including proliferation, differentiation, and gene and protein expression by both adult differentiated and stem cells (Wang, JH et. al, 2008).

#### 2.1.1 External Mechanical Forces

Two basic ways to apply tensile mechanical forces to cells a substrate such as silicon include uniaxial or biaxial. Uniaxial stretching is appropriate for application of mechanical forces to cells originating from tendons and ligaments; these cells are aligned with the long axis of the tissue and are primarily subjected to uniaxial stretching in vivo (Chen et. al, 1997). On the other hand, biaxial stretching is applied to cells that are subjected to tensile forces in all directions in vivo, such as dermal fibroblasts. Several biaxial stretching systems have been devised, which typically use circular elastic membranes to produce isotropic strains independent of stretching direction (Wang, J. et. al, 2010). In addition to tensile forces, compressive forces can also be applied to cells that are subjected to compression in vivo. A common means of applying compressive forces is through applications of hydrostatic pressure (Wang, J. et. al, 2010). Other techniques use direct platen abutment to apply compressive forces to cells. These types of loading systems include unconfined compression, in which constant or low-cycle intermittent loads are delivered by manually applying weights (Wang, J. et. al, 2010). These compressive loading systems can be used to investigate mechanobiological responses of cells in tissues primarily subjected to compression in vivo, such as articular cartilage.

#### 2.1.2 Internal Mechanical Forces

In the body, cells generate mechanical forces themselves; these forces are commonly referred to as intracellular tension. However, in non-muscle cells, intracellular tension is produced by the cross-bridging of actomyosin. This process is caused by ATP hydrolysis (Wang, J. et. al, 2010). These tensile forces are then transmitted to the ECM via focal adhesions, and the

forces acting on ECM are called cell traction forces (CTFs). CTFs play a critical function in cell mechanobiology, as they direct ECM assembly, control cell shape, permit cell movement, and uphold cellular tensional homeostasis (Balaban et. al, 2001). CTFs also deform the ECM network and cause stress and strain in the network, which in turn modulate cellular functions such as gene expression and protein secretion. Hence, CTFs are critical in many fundamental biological processes such as embryogenesis, angiogenesis, and wound healing (Wang, J. et. al, 2010).

However, mechanobiology studies rely on cell-substrate adhesions which transmit external mechanical forces to cells. This is because external forces acting on cells can alter their internal forces, thus affecting cellular mechanobiological responses. Interestingly, factors such as substrate stiffness also have a significant influence on cell behavior. For example substrate stiffness only can direct specific differentiation of human mesenchymal stem cells (hMSCs); soft substrates (0.1-1 kPa) mimicking brain tissues are neurogenic, whereas stiffer substrates (8-17 kPa) mimicking muscle are myogenic. Finally, even stiffer substrates (25- 40 kPa) resembling osteoid matrix can induce hMSCs to undergo osteogenic differentiation" (Wang, J. et. al, 2010).

#### 2.1.3 Mechanobiological Responses

Depending cell type and loading conditions, various mechanical forces on cells promotes a variety of cellular responses/functions, including cell proliferation, differentiation, gene expression and synthesis of ECM proteins, production of cytokines and growth factors (Wang, J. et. al, 2010). In a study conducted by Rehfeldt et al, human tendon fibroblasts were stretched at different magnitudes and showed increase in proliferation as well as gene expression and protein production of type I collagen (Rehfeldt et. al, 2007). Another study, conducted by Yang et al, cyclically stretched human tendon fibroblasts at a magnitude of 5% and a frequency of 1 Hz for

24 hours, resulted showed significant increases in cell proliferation. However, when the cells were stretched for 48 hours cell proliferations was inhibited, this of course indicating stretching-induced proliferation of tendon fibroblasts depends on stretching duration (Yang, G. et. al, 2004). Another test in human periodontal ligament fibroblasts, conducted at 10% cyclic equibiaxial compression test which decreased type I collagen mRNA expression and reduced synthesis of fibronectin as well as the amount of total protein; however, the same level of cyclic stretching increased type I collagen mRNA levels and total protein levels (He et. al, 2004).

These studies illustrate how tensile and compressive forces with the same magnitude induce differential cellular mechanobiological responses. Other than affecting cell proliferation and protein expression, mechanical forces also promote the expression and production of inflammatory mediators, including COX-2, PGE2, and LTB4, in a stretching magnitude-dependent fashion (Wang, J. et. al, 2010). These results suggest that when tissues such as tendons are injured, appropriate levels of exercise could be beneficial as it may reduce the inflammatory response. On the other hand, excessive loading of injured tendons, which may worsen tissue inflammation, could be detrimental. These various mechanical forces have noteworthy effects on cells, understanding these forces and their affects can eventually lead to a better understanding of cell function, homeostasis and equilibrium. Understanding these can lead to future success in various biomedical applications such as: tissue-engineering, wound healing, cancer treatment, gene therapy and tissue remodeling.

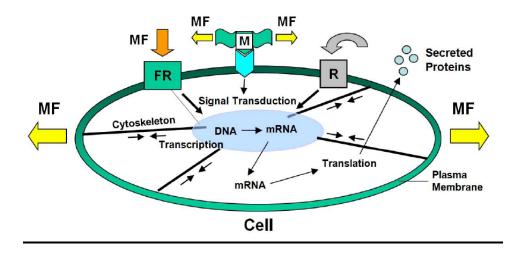


Figure 2.1: Schematic illustration of the "mechanical nature" of cellular mechanotransduction mechanisms (Taylor and Francis, 2008).

Mechanical forces (**MF**) can induce mechanotransduction by directly altering conformation of an extracellular matrix (**ECM**) protein and integrin configuration and transmitting forces to the cytoskeleton and nucleus, thus eventually affecting transcription and translation. Also, mechanical forces can unfold a domain of the extracellular protein (**M**) and expose a cryptic site that may serve as an activating ligand for a cell surface receptor, resulting in a series of signaling events. Also, when mechanical forces are applied to "force receptors" (**FR**), such as integrins and G proteins, they initiate signal transduction, resulting in transcription followed by translation. As a result, soluble factors are secreted into the ECM, which act on the receptor (**R**) and then initiate a cascade of signaling events. Note that double arrows indicate intracellular tensions in the actin filaments (Modified with permission from Wang and Thampatty, Taylor & Francis. 2008).

#### 2.1.4 Cellular Chemical Signaling (Mechanotransduction)

For cells to respond to mechanical forces, these forces must be transformed into chemical signals inside the cell to induce a cascade of cellular and molecular responses. The cellular process of converting mechanical stimulus into chemical signals is called cellular mechanotransduction. Figure 2.1 illustrates and describes this process above. The mechanism of cellular mechanotransduction is still not entirely clear, however, it is generally established that external mechanical forces are transferred to a cell through integrin-mediated adhesions in the ECM (Juliano et. al, 1993). Integrins hold large ECM domains responsible for binding substrates to cytoplasmic domains, and are the primary adhesive receptors and mechanotransducers which link the cytoskeleton to the ECM. Consequently this ECM-integrin-cytoskeleton network plays a key role in mechano-signaling processes. Presence of mechanical stress to integrins is capable of altering the cytoskeleton and activated gene expression in a stress-dependent manner.

#### 2.2 Cell Stretch Devices

In order to determine the various means to stretch cells, an extensive literature review of devices currently in existence was conducted. Each device has its own specific function to obtain specific stretching of cells, and its own specific way to obtain those desired results. The main methods of stretching cells can be broken down into systems using motors, vacuums, and a pressure force.

#### 2.2.1 Pressure to Achieve Stretch

The direct platen abutment approach of compressing the cells uses a manual application of a force directly downward on the membrane to stretch (Brown, 2000). It has been used widely

in collagen applications because it mimics in vivo conditions closely. A device that uses this method has been developed by Tanaka to target specifically the low strain regime in bone. This device uses an actuator to drive the force of the platen to compress the membrane, which in this case is a 3-D bone matrix (Tanaka, 1999). This device runs at a frequency of 100Hz and can produce cyclic loading and normal sinusoidal wave loading, but also non-typical loading (Tanaka, 1999).

Another device design uses pneumatics to compress the cells (Torzilli et al, 1997). Advantages of these devices are the ability to have cyclic loading which produces sinusoidal compression of the membrane or non-sinusoidal loading of the membrane. The forms of stretch it can attain include unconfined compression with no support of the materials from the sides as well as confined compression with support on the sides. This has been shown to be an effective form of stretching cartilage cells to determine metabolic responses (Torzilli et. al, 1996). This study involved stretching of cartilage cells with stresses ranging from 0.5-24MPa were loaded at 1Hz sinusoidal for a 2-24 hour period of time. It is also able to obtain homogeneous compression as in the hydrostatic pressurization mechanism. Disadvantages would be its ability to attain strip biaxial strain, as well as keep the cells alive during the stretch (Brown, 2000).

#### 2.2.2 Vacuum to Achieve Stretch

A widely used means of placing cells under compression is through the use of hydrostatic pressure (Brown, 2000). The pressure can use vacuum pressure which is negative or positive pressure to compress cells, and thus stretch them. There are several devices which have been designed using this basic method, and are useful for their given applications. Some specific advantages of this device are the ability to obtain homogeneous compression, which allows for

equibiaxial strain to be easily attainable. Also, the pressure does not prevent the cells from obtaining any nutrients from the cell media which allows them to remain alive during the stretch (Brown, 2000).

Some disadvantages of this device also come with the homogeneous compression in that it becomes that much more difficult for the cells to attain strip biaxial strain. Modifications to the compression chamber would have to be made for the stretch device to be able to attain strip and equibiaxial strain. Other disadvantages are the use of either pressurized O<sub>2</sub> or CO<sub>2</sub> can prove to be harmful to the cells under strain. In order to keep the cells alive, modifications need to be made to the cell media to compensate for that pressurized environment. Another disadvantage of hydrostatic pressurization as a cell compression method is that cells experience forces regularly that are different than just hydrostatic in nature, which therefore does not serve as a totally useful in vitro model (Brown, 2000).

This form of applying a mechanical load to cells includes circular substrates that have loads applied to them, and subsequently produce the strain on cells. There are four main ways to attain this strain on the cells. A curved platen may be used to provide the uniform strain on the cells (Williams et al, 1992). This study did show that surface strains due to bending, however, are not negligible and result in large strains near the clamped regions of the membrane (Williams et al, 1992). Another way to apply this load is through prong displacement, in which a collagen substrate is loaded and unloaded with a prong being moved by a stepper motor control (Vandenburgh, 1988). The third means of displacing the membrane would be though a fluid movement. The final option for movement of the membrane would be through vacuum displacement, in which the pressure is applied to the underside of the well and the substrate is

then stretched from the pressure (Brown, 2000). These four methods can be seen below in Figure 2.2.

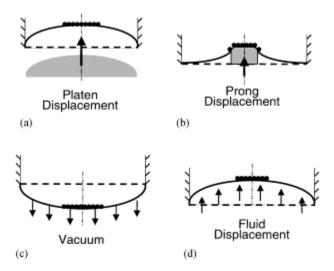


Figure 2.2: Methods of out-of-plane circular substrate distention (Brown, 2000).

There is one commercially available device that actually takes that concept of vacuum circular substrate distention, and that device is called the Flexcell® System. This system attempts to account for the various stresses experienced in the body, and mimic them as close to in vivo conditions as possible using a vacuum pump to push down on the substrate and apply strain (Flexcell International Corporation, 2004).

This device has been used extensively used in literature for various applications. In one application, it was used to test the development of tissue engineered tendons for anterior cruciate ligament replacement (Garvin et al, 2003). This used the Flexcell device to apply strain to the tendon to test the mechanical properties. With this device, the cells were alive with 5% CO<sub>2</sub> and 37°C by placing the base plate in an incubator while still applying the vacuum pressure for a period of 1.5 hours (Garvin et al, 2003). The cells are loaded in a circular six plate well with

anchors on both sides of nonwoven mesh, holes to accommodate the vacuum flow, and the rubber membrane from one side to the other. This can be seen below in Figure 2.3.

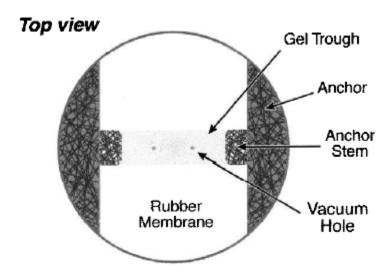


Figure 2.3: Top view of Flexcell well with substrate (Garvin et al, 2003).

The cells were placed in a six well dish with a Delrin Trough Loader insert that allows for the space beneath the culture to be entirely filled. The insert has holes that allow for air to escape from the top of the rubber membrane and push down on the loader (Garvin et al, 2003). Once the cells have been seeded properly on the gel, the trough loader is removed, leaving a space for the deformation of the substrate, which can be seen in Figure 2.3, which shows the side view of the substrate loading process.

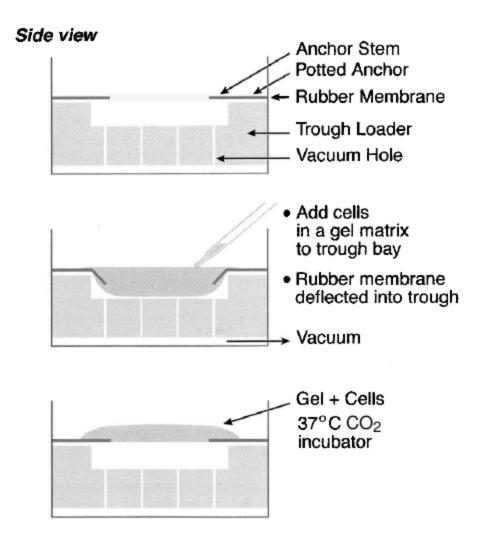


Figure 2.4: Flexcell side view of well with substrate (Garvin et al. 2003).

This design allows for uniaxial movement of the substrate, as well as equibiaxial strain applied to the substrate (Garvin et al, 2003). The archtangle shape of the rubber (rectangle with circular ends) allows for a nice long axis of a fixed substrate, and then provides the sides that are not fixed to move downward when the vacuum force is applied (Triantafillopoulos et al, 2004). Some limitations of this device are the overall uniform strain and the anisotropy of the strain profiles because cells and tissues do not always experience uniform strain (Brown, 2000).

The Flexcell device is in Professor Billiar's lab currently, and does a very good job of obtaining strip and equibiaxial stretch on a soft substrate. The main issue with this device is the inability to see the real time realignment of the cells in reaction to the stretch.

#### 2.2.3 Motor Driven Stretch Devices

Finally, this stretch can be achieved through biaxial strain which physically pulls the substrate outward in two directions perpendicular of one another (Brown, 2000). This can be seen in multiple devices. In a device that was used by Norton et al, a cross-shaped PTFE membrane was clamped to a two micrometer-driven mechanism with an accuracy of 5µm, to move the arms of the cross substrate and stretch the membrane (Norton et al, 1995). The topview of the device can be seen below in Figure 2.5.

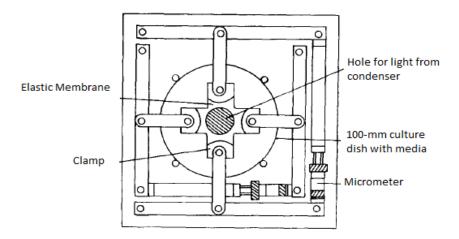


Figure 2.5: Micrometer-driven biaxial device (Adapted from Norton et al, 1995).

Other devices driven by motors include the Strex Cell Stretching System, which is made by B-Bridge International, Inc. The system can be seen below in Figure 2.6. It seeds cells on a soft-substrate silicon membrane that is shaped with a cross structure to reduce the Poisson effect created in biaxial stretch. The device allows for the biaxial stretch with a two motor system that

drive threaded rods to move the corners of the silicon well to stretch it, and the cells seeded on top of the well (B-Bridge International, Inc., 2009).



Figure 2.6: Strex Cell Stretch Device (B-Bridge International, Inc. 2009).

The motors are controlled by a computer driven program that will dictate the desired amount to stretch the membrane. To overcome the moment created by pull on any one axis, the arms are also attached to bearing slides that will keep the membrane in perpendicular alignment with the device and keep the stretch either strip biaxial if only one motor is running or equibiaxial if the two motors are running (B-Bridge International, Inc., 2009).

Various other devices currently used in labs are driven by four motors to obtain equibiaxial and strip biaxial strain. With each motor controlling one of the four axes of the membrane being stretched, it allows for more precise control to ensure the desired percentage of stretch is obtained at any given point in time due to the Poisson effect created. Currently in Professor Billiar's lab is a four motor driven device.

## 3 - Design Approach

Our goal is to design, build, and test a cell stretching device. The main objective was for the device to be able to stretch the cells biaxially, meaning both strip and equibiaxial stretching. The client expressed desire for the device to stretch the tissue for a minimum of six hours, but with the hopes of obtaining a stretching period closer to that of ten to twelve hours. Also, the device should be able to strain the tissue sample at a uniform rate between 2-20% at 1 Hz. The cells also need to be kept alive and at homeostatic conditions during testing.

After these objectives, other less critical, yet still important features were for the device to have a suitable cell viewing area of 1 square cm, have a stationary viewing area during the stretching process, and to have the device itself be inexpensive.

### 3.1 - Design Alternatives

In order to determine the client's priorities in the design, a pairwise comparison chart was constructed so that objectives could be ranked in importance. Once a final design was determined, this was a useful chart in the event that some aspects of the machine cannot be completed because the cost or time constraints cannot be met. From here, we were able to determine the top priorities of this machine should be stretching with strip and equibiaxial strain, keeping the cells alive during the stretch through the use of incubation, fitting the device under a microscope stage for real-time viewing of the cells during stretch, outfitting the device to stretch the pre-manufactured Strex well, and to keep the cost under \$1000. Secondary objectives include being able to adjust the strain rates and strain percentages of the motor and to minimize the cost. The smaller objectives are keeping the device inexpensive and maximizing the viewing area to see the cellular response of the stretch. The filled our Pairwise Comparison Chart from our advisor can be seen in Appendix A.

From the objectives, we were able to determine a list of functions the machine needed to perform in order to successfully meet the constraints and objectives of the project. From those functions, we were able to determine means to achieve them, which led into preliminary design options for our machine. In the functions means tree in Appendix B, we started with basic functions and brainstormed possible ways to make this happen based on previously reviewed stretching machines or other methods of stretching membranes that may not necessarily be associated with cells.

#### 3.1.1 - Four Motors System

A stretch system based off a clamped four motor platform has several advantages. This system can certainly achieve strip biaxial and equibiaxial strain and can do so without moving the center of the tissue sample in order to view its response to the stretch under the microscope. Each motor is responsible for moving one axis of the square well, and would be able to be controlled independently of each other. This allows for easy controlling for strip and equibiaxial strain. This type of system can also perform various strain percentages, strain rates, and duty cycles, though this is all based off the power, accuracy, and capabilities of the motor.

Another design option is a four motor system with linkages. This system is useful to create a stationary viewing area to see the cellular response to the stretch of the cells. Each motor is able to control the linkage movements and ensure strip and equibiaxial strain on the tissue.

This will be easy to add or remove strain on the cells and have a quick response to create variable duty cycles in the movement of the motors with proper programming.

A concern with a four motor system is the design space that is reduced in order to have enough room for all four motors to move and be properly vented to prevent overheating and keep within the size constraint of the microscope viewing area. This could be overcome by moving the

motors to a plane different from the membrane being stretched, however, this would involve more design of parts that would allow for the motors to be on a different plane that that of the stretched membrane. Each motor will also have to be equipped with a motor controller board to run it, which is expensive and could increase the cost of the device.

#### 3.1.2 - Two Motors Systems

The two motor threaded rod system is yet another design option. The two motors have an advantage over the four motor systems because they take up less room around the microscope stage; require less power than four motors do, and cost less money. Threaded rods are also an advantageous stretching mechanism because they are an easily obtainable system used in many applications. This aspect makes them easy to understand and work with because they are so commonly used. Using this method allows for both strip and equibiaxial stretching.

A commercially available version of this design is known as the Strex®. It involves two motors with threaded rods attached to guide rods that are connected to each corner of a square soft membrane. The bottom left corner remains stationary, but the other three corners move, stretching the plate in either one or two directions simultaneously. Reverse-engineering this is a design option, and a 3-D image model of it made in SolidWorks can be seen below.

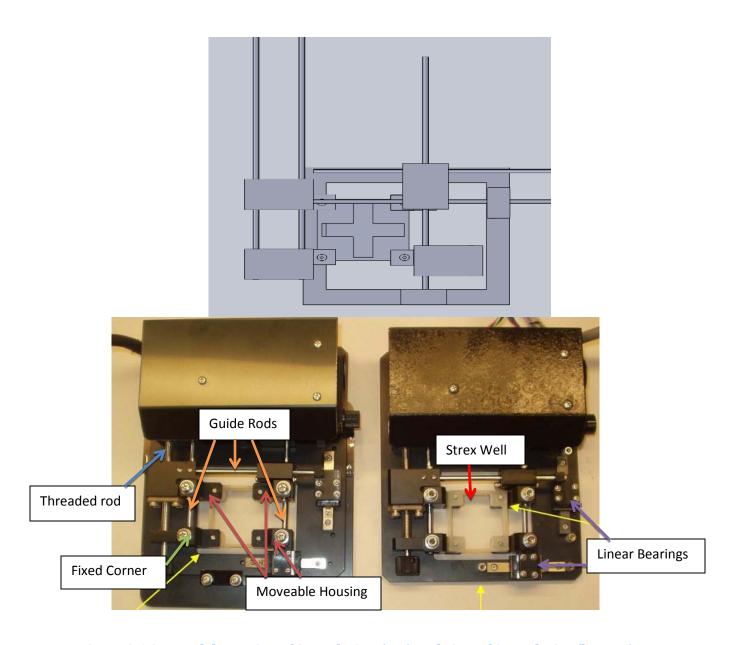


Figure 3.1 3-D Model Drawing of Strex design (top) and view of Strex device (bottom)

Each corner is attached to a moving carriage that is attached to a threaded rod driven by motor control and a guide rod that keeps the movement perfectly perpendicular to the threaded rod. To overcome the moment force created there are linear bearings attached to the rails which will allow for the motion to remain in the correct direction. The motor drives threaded rods to move the housings attached to each arm away from the original location of the Strex well it is attached to, and allows for the aligned movement of the well (B-Bridge International, 2010).

There are some disadvantages to this design, though. For example, the main concern with using the threaded rods to stretch the tissue sample is that the rods may not be able to move fast enough to strain at the desired varying duty cycles because of the threaded design. Another major concern is that when the tissue it being stretched, it does not stay centered over the viewing area, making it difficult to monitor cell response throughout the entire stretching process. One last issue is that two motors may not be able to generate enough power to move the machine at the desired rate.

Adding linkages is another variation to the two motor design. Two motors being the benefits of having a larger viewing area, a cheaper cost, and do not need as much power as four motors do to run the device. Linkages are also a useful tool for tissue stretching, being able to both biaxially stretch and provide means for varying duty cycles.

A major disadvantage of the linkage system, as touched upon before, is the possibility of unwanted stresses occurring on the edges where the linkages are attached to the tissue sample, which could potentially alter the collected data. Disadvantages of the two motor system include not having enough power to move the machine itself and not keeping the tissue in a centered position.

Another potential stretching design will resemble the four motor design however; the motors will be replaced by pistons driven by compressed air. A piston driven stretching device, much like the four motor design, will be able to achieve both strip biaxial and equibiaxial strain. An advantage to pistons is that there will be no need to learn how to then write complicated programs to run them, although the air pressure would still need to be controlled. There can be multiple compressors giving out different levels of compressed air in order to achieve both types

of strain. Compressed air can be easily dispersed at different rates to perform variable duty cycles, strain rates and strain levels.

Although there are many benefits to using pneumatic driven pistons, there are some drawbacks. Compressors can be big and bulky and can take up valuable space. There is limited space to deal with around a microscope and compressors could impede. Also, though the pistons would be faster than motors, in order to power the pistons air hoses need to be run to them which could become cumbersome when attempting to run experiments. The intended design of this stretch device with two sets of opposing motors may be large and may prove difficult to fit within the given space under the scope. An exterior structure may have to be built in order to support the system instead of mounting it directly to the microscope.

Using hydraulic motors to stretch the tissues is another design option. Hydraulics are useful because there is direct control of the motor via control valves, which can be useful to ensure pure strip biaxial and equibiaxial strain on the membrane. A noteworthy advantage of this option is the ability of hydraulic motors to apply significant power with very small tubes and hoses. This reduces the amount of parts around the microscope stage which can aid in the visibility of the tissue with the microscope.

This option, however, could prove difficult in programming the motors to work in synchronization, especially if four different hydraulic pumps are used. Another drawback could be in the variable duty cycles, because the pumps can quickly add or remove fluid to add or remove the force, but the acceleration of the piston at that speed may not be quick enough or slow enough for the needs of a fast or slow recovery in the stretch mechanisms.

#### 3.1.3 - Vacuum (Flexcell)

A vacuum based system works especially well for stretching cells equibiaxially. These systems are consistent, and exhibit good homogeneous stress on the cell samples. A system like the Flexcell would work very well, provide consistent data during testing, and culture plates would be easily available. The Flexcell device is existent in the lab at WPI, but its main issue is the inability to view the real time reaction to the cell stretch because of the closed top design of the vacuum chamber. Attempts have been made to modify the device however the results have not been able to view the results in real time. However, this device would not need any building or programming which can save large amounts of time in the design process. Ultimately this system would allow for more time testing samples and obtaining data.

#### 3.1.4 - One motor system

This design consists of one motor which applies downward pressure to the tissue sample to stretch it over stationary posts. Stretching the tissue over stationary posts will achieve strain over the tissue. One of the advantages of this design is first and foremost its simplicity. There is only one motor which moves upwards and downwards and can be programmed to achieve variable duty cycles and strain rates. The motor pushing the tissue is the only moving part of this design which also contributes to its simplicity. This design will be a relatively small structure so it will not be difficult to fit under the microscope stage, but will be difficult to find a way to view the cells during the stretch. Also due to the one motor, both strip biaxial and equibiaxial strain is more difficult to achieve, and would most likely require some movement of parts to make sure. Like with all motors, a drawback is that programs must be written in order for them to function correctly; these complicated programs will have to be learned and executed.

Another one motor design angles the motor at 45° in the upper right corner of the Strex well to obtain equibiaxial strain. Linear bearings are attached to guide rods which are then

attached to the four corners of the Strex plate. The motor attaches to an arm and flywheel and motor which then causes the movement of the upper right corner linear bearings which translate the motion to the other two corners of the membrane. Strip biaxial strain is attained by adjusting the angle at which the motor pulls to as close to zero as possible, and then accounting for some error in the Poisson effect which will be seen. This can be seen below in Figure 3.2.

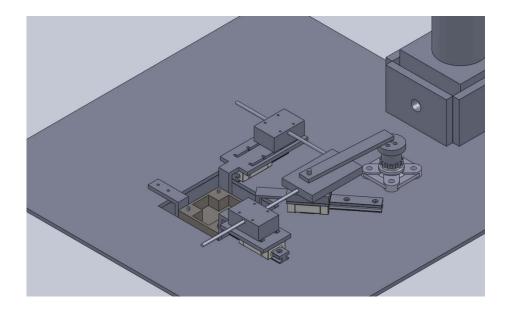


Figure 3.2: Linear Bearings One Motor design (equibiaxial set up)

This device allows for no programming, and the parts are easily available off the shelf. However, the adjustment of the motor and linear bearing system will leave room for error, so the precise angles at which the device needs to be to obtain specific strains will need to be very accurate.

## 3.2 - Design Process

In order to determine what design would meet the most objectives desired by the client and fulfill all the constraints, we developed a function and means chart to determine what ways these various objectives could be obtained in our designs.

We developed this chart, shown in Appendix B, in order to help give us a better visual idea of which design would be the most efficient at completing our intended tasks. On the Y-axis, we had both our constraints and objectives listed. On the X-axis, we listed our potential design ideas. Where the design ideas meet with the individual constraints is where we assigned values. These assigned values are on a scale of 0-100 with 0 meaning the design will not be able to perform the constraint/objective and 100 meaning the design will definitely be able to complete the constraint/objective. We assigned these values based upon our personal opinions gained through research on whether or not the constraints and objectives could be achieved.

After assigning values, we tallied the total points up per design and ranked them accordingly.

The two main designs that are being considered are reverse engineering the Strex design, and the one motor linear bearing design. A comparison of the two designs can be seen below in Table 3.1.

Table 3.1 Comparison of Strex and One motor design

Strex Design Reverse Engineer	<u>Functions</u>	One motor design
Capable of uniaxial stretch	-Uniaxial stretch	Capable of uniaxial stretch
Capable of equibiaxial stretch	-Equibiaxial stretch	Capable of equibiaxial stretch
The Strex based design would be	-Strain rates	This design would allow for variable
capable of achieving 2-20% strain		strain %, however it would most
with any strain % in-between.		likely be capable of a particular set
		of strain percentages:
		2,5,10,15,20% strain
Adjusting frequency will not be an	-Frequency	Adjusting frequency and varying

issue with this device, frequencies of (.5, 1, 1.5, and 2) Hz can easily be achieved. Also the frequency is not limited to these numbers it can go anywhere between and above.  This device will be capable of varying duty cycles.  These devices will both fit on the Zeiss stage this design will be slightly larger.  Compatible with Strex silicon well.	-Duty cycles  Size	frequency is possible, however the device will most likely be limited to several predetermined frequency setting that would be dependent of power source/ current.  Varying duty cycles with this device will require modifications; the user would have to change an electrical component (diode).  Slightly smaller/lighter than Strex design.  Compatible with Strex silicon well.
Companiole with Silex shicon well.	Well type	Companiole with Silex sincon wen.
Will require time consuming programming. Motor control boards or stepper motor drivers.  Programming is difficult to learn and small mistakes in program can cause un-noticed/unwanted results.	Programming	None is required. This system relies completely on mechanics and current for motion.
Will use stepper motors, however, the other components would most likely be custom made with some being premade. The fabrication will be more difficult in comparison to a model with mostly premade parts. Also the device will require a computer with expensive motor control boards to run the device.  Also the motors will have to change direction often which results in greater power consumption which may cause overheating. (High stress on motors)	Parts	Most parts will be premade from a distributor such as PIC -Design or small parts. These components are reliable, available in various sizes and finishes. Some components may be customized, however these components will include mostly simple parts which are cheaper to make from free scraps or cheap aluminum.  The motor in this system will turn in one direction, allowing for a smooth stretch and minimal power consumption. (Should be minimal stress on motor)
More intricate design, with more components. Both systems require near perfect assembly, however, with more components it increases difficulty of having whole system in sync. Though, if the device is made properly is should allow for smooth stretch with minimal friction.	Mechanics/Dynamics	With fewer components this device is dynamically and mechanically simpler. However, like the Strex design to prevent friction/jamming, the components need to be assembled with no error. Also with the motor never having to change direction it may provide a smoother more homogeneous stretch.
Greater cost due to the second motor, more components and more intricate mechanics. Also the motor control boards can become quite expensive if not available. Stepper motor drivers are much cheaper however adds cost as well.	Cost	Lower cost due to minimal components and no motor controlling hardware.
This system will allow for simple use once ready. The user can simply	<u>User Friendliness</u>	This device requires the user to manually adjust the device to change

command the device to perform	strain type, strain %, and duty
whichever stretch it wants and can	cycles. The user would carefully
run it.	have to read the manual to ensure
	the device was appropriately
	adjusted. (If user does not pay
	attention to detail it is possible for
	miss-setup)

## 3.3- Final Design

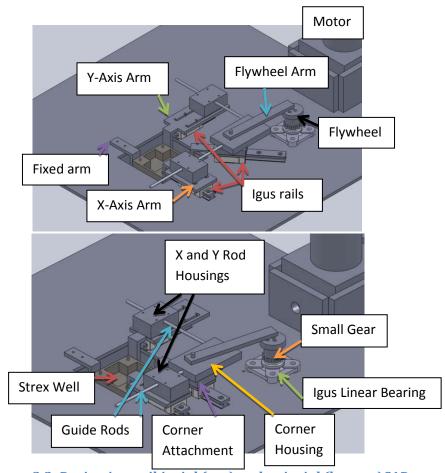


Figure 3.2: Design in equibiaxial (top) and uniaxial (bottom)CAD

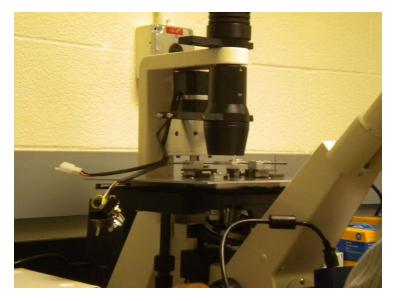




Figure 3.3: Our device on the Zeiss microscope (top) and close up with cells seeded on the well

## 3.4 - Device Description

Through careful deliberation, our team decided that a one-motor concept for a stretch device would give us the best opportunity to meet all objectives taken from our client statement. Having chosen the general design, we began to search for materials and parts that could be used to achieve major aspects of the device: the device must be powered by a single motor and must stretch a Strex silicone well both uniaxially and equibiaxially.

#### 3.4.1 - Motor Selection

Calculations for determining the force, torque and power needed to stretch the wells were accomplished for initial considerations regarding choices of motors. For this design a synchronous motor suits the device best, having a system that limits the stretch to its preconfigured state the motor simply has to turn at the specified frequency in a single direction. This is why stepper motors were not considered; they do offer precise motion due to the fact that as the drivetrain turns it does so in steps not in a smooth/fluid motion. The material properties of the ST-190-XY well from B-Bridge were not available in the specifications, however, the well itself is made from silicone elastomer comprised mainly of polydiethylsiloxane (B-Bridge International). In order to determine certain specifications needed for our motor, we set up an experiment on an Instron machine to stretch the Strex well and find its material properties. We know from calculations prior to the test that the max distance needed in travel at maximum strain (20%) for the well would be 6 mm or 0.006m. We programmed the Instron machine to stretch to this distance and record all exerted forces. At the completion of the test we found that the maximum force produced was approximately 19 Newtons. Using this number, along with the maximum distance needed in stretch for at maximum strain, we calculated that needed torque would equal 0.114 Nm. However, this calculation is solely for uniaxial stretch as the Instron machine could only pull in one axis. Therefore, we multiplied our initial uniaxial torque value by two in order to account for the two axes of stretching. This new value of 0.228 Nm would be used as the value or torque necessary to achieve stretch in the equibiaxial position. These two values depict bare minimum values of torque needed to successfully complete uniaxial and equibiaxial stretching of the Strex well. We decided to increase they torque due to certain aspects of our device's potential usage. Our device may potentially undergo experiments where it is continuously running for long periods of time which would place a large strain on the motor.

Extra torque would be beneficial in ensuring the motor is pulling at a steady pace for the extended period of the test. Also concerning to our group was the fact that we are utilizing a 1:3 gear ratio in order to increase speed to meet one of our objectives. However, using a 1:3 gear ratio exhibits its 1/3 fractional percentage on the motor's torque as well. To conclude the original torque of our chosen motor, when divided by 3, must be well over the 0.228 Nm torque needed for equibiaxial stretch to achieve this goal, as well as, to provide a factor of safety for the overall effect of friction throughout the device.

Taking all of these factors into consideration, our team was able to decide upon a motor that would be efficient for achieving all specifications necessary for our device as described above. The motor chosen is a brushless 12V DC Synchronous Gear Motor manufactured by Wonder Motor. Unlike stepper motors which provide movement in choppy steps, synchronous motors provide smooth operation. This will reduce vibrations which could disturb the seeded cells on the Strex well. This particular motor has a maximum output of 38 RPM while providing 8.4 Nm of torque. As stated above, the chosen motor's torque, when divided by 3, must well exceed the 0.228 Nm needed for equibiaxial stretch. If you take 8.4 Nm and divide by 3, you are still left with 2.8 Nm of torque. This value well exceeds 0.228 Nm which is excellent seeing as though we also need to account for all friction experienced throughout the device. Overall, the chosen motor will effectively stretch the well to our certain specifications and provide this stretch for extended periods of time without over exerting itself and failing.

To place this motor in the appropriate position on the device, it was necessary to devise a bracket system to contain the motor. The mounting bracket was designed in order to utilize the motor in an inverted position. This inversion allows us to have a direct belt drive design rather than run the belt off of numerous gears which would increase the net friction in the system.

### 3.4.2 - Gear and Pulley System

We are utilizing a 2:1 gear and pulley system to achieve desired rates of stretch as described in our project's objectives. Gears are necessary because we chose our motor based upon specifications needed for torque; rate of stretch was a secondary concern for us. The motor has a maximum speed of 38 RPM's; this equates to approximately 0.633 Hz. Referencing our objectives it is clear that we are aiming for a rate of stretch which can vary from anywhere between 1-2 Hz. In order to achieve the maximum rate of 2 Hz, we must utilize a 1:3 gear ratio. This 1:3 ratio essentially means that the gear coming off of the motor shaft will be three times the diameter and have three times as many gear teeth than the gear which will help directly stretch the well. So, by the time the gear on the motor completes one revolution, the gear directing stretch will have completed three revolutions which is where the increase of speed originates. The gears we are using are manufactured by W.M. Berg. Respectively on the motor and directing the stretch the diameters measure 47.24 mm with 36 gear teeth and 16.8 mm with 12 gar teeth. These are 3-d pulleys which entail they are turned by a belt with teeth that interlock with the teeth coming off of the gears. The belt being used to conjoin these two gears is a 3-d pulley also from W.M. Berg.

The gear directly causing stretch will do so by turning a crankshaft which will be placed within the inner bore of the gear. This assembly of gear and crankshaft will sit within a custom made bearing with the gear sitting above the top of the bearing to ensure it will spin freely and not be impeded. The crankshaft has a one inch diameter flat top surface with three pre-drilled holes. These holes are off center by 3mm, 4.5 mm, and 6 mm representing varying strain percents of 10, 15, and 20 percent. Depending on the type of stretch being tested, the location of the crankshaft will change. To achieve stretch, an arm will be positioned over one of these holes

and secured; the other end of the arm will be anchored to a housing which will be explained in the following section.

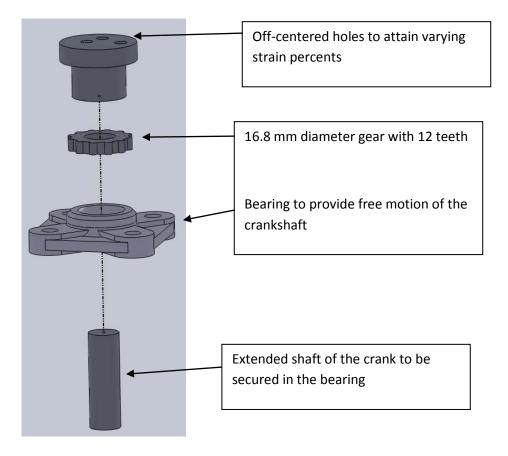


Figure 3.4 -Exploded View of Crank Shaft

# 3.4.3 - Linear Guide System to Directly Stretch Strex Well

This stretch device must achieve both uniaxial, as well as, equibiaxial stretch on a silicone Strex well. To ensure this stretch occurs as smoothly and linearly as physically possible, a rail system had to be installed to counter the Poisson Effect from the silicone material.

The platform which our entire device will sit upon is a 12"x12" piece of aluminum with a 6mmx6mm square milled out in the center for viewing of cells. Because our chosen microscope came equipped with a movable X-Y stage which added height, we also needed to install an adapter to clear this distance. The adapter plate was designed with two functional aspects in mind, one was to allow the device to be compatible to the microscope stage and second to allow the plate to clear the vertical distance between the of the movable X-Y stage equipped to the microscope platform. The adapter plate was designed to fit snugly in the X-Y stage, which is why no fasteners are necessary to secure the device in place. One can simply insert and remove the device with little trouble. To provide for smoother travel and to counter the moment the motor causes in the rear of the device, acrylic spacers were made. The acrylic spacers have a height of 9/16" this is because the adapter itself is 1/2" in height, though it is sitting on a 1/16" lip. The combination of both the adapter and spacers provide a level foundation for the device and allow smooth motion due to positive weight distribution. Also the acrylic can slide on the stage without scratching the coated surface of the microscope platform. Dimensions of both the adapter and the spacers can be viewed in Appendix H.

For stretch in both the X and Y axis linear rails were implemented to keep forces exerted in one direction with only negligible tolerance in the axis of stretch. A corner rail is used to apply direct strain to the upper right corner of the Strex well; stretch will be initiated from that corner. However, the type of stretch will depend on the angle of the rail. For uniaxial stretch, the rail acting upon the top right corner of the well will be parallel to the X-axis rail and be aligned so that the back edges of both the rail and Strex well line up. As the motor applies a force through the crankshaft, the Strex well will only deform in the X-axis through movement along only two rails. Conversely, if equibiaxial stretch is being tested, the rail will shift up 45 degrees towards

the Y-axis. The end of the rail will line up so it is centered with the top right corner of the Strex well. In this stretch set-up, when the motor applies force through the crankshaft, that force is applied at a 45 degree angle through the top right corner of the well. Exerting force at a 45 degree angle will evenly distribute along both the X and Y axis and be kept linear by the rigid rails. The rails used in our device are manufactured by the company Igus. They are composed of 6063-T6 Aluminum. The main reason for choosing these particular rails had to do with their extremely low profile in comparison to other similar products. They stand a mere centimeter tall which helped our team to decreases overall height of the device to fit well underneath our microscope.

Sitting directly atop the rails are cast zinc chromated carriages which will slide along the rail to provide a smooth means of stretch. Unlike most bearings, these carriages do not glide using a ball bearing system. Instead, the inner surface which interacts with the aluminum rail is coated with a wear-resistant iglide® J plastic sliding pad. This custom plastic created by Igus is self lubricating which does away with the need for ball bearings or constant application of lubrication to avoid binding. Each carriage has four screw holes for attachment to other pieces.

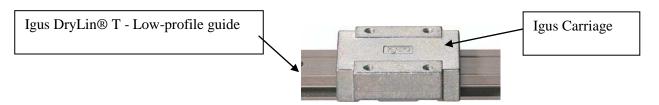


Figure 3.5 -Igus Linear Rail and Carriage System

The next component of the linear guide system which is being placed upon each carriage is an arm that will create a bracket to hold the Strex well in place close to the objective lens.

Each arm is made out of 1/8" oil hardened steel and has been cut and welded into a precise

position to hold the well perfectly. What is particularly innovative about both the X and Y axis arms is their multi-positional design. Each arm has channels and locking slots enabling the user a more simplistic means to change the stretch position between uniaxial and equibiaxial stretch. Finally, there is a stationary arm anchored to the device platform which will remain rigid to secure the bottom left corner of the Strex well.

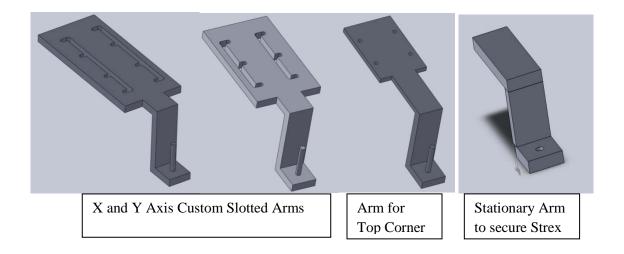


Figure 3.6 - Four Stretch Arms

The final component of the linear guide system consists of pieces which are the same for the X and Y axis rails, but slightly modified for the rail acting upon the top right corner of the well. For the X and Y axis the last component is simply a 3mmx2mm piece of ½" aluminum. The piece has four threaded holes which line up directly with the four holes on the carriage. The threads from this housing in combination with the carriage holes will create a very tight connection to reduce vibrations during operation. In addition to these four holes, there is also a clear hole through the 3mm long face of the housing. The clear hole will be 1/8" in diameter and sit in the center of that face; the center of the hole be 1/8" from the top edge. However, the housing sitting atop the arm on the rail applying force to the top corner of the well is a bit

different and consists of a two part system. The first part will consist of a 2.5cmx2cm piece of aluminum and have four screw holes that match up with those in the carriage, however, this piece will be only ¼" thick and have one additional hole. The additional hole will be centered 7.5 mm in from the 2cm side and 1 cm in from the 2.5 cm side. A screw will protrude through this hole from the bottom and be countersunk to maintain a level surface. The protruding screw is where the second part of this housing will be attached. The second part of the housing is a 3cmx5cm piece of ¼" aluminum. There will be a threaded hole 15mm in from both the 3cm side and 5cm side that will attach both pieces.

In order to further increase the linearity of the forces being distributed in each axis, stainless steel rod will be inserted into the 1/8" clear holes in the sides of each of the three housings.

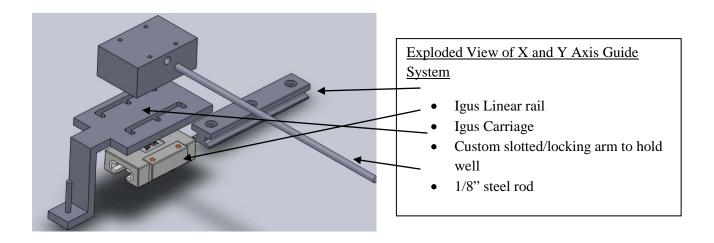


Figure 3.7 - Exploded View of X and Y Guided System

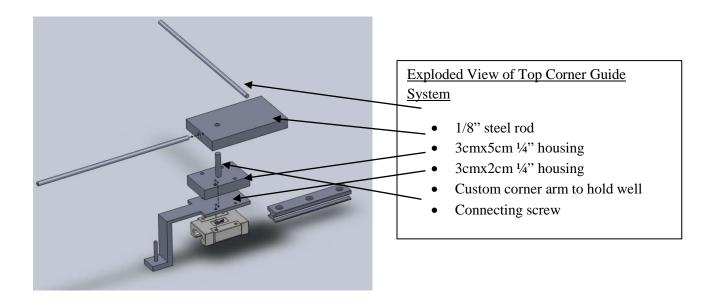


Figure 3.8 - Exploded View of Corner Assembly

# 3.4 - Microscopy Methods Currently Used

In order to choose a suitable microscope to use in combination with our device, our group first decided on objectives and constraints needed in a microscope for it to be considered sufficient. There are many microscopes at our team's disposal, however we needed to carefully research each one in order to find the best fit.

# **Objectives**

Although not quite as important to achieve as constraints, objectives (if met) will make our device much more professional and successful. The device should move fluidly around the platform of our chosen microscope to provide a smooth transition between viewing cells. To achieve that our group wants to keep the device as light as possible to keep torque down which would increase friction with the surface. Also the device was designed with protecting the microscope in mind, considerations are included in the Device Description.

### **Constraints**

One of our main concerns was the ability to view cells through the Strex well. We could not successfully complete our project if the cells that we are stretching could not be viewed. Due to its importance, the ability to view cells through the Strex well fell under constraints. Another concern of the group was the dimensions of available microscopes. We want to view the cells being stretched in real time; therefore our device has to fit within the dimensions of a microscope's platform. The most important dimension needed in a proper microscope is the distance been the objective and the seeded cells on the Strex device. Through trial and error, we found that a 60X objective lens would be necessary to view cell movements such as cytoskeleton realignment. With that in mind, it is key to know that as the magnification strength increases in an objective lens the closer it needs to be to what it is viewing in order to be effective.

In order to precisely move the stretched Strex well around to view individual cells, a microscope with this ability was necessary. Given our constrained supplies, the group decided a movable X-Y platform would suffice for Strex well movement. However, movable X-Y plates take away from the overall height that our device could be so the group had to take that into consideration upon designing our project.

### **Best-Fit Microscope**

Through careful examination of available microscopes and much deliberation, the group was able to decide upon one microscope which would be utilized for our project. That microscope is the Axiovert CFL40. This is an inverted microscope which allows us more height to design our device with. Also, because it is inverted the light source comes from above while the objective lens comes from below. This allows sufficient light for viewing and the objective lens is able to move quite close to the Strex well to compensate for the increased magnification

of the 60X lens. The dimensions of the microscope platform allow ample space for our device to be anchored upon. A movable X-Y plate is already installed on this microscope which is beneficial, however, our group will need to design around the fact that it decreases the available height between the top of our device and the light source.

# 3.5 - Incubation Methods Currently Used

Systems are needed to control isolated environments to maintain specific conditions for cells. Generally speaking these conditions are the same; maintaining appropriate temperature,  $CO_2 \& O_2$  levels, humidity and pH. These systems can be applied to microscopy by using microscope cage incubators, which have been shown to work as well as regular bench-top incubators. Most of these cage incubators use gentle streams of warm air to control the temperature. To monitor the temperature most systems use a thermocouple is inserted into a reference well to control the temperature as close as possible to the sample. Most systems aim to ensure a specimen temperature stability of  $\pm 0.1 \,^{\circ}$ C.

These cage incubators can be designed for virtually any inverted or upright microscope that is available. These models are generally customizable for specific needs and include various chambers and interchangeable plate adapters which allow for the use of any cell culture support (petri, glass slides, multiwell plates, and others). These cages also are compatible with manual and digital  $CO_2$  /  $O_2$  gas controllers. pH is generally controlled with buffers however, humidity is important to ensure water from the cell bath does not evaporate out of solution increasing the concentrations of potentially harmful chemicals. All these systems generally function at similar accuracies and specifications.

The amount of time the cells need to survive is related to the type of accuracy in a system one would look for. The longer a test needs to be run, the higher accuracy one needs to maintain appropriate temperatures for their samples.

# 4 – Device Validation

In order to ensure the device could work properly, validation had to be conducted on mechanical components as well as the ability of the device to view cells under the microscope during stretch. High density mapping was used to map strain fields in the well, and other tests were performed to validate the mechanical function of the device.

# 4.1 Test Trials - High Density Mapping

Precision and validation are key components to a successful project. Although the Strex manual states that the seeded cells are subject to the same strain as the well is being pulled to (B-Bridge International, 2009), we wanted to personally validate this fact on our own to confirm uniform strain across the well. To give our device supporting evidence, our group decided to use high density mapping, or HDM, as a validation tool. HDM is a technique that uses pixel displacement after a sample it stretched to map its strain (Kelly D et al., 2007). Kelly, along with other researchers, developed this technique to observe a heart over the course of a heartbeat. A camera capturing the sample's movement is connected to a computer, where the data is processed. Once the test is complete, the user can review the pictures side by side to see how far each area moved between frames. This information is exported to an excel spreadsheet to calculate the exact values (Kelly D et al, 2007). This tool is very useful for mapping strain at various points on the sample, to ensure the correct strain value is being used. This is exactly what we used the program for.

To validate the strain field, the Strex well was stretched to 20% strain. This value was chosen because it is the highest strain our device will stretch the well to. The area that was

focused on was a 20 mm<sup>2</sup> region, because this is where the cells will lie during stretching. This region is shown in the figure below.

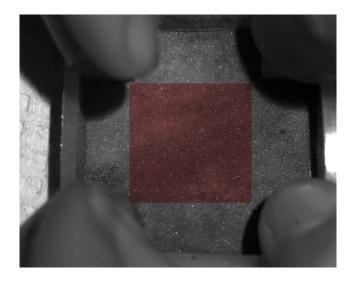


Figure 4.1: Hand stretched Strex well to 20%

As a reference, there are 31 pixels per mm. The HDM program was set to record 16 pixel shifts for 32 x 32 pixel boxes. After the high speed camera, a Photron Fastcam 1280 PCI, captured the well being pulled to this strain, the images were compiled using MATLAB. The data was also filtered using a Gaussian filter by inputting the function *fspecial('gaussian', hsize, sigma)*. The code for the compiling function is included in Appendix I. From there, the matrices for both the u and v displacement were totaled using the function *utotal=sum(udata\_s, 3)*. The s signifies that the data has been smoothed. Once the matrices were condensed, the strain was able to be calculated using the displacement of each point.

To calculate the strain, the displacement of one point was subtracted from the following point, and the total was divided by the pixel shift used. For example,  $(X_{1,2}-X_{1,1})/16$  would be the formula for the first strain point in the X direction.  $(Y_{2,1}-Y_{1,1})/16$  would be the formula for the first strain point in the Y direction.  $\{[(X_{2,1}-X_{1,1})/16]+[(Y_{1,2}-Y_{1,1})/16]\}/2$  would be the formula

for the first shear strain point. These calculations produced another set of matrices. MATLAB was then used to create a contour plot. The figures below show the X axis strain, the Y axis strain, and the shear strain respectively.

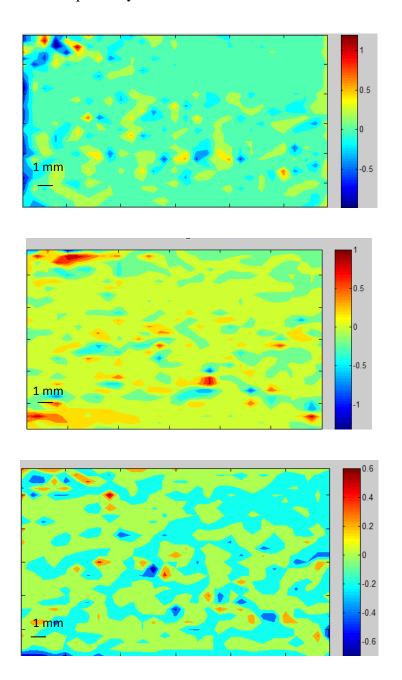


Figure 4.2: HDM results in x axis (top), y axis (middle) and shear strain (bottom)

The green color represents 20% strain, the warmer colors represent a higher strain, and the cooler colors represent a lower strain. As depicted in the pictures, the strain is fairly uniform across the selected section of the well. The abnormal pockets are most likely due to the method used to stretch the well. In the future, these strain fields can be used as a template to compare our device's strain percentages. This will ensure the wells are being homogeneously stretched to the correct strain.

### 4.2 Device Validation

• Arms reached well (well fit on device)

As seen earlier, the Strex well has holes 30mm apart that allow the device to mount to a stretch device. The arms of our device were obviously designed to work within this constraint. In both Equibiaxial and Uniaxial positions the arms reached the appropriate distance to allow the standoffs to slide into place.

### Speed Change

Using a stop watch and counting rotations change in rpms were validated. The device can accurately stretch between 0 and 1.9 Hz with increments of 0.5 Hz or 30 rpm.

### • Both strain types work

The device was simply tested in both equibiaxial and uniaxial setting, and in both cases the device functioned as designed. The

Can change % strain
 By utilizing the three off-centered holes on the top face of the crankshaft, strain
 percentages exhibited on the well can vary between 10, 15, and 20 percent.

### Stretched well

- The device was capable of stretching the well in both settings and the steel rods or linear bearings did not bind during the stretch.
- Device compatibility with Zeiss Microscope:

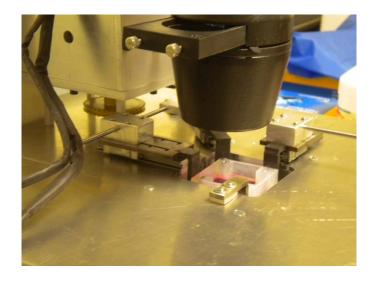


Figure 4.3: Stretch Device on Zeiss Microscope

The device was outfitted with an adapter below the 1/8" plate, the adapter had dimensions based of the X-Y stage of 5.25" x 3.5". The adapter was checked to ensure it would mount securely in place, then clearance was checked to ensure it cleared the back of the moving piece. After 6 cm<sup>2</sup> viewing area was milled in the center to allow the light to come down through the well. The hole also allows the arms to come down and allows enough space for them to move.

## Motion on scope:

Motion of the device and X-Y stage was simply validated by placing the entire device on the stage and checking the weight would not exceed the X-Y stages weight limit. Also in the rear of the device acrylic spacers were used to support the weight of the motor and avoid scratching the stage.

### o Viewing cells:

The arms on the linear guides drop to ensure the cells are close enough to the objective. The device was placed on the scope and the distance to the objective was checked. The objective could come up all the way to the bottom of the well using the coarse adjustment, using the fine adjustment the cells can be viewed in focus. A picture of the seeded cells was taken below on the Zeiss microscope to validate they in fact can be seen.

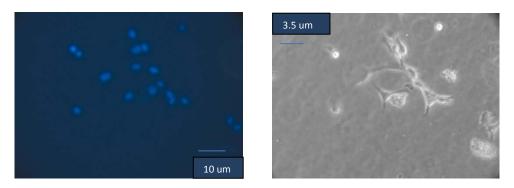


Figure 4.4: Rat Aorta Cells at 40X magnification with Hoescht and Phase Views

Additional Photos can be found in Appendix P.

### • Heat Test:

o The motor was ran for an extended period of time (4 hrs) at room temperature and the temperature was taken every half hour. The results of the test show the temperature never reached a level that can damage the motor or ruin any data.

Table 4.1: Heat Test Results

Heat Test Results		
Time (hrs)	Temperature (Celcius)	
0	25	
.5	26	
1	25	
1.5	27	
2	31	
2.5	32	
3	33	
3.5	33	
4	34	

Graphical analysis of the heat test shows a steady but slow increase in temperature. From (Figure x) using a best fit line the change in temperature attained was 2.5 degrees/hour, therefore for every hour the temperature increased only 2.5 degrees. If the test would have ran for another 4 hours the expected temperature would have been 45 degrees. This temperature is not very high nor is it anywhere near a temperature that could damage the motor or cause complications.

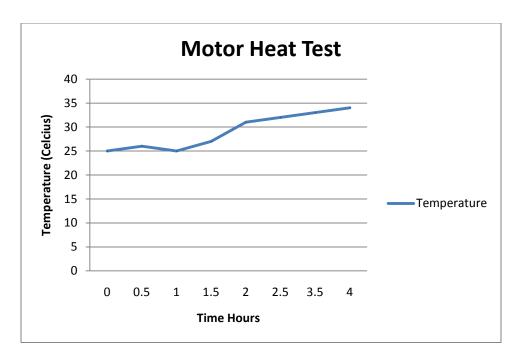


Figure 4.4: Graphical Results of Heat Test Over Time

# 5 - Conclusions and Recommendations

### 5.1 Conclusions

Our completed device was able to meet our constraints of achieving strip and equibiaxial stretch, however, the corner linear bearing needs to be moved to achieve this. The device can stretch at different rates by adjusting which hole on the flywheel the arm is attached to by simply changing the distance from the corner housing screw to the crank shaft hole or changing the hole distances on the arm. This change is described in Appendix E (Device General Use).

The device can be equipped to the Zeiss Axiovert CFL40 microscope and is capable of viewing cells during stretch, while placed on the microscope stage. The device fits in an incubator that fits around the Zeiss microscope to keep the cells alive during stretch. It can run for extended periods of time and the motor does not cause a change in temperature that will compromise the cells. The device is able to stretch cells at 1.9Hz at percent strains of ten, fifteen, and twenty percent strain.

The overall cost of the device was around \$880, including the cost of shipping of various parts. This came in well under the \$1000 budget. A large component of this cost was the \$250 spent to purchase Strex wells. The wells had to be purchased in bulk from B-Bridge, but this will be useful for many future tests. If the initial cost of the Strex wells were removed, the cost of the device would be \$630, which is significantly less than the \$1000 budget. A complete budget breakdown can be seen in Appendix C.

### 5.2 Recommendations

First, we would suggest allowing the device to include variable duty cycles. This aspect would allow the user to study how cells react to variable loading and unloading rates. A resistor

can be used to change the speed of the motor to attain these variable duty cycles. Because the body is subject to many non-uniform stretching rates, observing cell reaction during these variable cycles can be extremely useful.

We also recommend making the device more compact. Currently, our device fits well on the microscope stage and allows for motion on the X-Y stage. However, a smaller and lighter device would allow for a much more marketable device. The majority of the device's bulk is derived from the motor. Using a smaller/lighter motor would decrease the device's weight significantly. Also using different materials could significantly lower the device's mass, many strong but light plastics are available for these types of applications.

As discussed before, our device can be modified to stretch cells either uniaxially or equibiaxially. This transition between strain types requires the user to follow several steps to before successfully changing the device setting (full directions can be found in Appendix E). In the future, we suggest creating an easier and more user friendly transition. For example, creating a sliding arm mechanism with a locking pin could accomplish this. This would allow for a shorter setup time for experiments and a reduction in the chance for error in the process.

Creating an easier transition between strain percentages would also be a useful aspect.

Currently, there are multiple arms used for this transition (full directions can be found in the User Manual in Appendix E). Creating one arm that can transition between strain percentages would be much simpler. This can be accomplished by using Heim joints and a turnbuckle. This mechanism is shown in the figure below. Also by being able to lengthen the arm, the device could potentially be used to compress the wells.



Figure 5.1: Heim Joints (Lunsford, 2010).

Heim joints allow for smoother stretch, because they are designed for dynamic applications. The turnbuckle has opposing threads, so it can either bring rods closer together or farther apart. These threaded rods can be marked to indicate where the correct distance for each strain percentage is located. This design would allow for a much more user friendly transition.

Linear bearings on the inside of the housing on the X and Y linear guides would be a smart design alteration. They allow for more fluid motion and reduce the amount of friction in the system. Over time, this will be a better design because friction may cause the rods to bend, bind and possibly lead to other mechanical problems. Some linear bearings available through Igus can be seen below in Figure 5.2. The linear bearings come with plastic inserts to allow them to be run without lubrication, and will remove the metal on metal contact in the design. Another potential modification to minimize friction and increase fluidity of stretch would be to ball bearings at the point of contact between the arm and the crankshaft, as well as the arm and the corner housing.



Figure 5.2: Igus plain linear bearing (left) and Igus enclosed linear bearing (right) (Igus, 2010).

Another recommendation would be to find an easier method of determining what speed the motor is running at. Currently, the only way to determine this is through counting the number of rotations experienced over the course of a period of time. The best way to improve this issue would be through a tachometer, which can measure the revolutions per minute of the crank shaft. This would be an inexpensive purchase, but would just have to fit in the design space currently available.

One final recommendation is to use stainless steel arms instead of oil hardened arms.

Stainless steel rods have a higher wear resistance and are already heat treated, the oil hardened arms are much more brittle and often difficult to work with. They are also become very brittle when welded, unless allowed to anneal properly. Some of our steel arms broke due to cooling to quickly when being TIG welded.

Overall these improvements would increase the user-friendliness of our device and would reduce the time to change between strip and equibiaxial strain. This device is an innovative approach to stretching cells and with these improvements can reduce time spent in the lab setting up experiments.

### References

- Ananthakrishnan R, Ehrlicher A. (2007). The forces behind cell movement. Int J Biol Sci 3(5):303-317.
- B-Bridge International, Inc. (2009). *Life Sciences Products by STREX*. http://lifesciences.b-bridge.com/products/supplier/STREX/.
- Balaban NQ, Schwarz US, Riveline D, Goichberg P, Tzur G, Sabanay I, Mahalu D, Safran S, Bershadsky A, Addadi L (2001). Force and focal adhesion assembly: a close relationship studied using elastic micropatterned substrates. Nat Cell Biol 3(5):466-472.
- Bieler, F.H.; Ott, C.E.; Thompson, M.S.; Seidel, R.; Ahrens, S; Epauri, D.R.; Wilkening, U.; Schaser, K.D.; Mundlos, S.; Duda, G.N. (2009). Biaxial cell stimulation: a mechanical validation. *Journal of Biomechanics*, 42(11). doi:10.1016/j.biomech.2009.04.013.
- Brown, Thomas D. (2000). Techniques for mechanical stimulation of cells in vitro: a review. *Journal of Biomechanics*, *33*(1). Doi: 10.10/16/S0021-9290(99)00177-3.
- Chen, C. S., Mrksich, M., Huang, S., Whitesides, G.M., Ingber, D.E. (1997). Geometric control of cell life and death. *Science* 276. doi: 276:1425–1428
- Flexcell International Corporation, Initials. (2004). *Flexcell international corporation*. Retrieved from <a href="http://www.flexcellint.com/">http://www.flexcellint.com/</a>.
- Fredberg, U., Stengaard-Pedersen K. (2008). Chronic tendinopathy tissue pathology, pain mechanisms, and etiology with a special focus on inflammation. *Scandinavian Journal of Medical Science Sports* 18(1):doi: 18(1):3-15.
- Garvin, Joanne; Qi, Jie; Maloney, Melissa; Banes, Albert, PhD. (2003). Novel system for engineering bioartificial tendons and application of mechanical load. *Tissue Engineering* 9(5).
- Hasel, Cornelia, Dürr, Suzanne, Brüderlein, Silke, Melzner, Ingo, and Möller, Peter (2002). "A cell culture system for long-term maintenance of elevated hydrostatic pressure with the option of additional tension" <u>Journal of Biomechanics</u> 35:579-584.
- He Y, Macarak EJ, Korostoff JM, Howard PS. (2004). Compression and tension: differential effects on matrix accumulation by periodontal ligament fibroblasts in vitro. Connect Tissue Res 45(1):28-39.
- Hsu, Hui-Ju, Chin-Fu Lee, Andrea Locke, Susan Q. Vanderzyl, and Roland Kaunas. (2010). "Stretch-Induced Stress Fiber Remodeling and the Activations of JNK and ERK Depend on Mechanical Strain Rate, but Not FAK." *Plos One* 5.8 Print.

- Igus, (2010).Plastics for Longer Life. http://igus.com/wpck/default.aspx?PageNr=3586&CL=US-en&scroll=367&step=5
- Juliano RL, Haskill S. (1993). Signal transduction from the extracellular matrix. *Journal of Cell Biology* 120(3):577-585.
- Kelly DJ, Azelogu EU, Kochupura PV, Sharma GS, Gaudette GR (2010). Accuracy and reproducibility of a subpixel extended phase correlation method to determine micron level displacements in the heart. *Medical Engineering and Physics* 29 (2007) 154-162
- Lemmon CA, Chen CS, Romer LH. (2009) Cell traction forces direct fibronectin matrix assembly. Biophys J, 96(2):729-738.
- Lunsford Racing. (2010). Parts and Accessories. http://www.lunsfordracing.com/mm5/merchant.mvc.
- Norton, Louis A.; Andersen, Kim L.; Arenholt-Bindslev, Dorothe; Anderson, Lis, Melsen, Birte. (1995). A methodical approach of shape changes in human oral cells perturbed by a simulated orthodontic strain in vivo. *Archs oral Biology*, 40(9). doi:0003-9969/95.
- Rehfeldt F, Engler AJ, Eckhardt A, Ahmed F, Discher DE. (2007). Cell responses to the mechanochemical microenvironment--implications for regenerative medicine and drug delivery. Adv Drug Deliv Rev 59(13):1329-1339.
- Tanaka, Shiego M. (1999). A new mechanical stimulator for cultured bone cells using piezoelectric actuator. *Journal of Biomechanics* 32. 427-430.
- Torzilli, Peter A; Grigiene, Rita; Huang, Charles; Friedman, Steven M; Doty, Stephen B; Boskey, Adele L; Lust, George. (1996). Characterization of cartilage metabolic response to static and dynamic stress using a mechanical explant test system. *Journal of Biomechanics*. 30 (1). Doi: S0021-9290(96)00117-0.
- Triantafillopoulos, Ioannis K; Banes, Albert, MD; Bowman, Karl, PhD; Maloney, Melissa; Garrett, William E; Karas, Spero G. (2004). Nandrolone decanoate and load increase remodeling and strength in human supraspinatus bioartificial tendons. *The American Journal of Sports Medicine*, *32*(4). doi:10.1177/0363546503261700.
- Vandenburgh, Herman H. (1988). A computerized mechanical cell stimulator for tissue culture: effects on skeletal muscle organogenesis. *In Vitro Cellular & Developmental Biology* 24 (7). doi: 10.1007/BF02623597.
- Wang, D., Xie, Y., Yuan, B., Xu, J., Gong, P., Jiang, X. (2010). A stretching device for imaging real-time molecular dynamics of live cells adhering to elastic membranes on inverted microscopes during the entire process of the stretch. *Integrative Biology 5-6* doi: 10.1039/B920644B.

- Wang, H., Ip, W., Boissy, R., Grood, E.S. (1995). Cell orientation response to cyclically deformed substrates: experimental validation of a cell model. *Journal of Biomechanics* 28(12) doi:10.1016/0021-9290(95)00101-8.
- Wang, J., Li, B. (2010). Mechanics rules cell biology. *Sports Medicine, Arthroscopy, Rehabilitation, Therapy & Technology* 2(16) doi: 10.1186/1758-2555-2-16.
- Wang JH, Thampatty BP (2008). Mechanobiology of adult and stem cells. Int Rev Cell Mol Biol 271:301-346.).
- Wang J.H., Iosifidis, M.I., Fu, F.H. (2006). Biomechanical basis for tendinopathy. *Clinical Orthopaedics and Related Research* 433. doi: 10.1097/01.blo.0000195927.81845.46.
- Wang, J., Yang, G., Li, Z. (2005). Controlling cell responses to cyclic mechanical stretching. *Annals of Biomedical Engineering 33(3)*. doi: 10.1007/s10439-005-1736-8.
- Williams, J.L.; Chen, J.H.; Belloli, D.M. (1992). Strain fields on cell stressing devices employing clamped circular elastic diaphragms as substrates. *Journal of Biomechanical Engineering* 114(3) doi: 10.1115/1.2891398.
- WM Berg. (2010). MIN-E-PITCH 3D SERIES GEARS. http://www.wmberg.com/catalog/product.aspx.
- Yang GG, Crawford RC, Wang JHC(2004). Proliferation and collagen production of human patellar tendon fibroblasts in response to cyclic uniaxial stretching in serum-free conditions. Journal of Biomechanics 37(10):1543-1550.

# **6.1** – Appendix A – Pairwise Comparison Chart

This was filled out by Professor Billiar (the client) in order to determine the necessities of tasks to be completed for this design.

	Homeostatic	Biaxial	Inexpensive	Culture	Duration	Variable	Total	Importance
	Conditions	Stretch		Viewing	of	Duty		
				Area	Stretch	Cycles		
Homeostatic		0	1	1	0.5	0	2.5	2
Conditions								
Biaxial	1		1	1	1	1	4	1
Stretch								
Inexpensive	0	0		0.5	0	0	0.5	3
Culture	0	0	0.5		0	0	0.5	3
Viewing								
Area								
Duration of	0.5	0	1	1		0	2.5	2
Stretch								
Variable	1	0	1	1	0		4	1
Duty Cycles								

# 6.2 – Appendix B – Functions Means Tree


- 1. Strip Biaxial stretch 2-20% at different rates of loading/unloading
  - a. One motor with other side anchored (spreading anchor)
  - b. Pull at uniform strain rate
- 2. Equal Biaxial stretch To perform biaxial stretch
  - a. Two motors pulling at equal rates
  - b. One motor with linkage mechanism
  - c. One motor modified to pull in two opposing directions
- 3. Homogeneous strain area
  - a. Pull membrane in two directions
    - i. 4 motor system, all moving at the same rates
    - ii. 2 motor system in circuits
    - iii. 4 motor system with linkages
  - b. Possibly achievable with strip biaxial stretch
- 4. Control temperature and pH of the media to maintain homeostasis
  - a. Incubator
  - b. Cell humidifier
  - c. Hot air heater
  - d. Acrylic polymer cage
- 5. See visual cell response to stimuli
  - a. Microscope (inverted)
  - b. Photography
- 6. Variable Duty cycles
  - a. Stepper motor
  - b. Pneumatics
  - c. Hydraulics
- 7. Usability
  - a. Computer program to integrate all functions to perform when needed
  - b. Compatible with computer system

# 6.3 - Appendix C- Budget Breakdown


Date	Vendor	Desc.	Cost	Shipping	Total
1/19/2011	BBridge	Stretch Chamber	\$250.00	\$31.00	\$281.00
	McMaster-				
2/18/2011	Carr	Steel Rods & Sheets	\$64.21	\$9.75	\$73.96
2/18/2011	IGUS	IGUBAL Rod end	\$20.40	\$5.49	\$25.89
2/23/2011	IGUS	Rail & Float	\$57.71	\$5.49	\$63.20
3/21/2011	S&S Tech	Electric motor	\$129.99	\$10.95	\$140.94
4/11/2011	WM Berg	Pulleys & chains	\$65.76	\$50.00	\$115.76
		IGUBAL flange			
4/11/2011	IGUS	bearing	\$7.99	\$5.49	\$13.48
4/11/2011	Lyn-Tron	Standoffs (samples)	\$0.00	\$0.00	\$0.00
4/15/2011	Vangy Tool	Axis Arm	\$155.00	\$0.00	\$155.00
	Brierly				
4/15/2011	Lombard	M2 tap	\$14.00	\$0.00	\$14.00
Total					\$883.23
Budget					(\$624.00)

# 6.4 - Appendix D- Strain Parameter Examples

### **Examples of strain parameters**

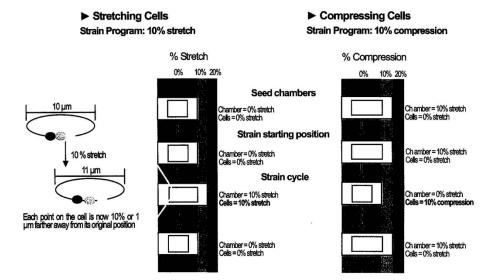

### ▶ 10% Stretch & 10% Compression of Cells Strain Program: 20% stretch or compression

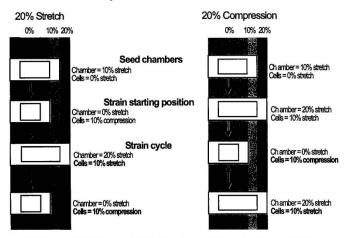

Use the Chamber Length Adjustment Knob to manually stretch the chamber. Use the knob to adjust for the seeding and start positions. Each 360° turn of the knob changes the total chamber length by 5%. or 1mm. ST-CH-04-XY chamber length is 20 mm.


V111405BC1 Rev103007

# 6.5 - Appendix E- General Use Manual

## 1. Strain Type Control

The arms in the X-axis & Y-axis have slots to control the type of strain, as the corner linear guide/crank shaft move from one angle to another the arm lengths should change as well. To move from equibiaxial strain to uniaxial strain as seen below four steps need to be taken:

- Loosening of the motor placement screw and switching from the shorter biaxial belt to the longer uniaxial belt.
  - The motor should be loosened and the belt should be removed first, however, the belt being put on should be placed and secured last.
- 2. Removal of crank shaft screw (4)
  - The crank shaft is also secured by four 4-40 screws, this must be done before the third step.
- 3. Loosening of X/Y-axis screws (8)
  - After loosening not removing the screws securing the x & y arms,
     slide and lock arms to desired position. These are M2x.4 screws and
     require a small flat screw driver.
- 4. Removal of linear guide screws (4)
  - Four M2x.4 screws secure the Igus linear guide in the top right corner
    of the device need to be removed to slide the guide to and from strain
    types.

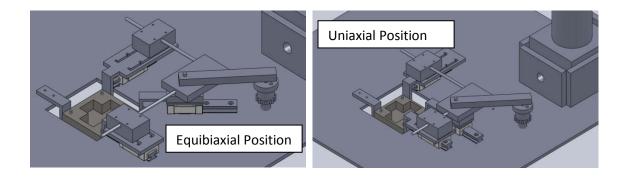

### 2. Strain Percent Control


Available strain percentages include 5, 10 and 15 percent strain. In the diagram below the arm is tagged by the red arrow, the housing connecting screw is indicated by the blue arrow and the strain percent positions are indicated by the black arrow (three 2-56 threaded holes), and the arm/crank shaft connecting point is indicated by the green arrow. Due to the closeness of the crank shaft holes the device will require separate arms for each strain percent. Changing strain percent is simple it requires two steps:

### 1. Removal of arm

• Two nuts and washers (green & blue positions) hold the arm down, they simply need to be removed to remove the arm.

### 2. Addition of new arm

• After removing the unwanted arm the new arm can be put in place (does not matter which way) then the washers and nuts should be securely re-tightened. Each arm should be labeled which strain percent it is for (5%,10%,and 15%).

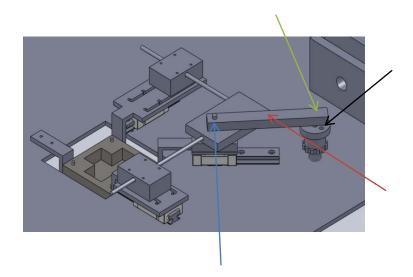

### 3. Motor Control

The motor being used puts out 8.4 Nm at 35rpm for torque which is more than enough to stretch the silicone wells. However, a 1:3 gear ratio is also being used, (this reduces the torque to 2.6Nm) which brings the top speed from 35 rpm to 114 rpm or 1.9 Hz. The motor is capable of turning clockwise or counterclockwise, and is also capable of variable speeds.

# 4. Running Device


When running the device to ensure wanted results the crank shaft must start at the zero strain point, this is to ensure the well is stretched and not compressed. Therefore before securing the belt in place the crank shaft should be already adjusted for whichever strain percentage is desired.

## 5. Strex Well Use

Adhesiveness and strength of chamber is not guaranteed for more than 3 uses. The chambers are durable enough to withstand approximately 900,000 stretches of 20%. Also the wells should not ever be stretched past 20%, because strain past 20% will significantly decrease the integrity of the silicone (B-Bridge International, Inc. 2009).

## 6. Mounting Strex Well

The Strex well has 4 pin holes that are 3mm in diameter, and these holes are used to mount the well to the stretch device. The holes are 30mm apart from their center, and 1/8" outer diameter standoffs will be used to secure the well. The 1/8" is slightly larger than 3mm, allowing the standoffs to provide a snug fit limiting vibration and motion in the Z-axis.

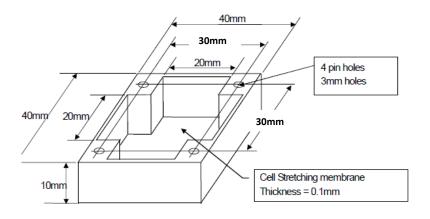

(B-Bridge International, Inc. 2009)

### 7. Autoclaving Strex Well

Sterilize chambers in an autoclave for 20 minutes at 121°C. The silicone chambers can withstand temperatures up to 180°C. Use of an autoclave is preferable. However, if an autoclave is not available, the chambers may be sterilized by

submerging them in 70% ethanol, rinsing with water, then drying in a sterile environment. Place the sterile silicone chambers in a Petri dish in preparation for coating (B-Bridge International, Inc. 2009).

- 8. Preparing Strex Well
- **Protein Coatings (Strex)** (B-Bridge International, Inc. 2009)

# **▶** Fibronectin Coating

Preparation of fibronectin solution:

- 1. Dilute human or bovine fibronectin to a final concentration of 50 to  $100 \,\mu g/ml$  in Phosphate Buffered Saline (PBS) Coating with fibronectin solution:
- Coating with Horonectin solution.
- 2. Pour 3-6 ml of the fibronectin solution into each strain chamber
- 3. Incubate at 37oC for more than 30 minutes.
- 4. Aspirate the fibronectin solution. If coating is successful, water will not be repelled after removing the fibronectin solution.
- 5. The liquid solution can be used to coat 3 or 4 chambers before discarding.

PBS (per liter):			
NaCl	8.00 g		
KCI	0.20 g		
Na₂HP₄ (anhyd.)	1.15 g		
KH <sub>2</sub> PO <sub>4</sub> (anhyd.)	0.20 g		
Note: Dulbecco's PBS in powder form for tissue culture applications is commercially available.			

(B-Bridge International, Inc. 2009)

## **▶** Gelatin Coating

Preparation of gelatin solution:

- 1. Add gelatin powder to PBS at a concentration of 2%
- 2. Autoclave the mixture to dissolve and sterilize

Coating with gelatin solution:

- 3. Pour 3-6 ml of the gelatin solution into each strain chamber
- 4. Incubate at 37°C for more than 30 minutes.
- 5. Aspirate the gelatin solution. If coating is successful, water will not be repelled after removing the gelatin solution.
- 6. The liquid solution can be used to coat 3 or 4 chambers before discarding.

## ► Collagen Coating (Cellmatrix 1-C, P, Type 3 or 4)

Preparation of collagen solution:

- 1. Combine 1 part collagen to 10 parts HCl, pH 3, in a sterile tube Coating with collagen solution:
- 2. Coat chamber with a thin layer
- 3. Aspirate excess
- 4. Dry in biological safety cabinet at 25°C or below. The chamber can be stored at the same temperature.
- 5. Wash the chamber twice with culture medium. If coating is successful, water will not be repelled.
  - **Cell Culture** (B-Bridge International, Inc. 2009)

Seed cells at the appropriate concentration in the freshly coated chamber. Do not over expose the cells to dissociation enzymes. Cells should be treated in the same manner (type and concentration of enzyme, temperature, and time for digestion) for all experiments.

Cells should not be cultured at a high cell density in the chambers. For example, epithelial cells often form a cell-sheet and the cell-cell adhesion seems to be stronger than a cell-surface adhesion. When this happens cells may detach from the chamber. Additionally, cultures that are grown over a week in the chambers may detach.

After overnight incubation, inspect cells with the microscope to ensure that they have adhered to the chamber.

### 9. Strex Well Issues

The PDMS (silicone) chambers are unfortunately very hydrophobic with two methyl-base on the surface. Cells adhere to surfaces coated with fibronectin or collagen through integrins. Compared to adhesion on plastic or glass wells Strex wells are slightly


different because they are not charged. If the cells on the Strex well are not attaching or are detaching during stretch, a higher concentration of either fibronectin or silicone may be necessary. These coatings of course must be applied before culturing the cells. (B-Bridge International, Inc. 2009)


### 10. Additional Information


- If there are wrinkles or bubbles on the bottom surface of the strain chamber when seeding cells this is because of the wells thin structure. The well is designed not to have any heterogeneous regions, though it is difficult to achieve. B-Bridge recommends using a small volume of ethanol in a Petri dish and lightly placing the chamber in the culture dish starting at one edge and moving toward the opposite edge of the chamber to remove air bubbles between the dish and chamber. The ethanol obviously must be evaporated before spreading a cell suspension on the chamber. (B-Bridge International, Inc. 2009)
- Obtaining protein or mRNA samples from the cells:
- 1. Proteins for Western blotting: Wash the cells once with PBS. Add SDS-PAGE sample loading dye directly into the chamber, and collect the cell extract by using a cell scraper (B-Bridge International, Inc. 2009).
- 2. Proteins for Immunoprecipitation: Wash the cells once with PBS. Add cell extract buffer directly into the chamber, and collect the cell extract by using a cell scraper (B-Bridge International, Inc. 2009).


3. RNA: Wash the cells once with PBS (for RNA preparation). Add RNA extraction buffer directly into the chamber, and collect the cell extract by using a cell scraper (B-Bridge International, Inc. 2009).


### 6.6 - Appendix F- Design Drawings

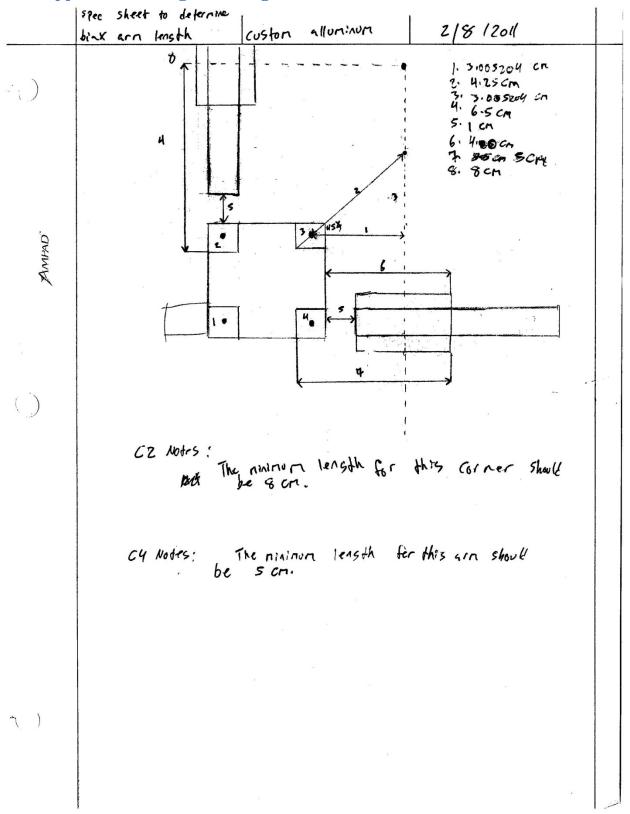

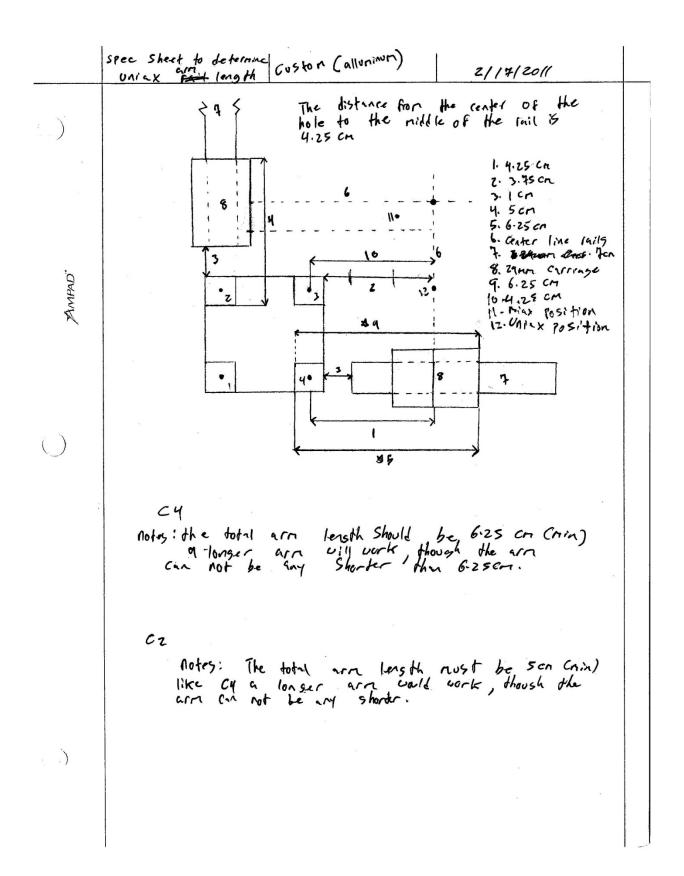

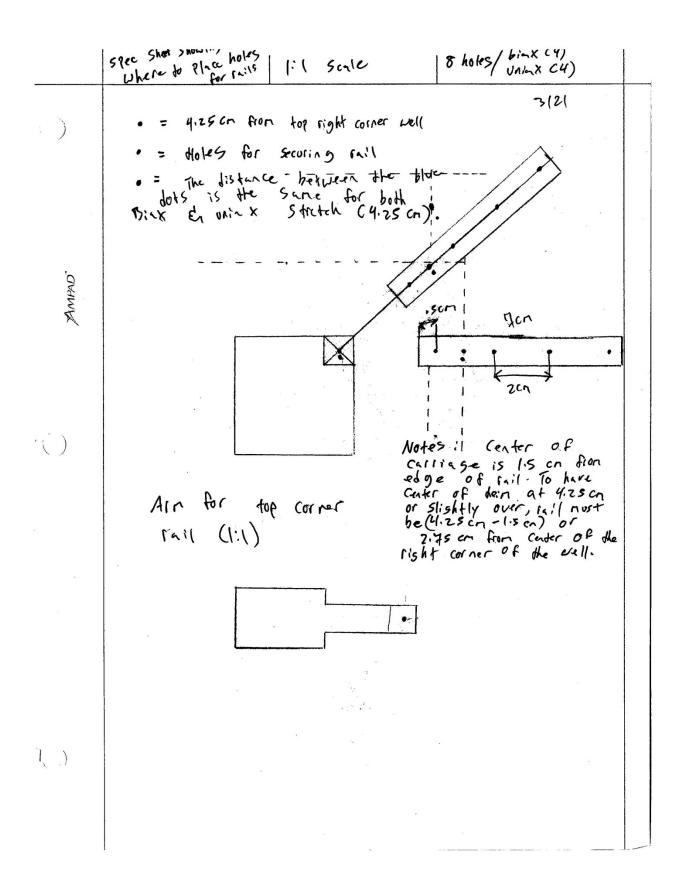


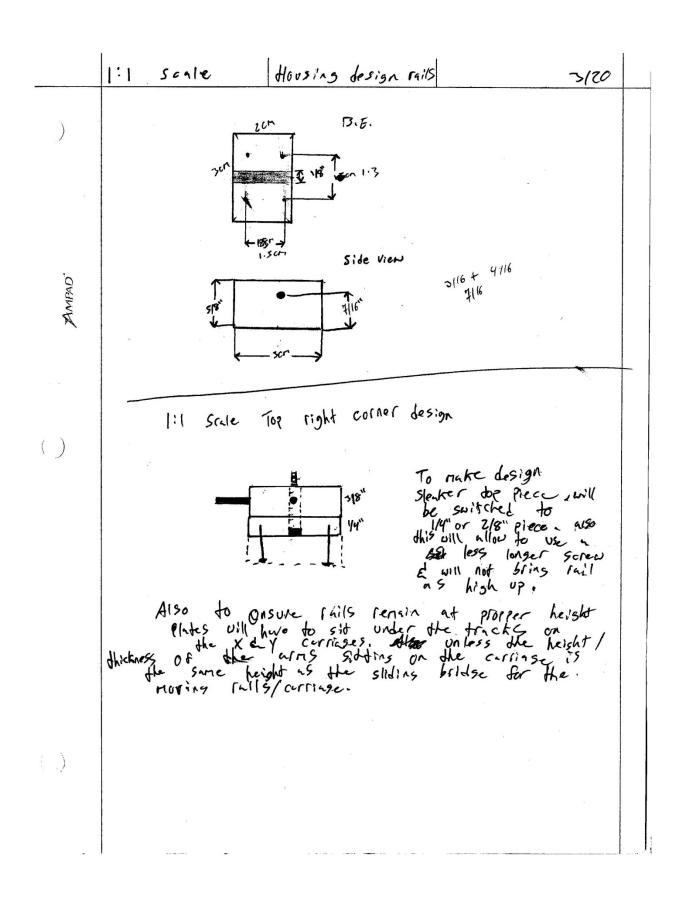

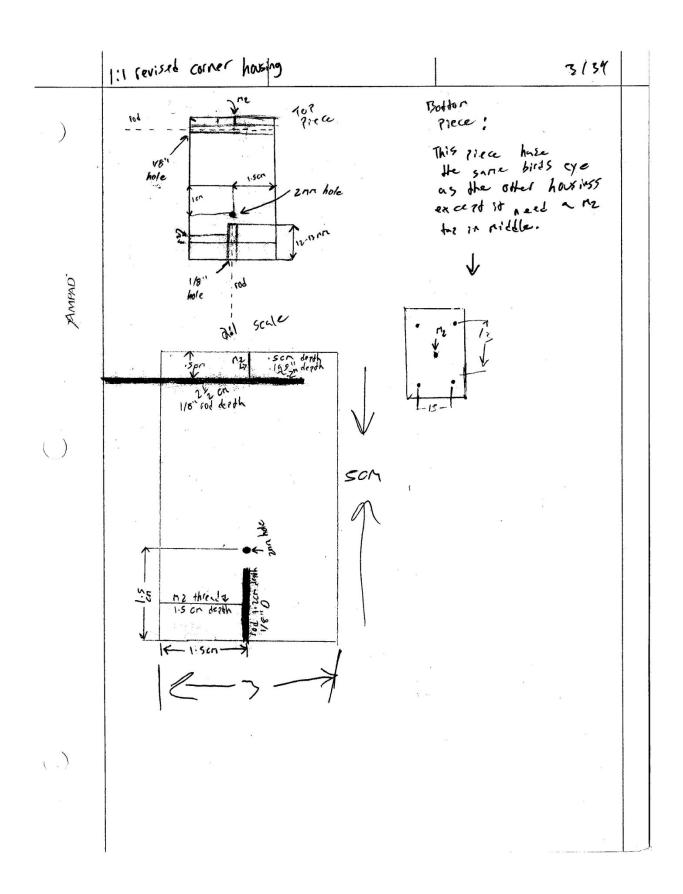


AMPAD

(, )

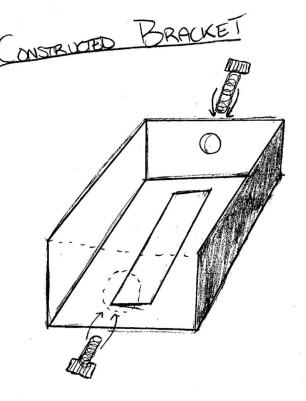



# LEG SUPPORTS

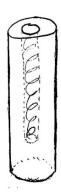


- Icm Dian
- 30mm height

  - height (estraints)
     should leave ample room for bett
- 6/32 screw
- Not located in each corner
  - · leave space for best position in equi/uni stretch

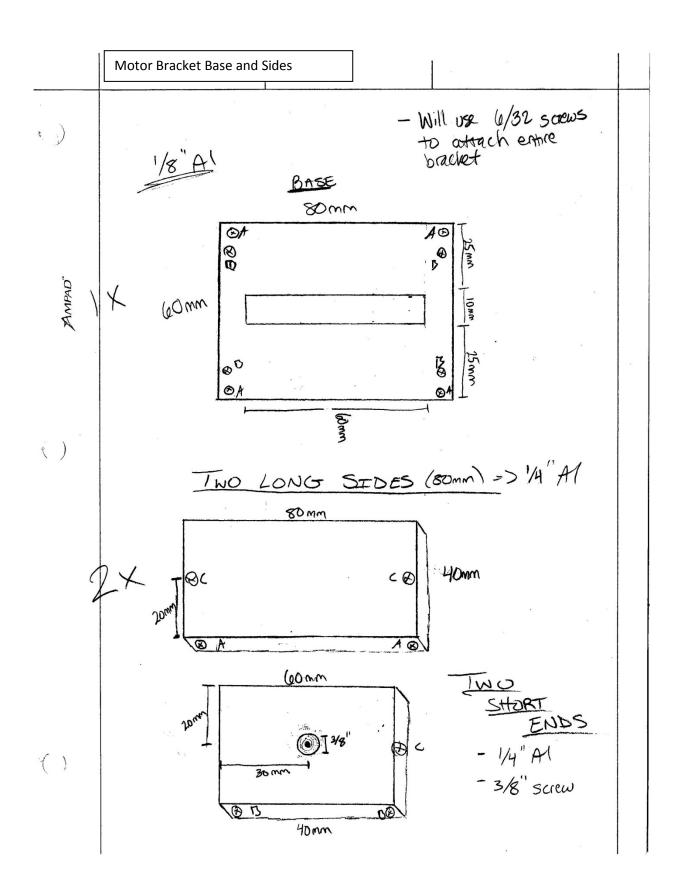


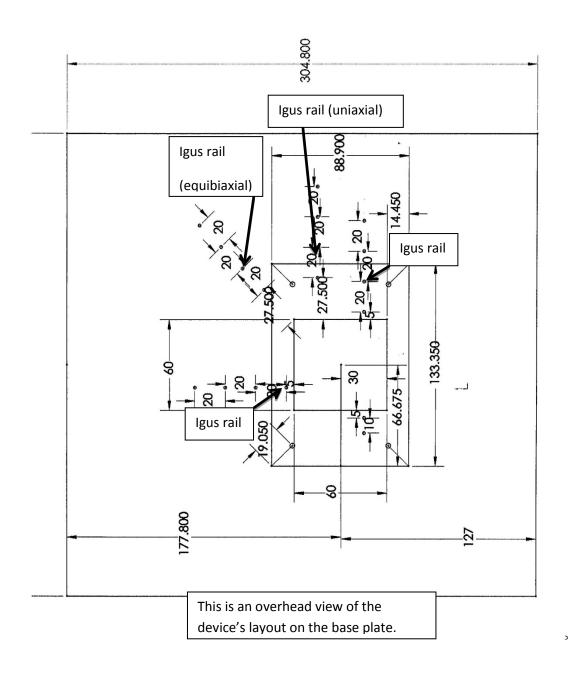

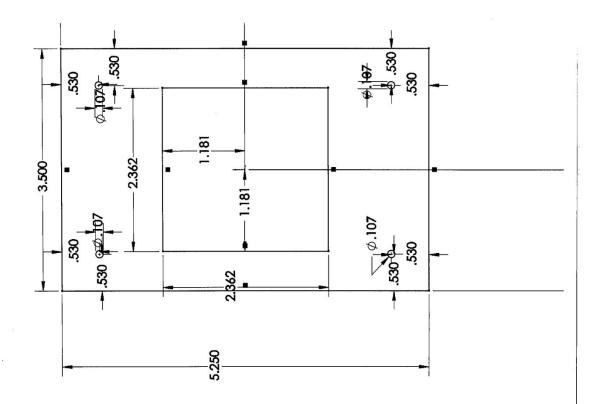

This is the X-Y stage adaptor that attaches below the base in order to maintain a level device while on the microscope stage.




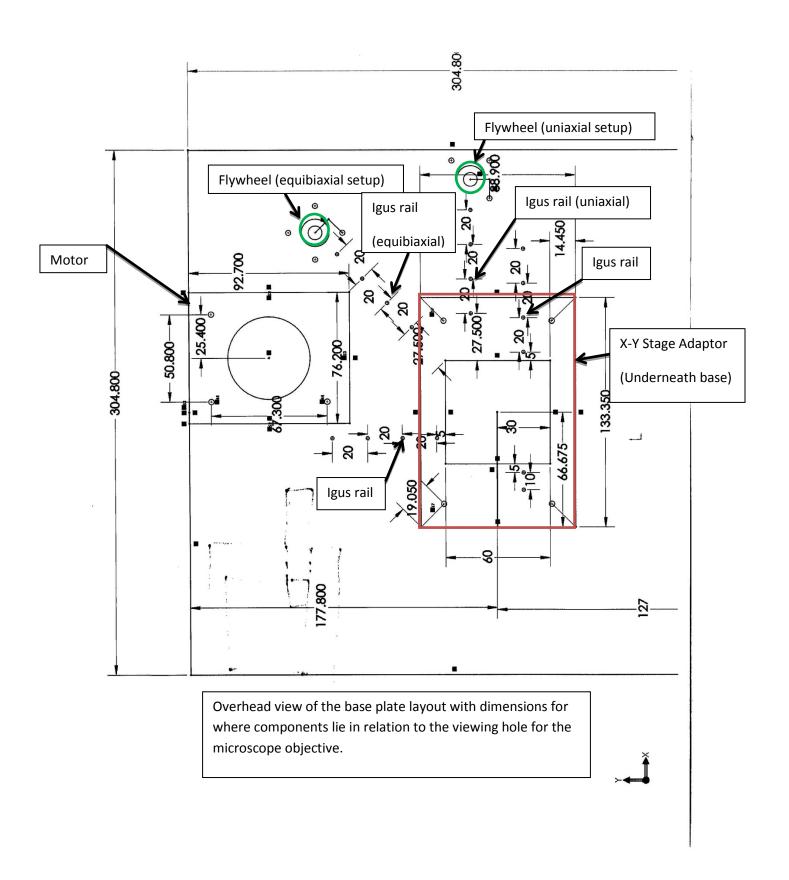

## 6.7 - Appendix G- Ordered Parts

## 3-D PULLEYS => 3 day lead time

CIR. PITCH	BORE	STYLE	MATERIALS	BERG'S ® NAME	BELT
4mm	Ø8	PIN HUB	Aluminum per DIN 3.1355 Anodized	Min-E-Pitch® Three D Drive	Operates with 3TDF Series

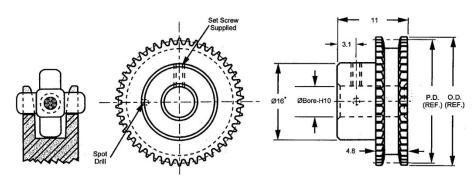

STOCK	NO. OF	PITCH	OUTSIDE
NO.	TEETH	DIA.	DIA.
3TDP263A-12	12•*	15.22	16.80
3TDP263A-13	13•*	16.49	18.07
3TDP263A-14	14.*	17.76	19.34
3TDP263A-15	15•*	19.03	20.61
3TDP263A-16	16•	20.29	21.87
3TDP263A-17	17	21.56	23.14
3TDP263A-18	18	22.83	24.41
3TDP263A-19	19	24.10	25.68
3TDP263A-20	20	25.37	26.95
3TDP263A-22	22	27.90	29.48
3TDP263A-24	24	30.44	32.02
3TDP263A-25	25	31.71	33.26
3TDP263A-26	26	32.98	34.56
3TDP263A-28	28	35.51	37.09
3TDP263A-30	30	38.05	39.63
3TDP263A-32	32	40.59	42.17
3TDP263A-35	35	44.39	45.97
3TDP263A-36	36	45.66	47.24
3TDP263A-40	40	50.47	52.32
3TDP263A-45	45	57.10	58.66
3TDP263A-50	50	63.42	65.00
3TDP263A-55	55	69.76	71.34
3TDP263A-60	60	76.10	77.68
3TDP263A-65	65	82.45	84.02
3TDP263A-70	70	88.79	90.37
3TDP263A-75	75	95.13	96.71
3TDP263A-80	80	101.47	103.05




<sup>\*</sup> Sprockets with Ø19.0mm P.D. and smaller are recommended for idler use only.
For 12-16 teeth, hub diameter equals 10.0.
Other numbers of teeth are available on request.
Stainless Steel equivalent available.
Teeth could be anodized.

## 3-D BELT =) I day lead time



4mm	Polyurethane (Green) 8mm Dia. Stainless Steel Cable*	Min-E-Pitch®	Operates with 3TDP, 3TF, and 3MTB series.
GIRCULAR PITCH	MATERIALS	BERG'S ® NAME	PULLEY

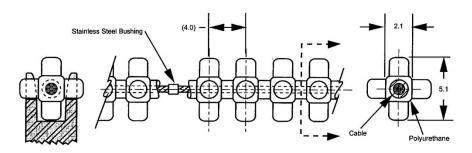

STOCK NO.	NO. OF PITCHES	LENGTH (Ref.)
3TDF-30-E 3TDF-35-E	30 36	120.0 144.0
3TDF-40-E 3TDF-45-E	40 46	160.0 184.0
3TDF-50-E	50	200.0
3TDF-55-E 3TDF-60-E	56 60	224.0 240.0
3TDF-70-E	70	280.0
3TDF-80-E	80	320.6
3TDF-90-E	90	360.0
3TDF-100-E	100	400.0
3TDF-110-E	110	440.0
3TDF-120-E	120	480.0
3TDF-130-E	130	520.0
3TDF-140-E	140	560.0
3TDF-150-E	150	600.0
3TDF-160-E	160	640.0 680.0
3TDF-170-E	170	
3TDF-180-E 3TDF-190-E	180 190	720.0 760.0
3TDF-190-E	200	800.0
3TDF-200-E	210	840.0
3TDF-210-E	220	880.0
3TDF-230-E	230	920.0
3TDF-240-E	240	960.0
3TDF-250-E	250	1000.0
3TDF-260-E	260	1040.0
3TDF-270-E	270	1080.0
3TDF-280-E	280	1120.0
3TDF-290-E	290	1160.0
3TDF-300-E	300	1200.0
3TDF-310-E	310	1240.0
3TDF-320-E	320	1280.0
3TDF-330-E	330	1320.0
3TDF-340-E	340	1360.0
3TDF-350-E	350	1400.0
3TDF-370-E	360	1440.0
3TDF-400-E	400	1600.0
3TDF-440-E	440	1760.0

- For Field Splice Kit order 3CCR-7.
- Available with Aramid Core. See page A-2 for alternate belt construction and reverse bending applications.
   Reduce Operating Load Rating by 50% for field splices.

- Three Dimensional Drive
  Zero Backlash
  Silent Drive
  Infinite Lengths
  90A Durometer
  No Lubrication
  Mass11 Grams/Meter
  Positive Drive
  Ultimate Tensile Strength 222N 40N/PIN
  When belt is used in open loop configuration
  (without splice, no crimp bushing).
  Temperature Range + 82°C to -26°C
  Recommended Operating Load 53N
  Recommended Max. Operating Speed 1.91 M/S




BULK FOOTAGE - NOT SPLICED							
STOCK							
NO.	LENGTH						
3TDF-1.5M	1.5M						
3TDF-3.0M	3.0M						
3TDF-7.5M	7.5M						
3TDF-15.0M	15.0M						
STDE SO OM	30.0M						

WWW.WMBERG.COM

**MA 35** 


1-800-232-BERG (U.S.A ONLY)

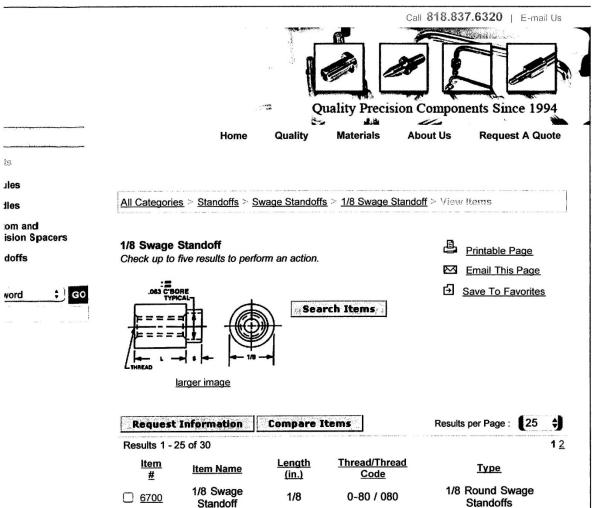

	. 100 1 10 10 10 10 10 10 10 10 10 10 10		- Alexandria de la compansión de la comp	
Results 1 - 2	25 of 30	CONTRACTOR OF STREET	THE RESIDENCE OF THE PARTY OF T	12
ltem #	Item Name	Length (in.)	Thread/Thread Code	<u>Type</u>
<u>6700</u>	1/8 Swage Standoff	1/8	0-80 / 080	1/8 Round Swage Standoffs
□ <u>6701</u>	1/8 Swage Standoff	3/16	0-80 / 080	1/8 Round Swage Standoffs
☐ <u>6702</u>	1/8 Swage Standoff	1/4	0-80 / 080	1/8 Round Swage Standoffs
□ <u>6703</u>	1/8 Swage Standoff	5/16	0-80 / 080	1/8 Round Swage Standoffs
□ <u>6704</u>	1/8 Swage Standoff	3/8	0-80 / 080	1/8 Round Swage Standoffs
☐ <u>6705</u>	1/8 Swage Standoff	7/16	0-80 / 080	1/8 Round Swage Standoffs
<u>6706</u>	1/8 Swage Standoff	1/2	0-80 / 080	1/8 Round Swage Standoffs
	Otaliaon			


 $http://vmgengin\,eering. tho mass net. com/vlew items/standoffs-swage-standoffs/1-8-s$ 

Page 1 of 2

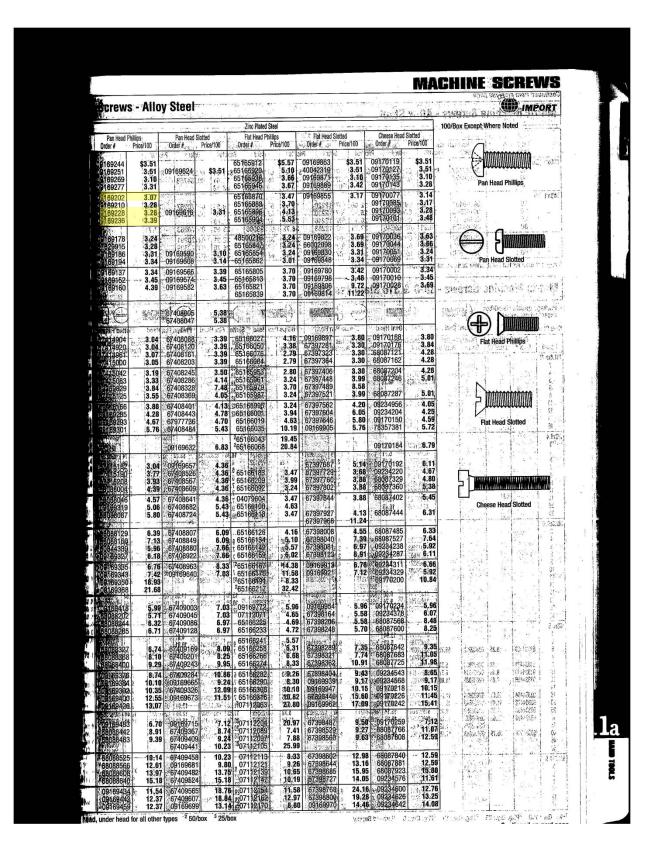


	ic Ma	inine 	terna kin	loo:	athors is not seen as	stíú Í	1.45	ated Steel	got2 ho x 10 or 17	11 1	Choose Bood SI	otted	SARCO SARC
V16x1 (C	Length (mm), ont'd)	llyddi.	Pan Her Order # 0916946	nd Phillips Price/100	Order#	ead Slotted Price/100	Order#	Price/100	Order #	Slotted Price/100	09234659	rice/100 \$14.97	8
M8x1.25	60 80	ecol e	0916947	17.63	6740964 0916970	3.17 99.70	30711219 30711221	816	67398842 09169988 09169996	33.18	09234758	19.63	37.00 7.00 20.00 80.00
 (1)	16 20 25 30		0916949 0916950 0916951 0916952	9 23.00 7 24.74	09 697 674096 091697 091697	26.14	20711222 20711223 20711224 30711225	8 19.45 6 23.16	6739888 67398925 67398966 67399000	24.83	09170267 09170275	22.24 24.74 27.71	97.26
(fi)	35.4		0916953 0916954 0916955	26.56 1 29.28	091697 674096 091697	29:73 2 31:27	20711226 20711227	28.59 30.10	67399046 6739908 267399121		09170291 09170309 09170317	29.93 31.59	6.7 <b>6</b> a10 <b>6</b> a-1 <b>6</b>
1.5	verall for fi	-	under head	for all other	types <sup>2</sup> ,50	0/box <sup>3</sup> 25/b	ox - 04.	3 721 03 M 3 721 03 157 8	-2042 -2042 -2149	26 G	9565 20109574 20109574 20109572	3.9.6 3.9.6 4.39	
Packag	e quantity	is 50/bo	Margaret 1	ws - 18°	8 & 31	6 Stain	ess Ste	<b>(6)</b> 0   37.8	1585 1589	Sporting	<b>(P)</b>		
except	where not less Steel Slotted O		China		hillips Oval	at an har	otted Pan Hea	id Phillips	Pan Head	Slotted Flat He	ead Phillips F	Pozi Flat	Slotted Fill
L (mm) 1.6x0.35	Order #	Price/100		Price/100 0r	ier# Price	Oluei #	1.387	rder# Price/10 EETal ers EETal	Order#C	Price/100 COrd	er#   Price/100 160   Price/3	3.05 1 3.19	Price/100
5 6 8	andani	e i i i Nonadi	880050 (58021	1801 1801 5.01	997629	5.92 8.20	3 <b>19</b> 3 <b>19</b> 3 <b>19</b>	6739 6739 24 6739	86725058 86725056 86725074 86725082 86725090	69.86 7 E9.86 4 E9.86	233, 06530a 05630a 93060ay	3.33 3.84 3.55	6.0
10 12 14	Detro		τ 330. 8	4.05 7.25 4.59 5.72	204 40t	.20 08224 .05 06224 .80 08170 .76 78357	645	6739 8.94 6739 4.63 6739 8.19 (0916)	86725108 86725116 1910a	42.69 4 642.69 4	150 .0-8:11 150 E. 151 E. 213 .55:14	3.85 6 4.28 6 4.67 6 5.76 6	000 m
16 M2x0.4 '3 4	73291 4	sales Tarana	5891271 089166	\$879 IR	997637	5 95 N		29E 25.E	8043 8008 7 86725124	100 96 100 83 <b>8</b> 1004 853	999. 829. 52098.0	02627016	\$4.74
5 6 8	reletie A	en 1999	ieżd, dedi	F F 88 7887 1880	997652 997660 997678	5.51   185601102 5.15   85601100 6.85   185601128 6.89   185601128	2.81 0 2.86 0 2.95 0 3.34 0	633709 7.3 633717 4. 633725 6. 633733 5.	8 186725132 1 186725140 5 186725157	4,30 853 6,15 853 6,86 853 7,06 853	10761 2.63 10779 2.42 10787 3.20 10795 9.37 10803 9.80	02637024 02627032 02627040 02627065 02627065	3.86 5.07 7.18 5.85
10 12 12 16 20 20	Slotted	hasH es	Che	Br 2. Kd	997694 997702 SO4	9.48   85601451 0.90   3   8   8   8   8   8   8   1   1   1   1	3.67	833741 X	186725173 186725181 186729193	7.06 858 7.85 9.72 6 10.45 a	10803 9.80 Sensi	02627065 1 G. A 0 00.6 0 08.7	6.29
25 M2.5x0.45 3 4		1	1/4	SE.ii	\$5.7 857700 527 857700	780304 EE.	300	4.15 6730	6126	13 200	1.029047 1.029047	0   es.a 002627073 002627687	5. <b>9</b> 0
5 6 8	a s: VCs	al ar	e de Refer Arker des	\$968 £106	012368 316657 995221	8.21 9560 1777 3.52 98560 1786 3.84 48560 2193	4.46.130	2633766 47 2633774 5: 2638762 2638762	19 186725223 19 18672523 14 12/86725249	7.85 (2053	10829 4.87 10837 3.34	02627686 02627699 02627107 02627107 02627115	6.00 6.71 6.71
10 12 14 16	() <u>4</u> (())		say Fastenar	1854 18 ±	997736	5.40   85601201 8.70   85601219   85601227   9.86   85601235	5.47 0 5.42 5.91 0	2633808 5. 2633816	186725266 5-1-5-3 18 - 186725272	10.20 853	10852 8.25 10860 11.30	02627028 86.1 02627131	9.54 9.54
20 25 30		3 e 1	log tystener Any rasienar i Strengto Ti unny Liucis	70.	997769	9.74 185601243 12.72 10 86 18 86	310	2633824 8. 2633832 17.	- 43000	7.100.51	10878 500-7016 21-1601 21-1601 21-1601	02627149 002627158 002627158 0 026 A	9.54 9.74
M3x0.5 4 5	93334423	\$8.25	en Under W en Under W 102632792	\$5.73 3.66	0077777	13.94 5.63   68019165 9.74   68019207 6.53   68019249	a come or bear his	2633840 4. 2633840 83.5	22 8672530 14 8672531	9.23 680 7.27 680	10886 4.01 18001 10.01 18048 8.83	0263) 663	- GÉ
7 8 10 12	93334431 93334449 93334456	8.13 9.39 8.83	02632818	8.77 COS	997827 893	: U AX EXIGNITIVE AND	N 5.95-1740	2633865 E 10 2638873 O 2 69 2638881 O 8 8	14 10.8672532 53 10.8672533	4.93 1680 1 3 6.93 1680	18126 -8: 11.29 18167 -3: 11.27	02631679 002631687 002631696	4.35 4.35 9.74
14 16 20 25	93334464 93334472 93334480	10.67 11.00 11.58	02632834 02632842	815	997835 997843 997850	11.19   68019322 11.30   0   6 f. 11.58   68019863 12.22   68019405 15.14   68019421		2688899   0.110 2688997   9.8 7 2688915   0.8 4	18 86,25355 55 86,2586	7.82 1680	10894 ) 6.23 18241 12.99 18282 13.91 10902 12.68	002631708 002631715 02631029 02631737	7,67 9,74 9,96 10,44
30 35 40	93334498 100000 293334506	22.60 55.79	02632867 02632875	12.29 548 19.00 148 9.39 8 11.22 648 11.12 708 11.48 628	997876 997884 997892	12.22 1880 9401 15.14 1680 9421 16.35 8560 256 21.82 8560 268 25.24 8560 263	11.30 10 7.07 8.10 ac	3633931 - 11	58 8672539 8672540	20.91 853 (20.91 853 (16.06 853	10910 15.01 10928 18.24 10936 18.59	002631746 6 re.g	15.49
45 50 M3.5x0.6	Squant issaur Selecti	100	Prije Sealani Prije Sealani Prije Sealani Nije and Hiyo	04.		19.04 10 20 20.04 10 20 20.04 10 20		THE PARTY OF THE P	32 8870F49	9 0.58 07 0 10 8 07 0	7:09441 163681	0 15.0 0.263 U.55 0.263 U.55 0.263 1760 0.263 1778 0.263 1778	- 4
6		120		188 188	997918 1997926	18.01 0 20 38.21 0 20 40.28 8560128 35.48 8560129	i i	2633949   10 2633956   50 2633964   49 2633972   6 g	32 8672543 86 8672544 82 8672545 39 8672646 28 8672847	27.99 27.99 27.99 27.99 853 29.14	400462 44.6 440018	02631760 -02631778	35.46 4.56 8.38 11.48 9.52

		WS -	18-8 &	k 316 Pan Head	Tree	less &												
		Price/100		Price/100		Price/100	Pozi Po Order#	(Price/100		Flat Head Price/100		Flat Head . Price/100	Pozi F Order #	lat Head Price/100		llister Head Price/190	Slotted Order#	Truss Hea Price/
			81997967 81997975 81997983 81997991 82432519	\$18.47 13.87 42.45 19.84 22.24	4	Litus Lit. 15 Lit. 15 Lite! 16	1,82816 1,0 <b>2634012</b> 1,82818 1,81819 1,81818 1,81818 1,81818	\$39.46	86725496 86725504 86725514 86725520 86806049	33.75 13.87 15.83			02631810 02631828 02631836 02631836	12.16		\$16.64	4/2 10° 5-36/60° 5-36/60°	1
		\$14.77, 5.58 5.33, 15.43	82432527 82432535 82432543 82432550 68013002 68013044 82432568 68013086	26.26 12.62 10.75 11.65 13.18 18.96 19.95 18.67	68019488 680195201 680195611	3 cr14.09 1.89:10.46 1.86:12.36	02634046 02634053 02634061	14.22 8.02 8.70 8.13 9.96 5.84 11.12	85806056 86806064 86806072 86806098 86806098	12.53 12.53 11.28 11.85 14.22 13.89	85310969 168018324 168018365 168018407 168018449	13.22 10.56 11.03 10.60	102631851 002631851 002631869 02631885 02631893 02631991 02631919	8.25 5.84 9.52 7.07 13.33 10.32	02623791	10.35	02626083 02626091 02626109	19 19
	7. 10. 20.	20.89 25.56 14.13 18.36 25.36	68013127 68013168 82432576 82432584 82432592 82432600 82432618 82432626	20.06 20.63 30.64 23.28 35.09 32.93 38.33 65.53	68019645 68019660 680196863 85601326 680197021	20.63 18.19 26.36 21.11 21.28.88 20.31	):92634103 (02634111 (02634129 28387827 (02634145	22.48 22.48 26.16 31.34	86806114 86806122 86806130 86806148 86806155 86806163 86806189 86806189	13.94 15.87 15.58 15.58	68018522 68018563 68018605 85310977 85310985 85310993	13.01 18.80 19.77 22.08 23.89	001 689 02631935 02631943 02631943 02631968 02631968 02631976 02631984	8.83 17.79 22.72 27.32 30.88 33.26 40.97	02623858 02623866 02623874 02623882 02623890 02623908 02623916 02623924	19.72 22.08 22.31 23.89 24.61 27.45 34.66	02626125 02626133 02626141	
	0006 0014 8022 017 8030	21.45 11.61 21.16 22.92	82432634 68013366 82432642 68013440 82432659 82432667	29.14 17.35 25.04 22.75 28.68	168019744£ 168019769 168019785£ 168019801£	24.66 21.45 21.06 27.93 21.53	035878* 002634152 002634657 002634665 002634673 002634681 002634699	29.03 22.60 17.64 15.95 29.94 24.91	86806205 86806213 86806221 86806239 86806247 86806254	23.51 17.38 15.20 18.89 26.36 16.06	85311009 168018647 168018688 168018720	26.16 14.27 14.27 13.78 15.74	02631992 .02632008 02632016 02632024 02632032	13.64 17.79 9.39 25.92 13.33	02623932 02623940 02623957 02623965 02623973 02623981 02623999	16.60 18.59 17.02 17.38 27.06 23.89	02626679 02626687 02626695	25. 24. 26.
	3048 8055 263063 1633071 663089 4417 683097	21.53 40.42 22.72 33.41 38.44 38.56	82432675 82432683 82432691 82432709 82432717 82432725 82432733	35.91 36.49 62.33 63.36 56.83		31.35 32.74 39.64 29.88 38.21	02634707 02634723 02634731 02634749 02634749	22.37 39.46 41.20 38.91 49.56 52.08	86806262 86806288 86806288 86806296 86806304 86806312 86806320 86806338	23.05 16.27 28.56 35.91 45.00 45.00 45.00	68018803 68018845 68018886 85311017 85311025 85311033	15.98 26.67 33.16 23.89 31.82	02632040 02632057 02632065 02632073 02632081	10.20 23.97 31.10 45.20 43.27	02624005 02624013 02624021 02624039 02624047 02624054 02624062	23.89 31.05 32.11 36.23 43.72 52.44 70.11	02626711 02626729 02626737 02626745 02626752 02626760 02626778	38. 41. 42. 31. 37. 42. 79.
10. 10.	0.683105 Cao 0.167.19 0.167.19	39.72	82432741 82432758 282432766 282432774-282432782 82432782 82432782 82432790 682432806 682432816	123.51 69.20 78.74 81.14 108.47 120.30 136.32	10810968 101 101 1041700 1041700	(16)0 long	02634772   90   8   6   6   7	100	86806346 86806353 811 191211117 86806361	69.47	18 <b>53 1047</b> 1 Head 1 Head 1 Prige/100	51.88 s	<b>02632907</b>	45.91	02624070 002624088 02624096 02624104 02624112	70.11 70.11	1 02626786 19312 204 19312 204 19312 4 19312 19312	(0) 14(1) 100.43 E (0) 1
	0283313 02833126 02833139	21:26 S 22:76 T	6. F f 82432824 £ 82432832 d68013846 d 82432840 b 82432857	25.01 25.01 26.24.20 23.61.0 33.49 36.27.45	08.8 85601409 68019967 68020007 85601417 68020049:	24.32 38.06 29.93 31.66 29.84	02634780: 02634798 02634806 02634814 02634822 8.50	35.78 1 32.48 9 18.36 0 43.37 S 26.07 9	\$ 7.80 2868062870 1868062955 \$86806403 586806413 \$86806429 \$ 85	29.59 20.89 25.45 30.53 23.31	68018928 68018969 E.DE	23.76 20.71 21.76	02632115 02632123 02632£31	22.60 19.84 26.07 20.09	02624120 02624138 02624146 02624153 02624658 02624666	33.04	02626794 02626802 02626810 02626828	41.6 40.4 40.0 43.6
17. 1970 -21. 23.	02633659 02633659 02633659 0263367 0.) 10065 02633675	49.46 S	282432931	26.26 60.14 53.24 62.33 63.22 90.90	19.43	40.70 42.76 48.70 53.32 01.01 S 54.74 S	02634830   02634848   02634855   02634863   02634874   02634889   02634897   8 SO I	43.37 51.54 70.11 66.90 V 91.35 8 71.16 0	86806445 86806452 86806452 86806460 266806478 86806478 86806494 1 6 6	25.36 37.32 53.49 53.49 63.49 7	86311082	24.70 32.17 41.03 33.62 39.19 51.56	02632156 02632651 02632669 02632677 02632685 0 50 1	16.64 25.72 41.20 50.27 50.97 6 103.6	02624674 02624682 02624690 02624708 02624716 02624724 02624732	31.43 35.24 36.23 42.39 72.32 72.45	02626836 02626844 02626851 02626869 02626877 02626885	52,7 055.4 52.3 53.7 48.2
* 8.0	938 2 944 <b>4</b> 954 <b>2</b>		282432949 282432956 282432964 282432972 282432980 282432988 282433004	97.55 146.14	888.6088 9.08 17.58 90.78 20.70 77.28	64.29 S 180 S 280 S 20 S 20 S	02634905 6250	90.69 8 90 7 64 9 9 1 2 32	868065194 2868065194 2868065284 287828612 87828620 87828638 87828638	7 81.819 2 ep.011 29.48 7 80.79 4 19.811	8801086 7.0.3	65.81	1.2.7	92	02624740 02624757 02624757 10600 02624765 02624773 02624781	84.01 84.01 84.01 106.84	02626898 	57.10 10 10 10 10 10 10 10 10 10 10 10 10 1
13.28 13.88 12.56 21.46 25.71 43.27 10.50 43.27 14.33	2 585	(F.SC)   M (G.SC)   M	285180003 285130011 285130029 285130037 285130045 285130052	80.58 F.8	86.85 88.69 5601466 5601474	74052 \$20654 \$3,000 \$0,000 \$0,000 \$0,000 \$0,000 \$0,000 \$0,	2634913 2634913 2634947 2634947 2634954	87.47 8 87.47 8 154.85 157.93 01 162.99 21	87828653 87828661 87828679 87828687 87828687 87828695 87828703	98.83 C 70.27 Of 41.43 43.61 F 62.00 C 80.80 80		10 10 10 10 10 10 10 10 10 10 10 10 10 1	2632749 2632727 2632727 2632736 2632743 2632760	59.20 59.33 89.63 111.24 a	02624799 02624807 02624807 02624815 02624823 02624823	54.06 50.93 66.57 65.81 68.70	02626901 02626919 02626927 02626935,	92.96 112.84 122.36 126.26
15.88 19.45 18.45 46.13 48.60	200 2 017 <b>3</b> 125 <b>3</b>	200 <b>3</b> 200 <b>3</b> 200 <b>3</b>	85130060 85130078 85130086 85130094 85130102 85130110	149.36 128 190.10	36.52 284888 31.36 34.74	214.42 St	2634962 2634970 2634988 2634996	169.20 221.21 @ 114.91 1 284.00	87828711 87828729 878287378 87828745	123.39 106.46 81 106.40 106.40	510 810	) 1833 - 183 193	2632768 26327764 41.61 2632784	146.48 01 102.97 3 150.61 3	02624856	87.68 118.22 132.91 146.36	02626943 02626950 02626968 02626976	115.14 151.29 164.72

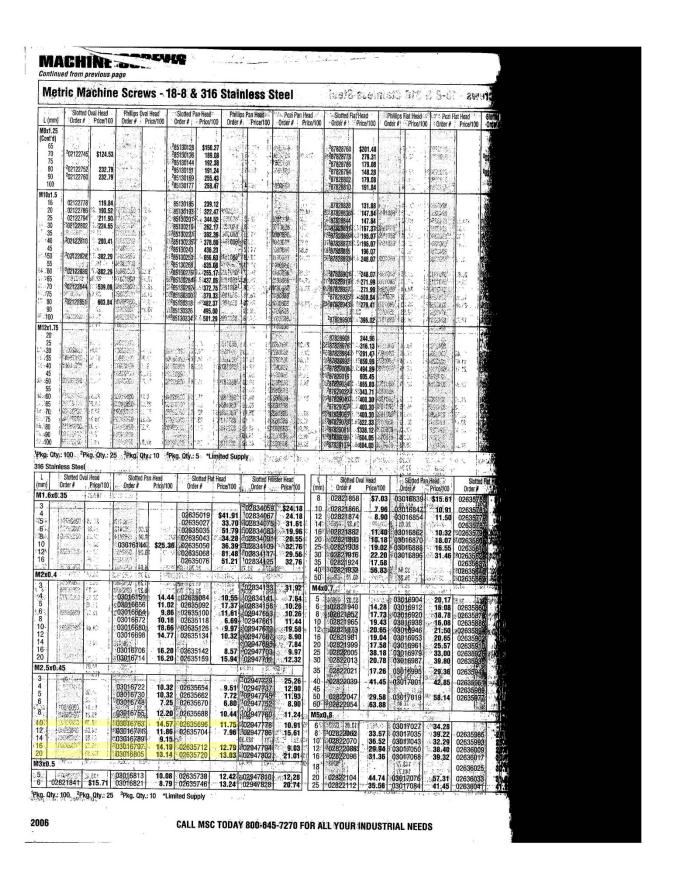

#### Screw Nuts - Military Specification 35649 / 35650

Price/100

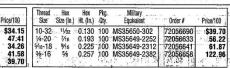

Material: steel Grade: 60M Finish: cadmium yellow plated



(III) IMPORT

- Charles	EJRO.	15	157	1		oc.	1		300	200	80.	 Styric's	artist.	521
j	Triu.	* 1	1.7		- 1	100	**	10			4	55.	;	

ALABATATOT LOS INTERNAS LOS SERVICIOS AS A SECULIAR AS A S

Çŧ .	c Plated Price/100	Class & Pla Order #	ain Steel Price/100	Class 10 P	fain Steel Price/100	18-8 Stain Order #	ess Steel Price/100	316 Stainless Steel Order # Price/100		
550 50 90	\$2.72 2.97 1.85 3.07	09170424 09170432 67476960 09170457	\$2.62 2.83 1.21 2.89	[ ] [ ] [ ] [ ] [ ] [ ] [ ] [ ] [ ] [ ]		68024009 68024041 68024082	\$6.12 7.30 4.93	<sup>2</sup> 62323555 <sup>2</sup> 62323563 <sup>2</sup> 65093940 <sup>2</sup> 65093957	\$11.07 11.12 7.60 14.56	
81 84 84	1.86 1.85 2.11 4.39	67477000 67477042 67477083 9 01010693	1.39 1.63 1.86 4.25	01010339 67477406	\$4.59 5.38	68024124 68024165 68024207 381707804	6.68 6.91 8.87 30.67	65093965 65093973 65093981	8.87 9.31 10.27	
48 89 21	4.74 8.17 12.90	67477125 67477166 67477208	3.72 6.84 9.59	67477448 67477489 67477521	9.42 18.83 26.89	68024249 68024280 68024322	23.85 43.10 60.77	65093999 65094005 65094013	28.83 55.19 80.19	
-	Price Ea.		Price Ea.	-44	Price Ea.	127	Price Ea.		Price Ea.	
82 04 46 87	0.25 0.22 0.49 0.58	67477240 67477281 67477323	0.15 0.39 0.58	67477562 67477604 67477646 67477687	0.39 0.52 1.25 0.96	68024363 68024405 68024447 68024488	1.03 1.28 2.90 2.93	65094021 65094039 65094047 65094054	2.07 2.12 3.12 3.32	
73 19 17	0.92 1.21 1.99 2.31	09170440 67477364 81195430 08018848	0.65 1.17 1.87 2.18	09170515 67477729 09170523 07366826	0.85 1.95 2.34 2.85	81707887 68024520 81707903 81707911	3.73 4.77 15.69 13.34	65094062 65094070 65094088 65094096	3.68 7.2 13.61 22.09	
		09170465 05736798	2.49 2.86	09170531 09170556	3.16 3.42	81707929 81707937	34,25 23.11	65094104 65094112 65094120 65094138	25,96 36,69 43,11 44,44	
	100 EV.	S10586 5	E de E	10000 3740835 i	Comments of	4)660 - 388448 - 388448	10 A 4 A 7 A 1	65094146 65094153 65094161	45.09 .96.02 .117.17	



Screw W. Across Ht.	Screw W. Across Ht.
Size Flats (mm) (mm	Size Flats (mm) (mm)
M2 4.0 1.6	M12 19.0 10.0
M2.5 5.0 2.0	M14 22.0 11.0
M3 5.5 2.4	M16 24.0 13.0
M4 7.0 3.2	M18 27.0 15.0
M5 8.0 4.0 M6 10.0 5.0 M7 11.0 5.5 M8 13.0 6.5 M10 17.0 8.0	M20 30.0 16.0 M24 36.0 19.0 M30 45.0 25.0 M36 54.0 30.0

Thread Size Pkg. Qty.	3 (3 Order #	Price/100
M8x1 100	67488643	\$15,57
M10x1.25 100	67488684	29.45
M12x1.25 100	67488726	41.44
This is a company B.	Stafnless Stol	Price Ea!
M14x1.5 1	67488767	0.67
M16x1.5 1	67488809	0.77
M18x1.5 8	67488841	0.85
M20x1.5 1	67488882	1.07

(III) IMPORT

. Military. Equivalent

MS35649-222 MS35649-242 MS35649-262 MS35649-282 MS35649-202

13 1 14 16 ST

89,46 52,30

teel with a bright finish of zinc plating for rust and corrosion resistance.

gth, for use with Grade 2 fasteners,
100se grade of low carbon steel with no finish unless a grade of nuts is
2, 5 & 18.

101 Element of the finish of zinc plating for rust and corrosion resistance to cracking and metal fatigue. Corrosion-resistant gold-colored coating.

102 Farage 8 Steel - High strength, use with Grade 8 steelers.

103 Farage 8 Steeler - High strength, use with Grade 9 screws and washers for the highest available strength and resistance to cracking and metal fatigue. Corrosion-resistant gold-colored coating.

103 Farage 8 Steeler - High strength ruse with Grade 9 screws and washers for the highest available strength and resistance to cracking and metal fatigue. Corrosion-resistant gold-colored coating.

103 Farage 8 Steeler - High strength ruse with Grade 9 screws and washers for the highest available strength and resistance to cracking and metal fatigue. Corrosion-resistant gold-colored coating.

103 Farage 8 Steeler - High strength ruse with Grade 9 screws and washers for the highest available strength and resistance to cracking and metal fatigue. Corrosion-resistant gold-colored coating.

104 Farage 8 Steeler - High strength ruse with Grade 9 screws and washers for the highest available strength and resistance to cracking and metal fatigue. Corrosion-resistant gold-colored coating.

Bold in Pkg. Qty. Only

no Plated Price/100	Grade 2 Plai Order#	n Steel Price/100	Left Hand Pla Order#	in Steel Price/100	Grade 5 Zini Order#	c Plated Price/100	Grade 5 Pla Order #	in Steel Price/100	Grade 8 Yell Dichron Order #	ow Zinc (182 nate Price/100	Grade 8 P Order #	lain Steel Price/100	L9 Yellow Zinc C Order #	Dichromate Price/100
\$1.70 1.91 2.60 2.67	67473728 67474601 67473769 67474643	\$1.61 1.77 2.60 2.75	67475442 67475962 67475483 67476002	9.23	67471086	\$2.51 2.84 3.82 3.87	52593324 52593332 52593340 52593357	\$1.56 2.45 2.97 3.74	05724406 85724409 05724414 85724417	\$3.04 4.01 4.13 4.50	67470245 67472126 67470286 67472167	4.08 5.07	67512905 03835923 67512913 03835931	
3.59 3.93 7.19 7.47	67473801 67474684 67473843 67474726	3.58 3.79 6.07 6.88	67475525 67476044 -67475566 <sup>3</sup> 67476085	14.30 15.85 30.11 30.12	67471128 67473009 67471169 67473041	5.06 5.57 8.43 9.71	52593365 52593373 52593381 352593399	3.05 4.90 5.37 9.43	05724422 85724425 05724430 85724433	6.09 6.80 11.69 14.02	67470328 67472209 67470369 367472241	9.37 7.69 12.59 17.98		36.36 37.18 53.91 54.46
0.03 10.36 17.12 17.36	<sup>3</sup> 67473884 <sup>3</sup> 67474767 67473926 -67474809	6.41 7.97 11.66 13.61	367475608 367476127 67476640 167476168	75.13		12.43 12.82 20.20 22.79	352593407 352593415 52593423 52593431	7,27 5.80 12.41 28.26	305724448 385724441 05724455 85724458	15:13 30.91	367470401 967472282 67470443 167472324		267513002 267513366 67513044 367513408	61.17 62.11 73.72 74.01
0.19 0.20 10.28	01 67473967 67474841 67474007 67474882	0.18	167475681 167476200 267476723 267476242	3 0.80	67471284 67473165 67471026 67473207	Price Ea. 0.21 0.24 0.34 0.36	52593449 52593456 52593464 52593472	Price Ea. 0.13 0.20 0.38 0.40	05724463 85724466 05724471 85724474	0.36 0.43	\$67470484 \$67472365 \$67470526 \$67472407	0.39 0.53	007 67513085 67513440 67513127 67513481	1.24 1.28 2.13 2.26
0.47 0.57 0.78 0.84	67474049 67474924 67474080 67474965	0.38 0.68	67475764 267476283 267475806 267476325	1.99 2.63		0.55 0.63 0.97 0.97	452593480 52593498 252593225 52593233	0.94 0.42	467470807 67472688 67470849 67472720	0.81 1.14	067470567 967472449 267470609 267472480	0.88 1.47	67613168 67513523 267613200 67613564	4.32 4.40 6.71 6.82
1.13 1.39	67474122 67475004	1.04 1.11	67475848 567476366	9.80 7.12	<sup>4</sup> 67471441 52592276	3.51 1.40	452593241 52593258	0.98	67470880	199-2.26	267470641	2.12	67513242	20.09 20.09
	CONTRACTOR STATE		°67476366 Pkg. Qty. 1 <sup>6</sup> P				1.0		67472761	Jak .	7387	Sugar		31-65

READY TO ORDER? - mscdirect.com

2023

#### WASHERS \* Flat Washers - Military Specification - are rest Stainless Steel Isels before more USA Manufactured from:18-8 stainless steel, these washers are burr-free, and passivated to meet stringent military specifications. Conform to MS, AN or NAS chemical and; physical specifications. MS washers have larger ID and OD's and are thicker than AN, and MAS. MAS washers - The "L" in the Dash Number column refers to the light (thinner) series of washers - Tolerance: +0081/-005" Screw Size Dash Number ISS. Min. Thick. (In.) (In.) (in.) MS15795 Military Specification 0 801 0.078 0.188 2 802 0.094 0.250 0.025 0,025 0,028 0.016 0.016 0.017 0.025 0.250 0.040 805 0.156 0.156 0.188 0.188 0.312 0.027 0.048 0.065 0.065 806 807 841 0.036 0.382 0.036

0.445 0.507 0.579 0.632

0.739

0.693

0.897

1.022

1.084

1.397 1.334 1.772 1.491

2.022 1.772 2.272

2.022

0.262

0.262

0.330

0.387

0.387 0.387 0.450 0.450

0.512 0.512

0.574 0.574

0.637 0.637 0.762 0.762

0.895 0.895

1.074

1,324 1,512 1,762

0.187

0.245

0.282

0.282 0.311 0.311

0.361

0.361

0.475

Aeronautics Standard 0.065 0.106 0.091 0.156 -0.104 0.187

0.036

0.036 0.051 0.051

0.051 0.051

0:064

0:064 0.051 0.064 0.074

0,086

0.122 0.108 0.136

0.108

0.014 0.028 0.014 0.028

0.014

0.014

0.014 0.028 0.028 0.057

0.028

0.057

0.057

0.028

0.028

0.028 0.057 0.028 0.057

0.082

0.014 0.014 0.014 0.029

0.014

0.014

0.014

0.029 0.0585 0.029 0.0585

0.219

0.219 0.257 0.296

0.327

0.351

0.382

0.445 0.476 0.507

0.538

0:569

0.678

0.834

0.834 0.960 0.960

1.084

0.127

0.151 0.151

0.176 0.176 0.205 0.205

0.267

0.330

0.392 0.402 0.455 0.455

0.525 0.525 0.580 0.642

0.767 0.892 1.017

0.117

0.130

0.145

0.145 0.171 0.171

0.197 0.197

0.257

815 816 817

820 821

822

823 824 825

826 827

C4L 0

C6L C6 C8L C8 C10L C10 8L 8

C416L C416 C516L C516

C616L C616 C716L C716

C816L C816

C916 C1016

C1216 C1416 C1616

NAS620 National

C0 C2 C3L C3

C4L C4 C5L C5

C6L C6

C10L C10

C416L C416

2036 /= New Item!

1/4

5/16L 5/16L

3/8L 3/8 7/16L 7/16

1/2 L 1/2 9/16

3/4

31

4L 5L 5

6L 8L 8

10L 10

1/4L 1/4

AN960 Aeronautics Specifications
2L C2L 0.101 0.266
2 0.20 0.101 0.266
3L 63L 0.107 0.266
3 63 0.107 0.266

5/16 5/16 3/8 812 813 814

3/8 7/16 7/16 1/2 818

1/2 5/8 5/8 3/4 819

3/4 7/8 7/8



F436 Structura	al Wash	ers		
Made to ASTM F436 spi Designed for use with A				
Screw OD ID Thick. (In	n.) Pkg.	Plain Steel	क्षान्त्र	Galvar
the state of the s	80 100 078	12142	\$5.15 470	8415 8416
3/8 13/16 13/32 051 .0 1/2 11/16 17/32 097 .1	80 100 078		7.38 770	8417 8418
5/8 15/16 11/16 122 :1 3/4 115/32 13/16 122 :1 3/8 13/4 15/16 136 :1	77 25 078	12209	29.10 7702	8419 8420
1 2 11/8 .136 .1	77 10 078	12126	53.09 7702	
11/8 21/4 11/4 13/6 .1 11/4 21/2 13/6 13/6 .1 11/2 3 15/8 .13/6 .1	77 10 078 77 5 078 77 5 078	12175	91.48 7702 94.19 7702 82.04 7702	8424
12 3 178 1130 11	1181 107 1181 107 1181 178	05 110.		.04202
Flat Washers	Extra	Thick	0.620	ċ2;
Bolt Size Pkg. (In.) Qty. Order#		It Size Pkg. (In.) Oty.	Order.#	Price
100 67492447 5/16 100 67492488	\$19.25	% 50 % 50	6749264 6749268	\$98
% 100 67492520 % 100 67492561	33.29 35.12	¾ 50 % 25	6749272 6749276	8 239 408
100 <b>67492603</b>	61.36 01a	1 25 <sub>0c</sub> 1 0 0 0	6749280	5 <b>82</b> 088
Metric Flat Wa		galili Kunio		4
Corrosion Res			and Latination	
High visibility blue color Metric Blue coating is 40 Bolt 0 00	)x more corro	sion resis	stant than s	andar
Bott ID QD Size (mm) (mm) M6 7.4 12.0	(mm) Oth	. TC On	ler# Pri	ce/f00/4 28.70
M8 9.6 17.0 M10 12.0 2105	3.50 10 4.50 10	0 -792	07726 07734 08028	29:85 36.88
M12 15.0 27.0 M16 18:4 34.0	4.60 10	798	08036	53,7 99,28
M20 22.5 42,5 M24 26,5 50,0 M30 33.6 60.0	4,60 50 4,60 25 4,60 20	798	08944 08051 08069	03.0 29.0 96.4
17.20   1881:73.10)	2.037-0.07 2.037-0.07 1.038-6.60	341	60.37 3	300
Metric Flat Wa	projection and thereto and in all	200 Pc	9 19.10 2 24.57 2 24.93	
50 1775511 48.31	rue sulo rue sulo	1	2017	3,50
50 407 16:201   43:22 	673-0 Act 9 2305-4864 676-4-686	((0	) 1600	22 37
7 - DE TOWN THE OF	976.0 320.0 66.0-760.0 66.0-772.5	IK-/X		100
Bolt ID OO Thick Pkg. Size (mm) (mm) (mm) Qty.	Order #	Price/100	18-8 Stal	
M2 2.2 5.0 0.3 100 M2.5 2.7 6.0 0.5 100	09063496 09170358 67492843	\$1.52 1.56	6802492 6802496	1
M3.5 3,7 8.0 0.5 100	09170882	0.91	6802500 90355081	2
M4 43 9.0 0.8 100 M5 5.3 10.0 1.0 100 M6 6.4 12.5 1.6 400	67492884 67492926 67492967	1.16 1.35	6802504) =6802508 ::6802512	V
M7 7.4 14.0 1.6 100 M8 8.4 16.0 1.6 100	67493007 Unlist 9	2.58	90917034 11 19110 31	12
M8 8.4 17.0 1.6 100 M10 10.5 20.0 2.0 100	67493049 019	2.74	6802516 10	1
M10 10.5 21.0 2.0 100 M12 > 13.0 24.0 2.5 100	67493080 67493122	4,36 8.22	6802520 6802524	000
M14 15.0 28.0 2.5 100 M16 17.0 30.0 3.0 100 M18 19.0 34.0 3.0 100	67498163 67498205 67498207	10.45	6802628 6802638	O O
	67493288	20.01	6802536 6802540	no n
M20 21,0 37,0 3.0 50 M22 23.0 39,00 3.0 50 M24 25,0 44,0 4.0 25 M27 28.0 50,00 4.0 25	67498320	23.23 25.55 59.26	0355092 168025 <b>44</b> 103550944	500 600 600
M30 31.0 56.0 4.0 25 M33 34.0 60.00 5.0 10-	67493361	72.04	503550951 -08550986	1
M36 37.0 66.00 5.0 5 M39 40.0 72.00 6.0 10	09170390	117.35	*03550986 *03550977 *03550986	.4
M45 46.0 85.0 7.0 5	A Second Trust	CAMPS 11		187


Flat Head Slotted Flat Head Phillins 1 Round Head College Box QUANTITIES ONLY - NOT SOLD IN BROKEN PACKAGES	erclassic wars beyer
Order #   100   Order #   10	Fillister Head Slotted Truss Head Slotted Thomas
40915423 \$19.05 40914632 \$8.90 40918443 \$4.80 appropriate 40000000 40000000 400000000 400000000 4000000	Order # 100 Order # 100 Min Africa
9317 3238 6.25 40914657 6.20 87909503 3.60 07112980 8.33 67409680 5.00 67413328 6.21 40915449 5.60 4091649 5.	x87419242 1 10.11 chen muni
40914681 5.44 40918476 4.89 7.13003 5.74 67409768 4.70 67413401 3.81 40915456 5.02 40914699 5.52 87909545 3.80 67413901 67409805 7.04 67413403 15.17	67.419366 10.01 (# 50ptor : 5 - 5 - 5 - 5 - 5 - 5 - 5 - 5 - 5 - 5
40915472 8.24 40914715 5.96 87909560 3.81 07113020 5.29 67409828 3.81 07413894 5.97 UNC	67419449 16.81 stp. it it it is in a 160 state of 160 sta
67391961 3.88 40914723 8.85 40916492 5.11 00 mag 40920050 4.65 4092699 5.85 7399901 7.81 40914731 7.53 87909586 5.89 0713090 40920050 4.65 4092699 5.85 73999586 5.89 0713090 40920050 4.65 4092699 5.86 73999586 5.89 0713090 40920050 4.65 4092699 5.86 73999586 5.89 0713090 40920050 4.65 4092699 5.86 73999586 5.89 0713090 40920050 4.65 4092699 5.86 73999586 5.89 0713090 40920050 4.65 4092699 5.86 73999586 5.89 0713090 40920050 4.65 4092699 5.86 73999586 5.89 0713090 40920050 4.65 4092699 5.86 73999586 5.89 0713090 40920050 4.65 4092699 5.86 73999586 5.80 0713090 40920050 4.65 4092699 5.86 73999586 5.80 0713090 40920050 4.65 4092699 5.86 73999586 5.80 0713090 40920050 4.65 40920000 4.65 40920000 4.65 40920000 4.65 40920000 4.65 40920000 4.65 40920000 4.65 40920000 4.65 40920000 4.65 409200000000000000000000000000000000000	6/419623 10.11
67392084 4.57 40914756; 7.85 187909594 5.28 40918567 6.39 40920075 5.01 40920085 67392126 4.85 40914764 6.08 8790950 5.44 67302126 4.55 40914764 6.08 8790950 5.11 6740906 674	10923633 6.26 10923641 5.43
67392209 611 872 872 87909628 4.68 071 3046 4.49 674 0001 5.83 40920860 5.49 67392209 611 879 096966 514 071 3046 4.49 674 0001 5.83 40920876 6.05 44	10023666 6.26 34.573 3555 400
0.5322282 8.19 37303651 9.25 0713061 7.39 00920994 9.59 75392324 5.23 00920994 9.59 75 00920996 5.37 00920996 5.50 75 00920999 5.50 75 0092099 5.50 75 0092099 5.50 75 0092099 5.50 75 0092099 5.50 75 0092099 5.50 75 0092099 5.50 75 0092099 5.50 75 0092099 5.50 75 009200 5.50 75 009200 5.50 7	25 SA CO. 7
7/392465 9.60 9.60 9.60 9.60 9.60 9.60 9.60 9.60	to receive a superior and a superior
7392409 278 67399462 335 5690541 129 40918623 346 674 0084 238 674 308 232 447 7392400 456 67390204 327 6790593 238 674 0084 238 674 0086 158 674 00	1788020 6.98 1419664 1.58 1419666 5.54 08704702 \$6.01 674
7392522 8.16 40914798 5.37 87909749 3.85 87913927 3.75 67410209 1.55 6791330 1.56 6791300 1.55 6791330 1.56 6791300 1.55 6	173188 6.99 419689 - 5.99 40923369 - 6.34 674
7911/31 4.1 67399369 7.51 87909750 1.82 071.13087 4.35 67410281 1.75 87913331 1.87 67392563 8.53 67399409 8.3 67909768 2.11 87913968 4.8 67410282 1.72 87913399 1.98 678	419/21 3.88 40923377 6.96 6741 419/62 5.54 6741
392605 7.00 40914806 7.80 87909784 2.34 0713403 6.00 67410407 216 87913363 3.34 674 392647 6.76 0711263 7.87 87909782 5.22 0711317 13.71 67410408 6.00 67410407 2.66 87913372 2.64 674	19946 12.34 13 88919 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1
392688 28.97 2.01 30920118 15.38 09173881 11.49 328	23682 14.22 52287 18.16 20324 20324 20324 20326 20326
392761 5.25 40914814 7.15 67401967 4.27 03958535 5.21 40920126 5.25 40920936 5.64 4093863 5.14 40914862 7.23 67401968 7.33 04412465 6.49 67419480 2.43 40920944 7.71 674	23690 9.23 4092 <b>943</b>
92699 645 99173329 6.12 67402008 5.84 67402040 9.08 67410522 6.80 40920951 6.94 67410523 7.09 40920969 7.44 6742	409231000
92969 4.75 0,402081 4.56 03958584 5.88 67410605 6.80 40920985 10.95 6742 4.83 4002093 6.80 40920985 10.95 6742	0125 32.40 00 32.4 33.4
33082 7.96 4094838 431 67402248 4.39 07113137 18.20 40920167 8.3 40921075 5.37 6.71	0166 12.48 interest states and the other states are states are states and the other states are states are states are states are states are states are states and the other states are s
9246 12.49 67402321 7.99 40918625 12.06 40920183 12.06 40921041 10.85 horses	2°78/
3207 2.80 365165615 9.80 87909925 4.66 40920191 9.86 40920392 9.7009000	Technology Control
3280 2.22 67399527 3.21 87909958 1.77 87913984 3.66 674/0720 1.89 674/20 3322 1.94 67399528 3.48 87909958 1.77 87913992 3.66 674/0720 2.13 87913982 2.12 674/20	240 14.45 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1
305 2.13 67399600 7.22 87909974 1.12 87914004 2.17 67410761 1.87 87913414 2.55 6874206 3.18 6739962 2.17 87909982 1.89 87914024 2.18 674208 2.54 87913422 2.88 674208 2.54 8791342 2.54 879134 2.54 8791342 2.54 879134 2.54 879134 2.54 879134 2.54 879134 2.54 879134 2	507 3.56 67416644 2.77 6741866 17 233 2.86 67416685 1.48 674184 7 364 7.39 67416727 2.83 2.74 2.74 2.74 2.74 2.74 2.74 2.74 2.74
822 2.15 67399766 -10.11 87910014 2.57 87914032 2.76 6741086 2.09 87913430 2.6 674204	06 4.78 1.67416768 -2.78 167415768 47 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1
355 3.10 67399840 4.35 87910022 2.34 87914065 3.90 67411099 2.50 87913463 3.34 674205 3.34 6741041 2.51 87913463 3.34 674205	21 12 48 87 44 0044 34 0 044 0 0 0 0 0 0
321         67399881         5.9         67910048         3.85         67910073         5.8         67411041         2.85         67413724         3.55         674206           5.55         40914848         13.71         67910055         3.85         87914095         14.24         67411922         3.23         67413724         3.55         674206           9.95         3.85         86809078         4.24         87910063         4.75         40918641         14.24         67411324         6.11         40924044         10.35         69143124         10.35         67411165         4.89         108047862         4.75         40918641         13.53         67411165         4.89         108047862         3.56         674206           7         5.31         76824434         10.11         67402446         5.81         40918658         27.35         409202271         12.9         40921082         0.02         0.02           8         6.43         6740246         5.81         40918658         27.35         40920217         12.9         40921082         0.02         0.02	62 6.86 67416925 3.70 67415986 4.0923385 9.89 40923245 7.00 67415886 4.0923385 9.89 40923245 7.00 67415886 4.0923385 9.89 40923245 7.00 67415886 4.00 674158
00 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0	24.8
20 8.98 78624616 19.98 67402529 8.94 480913666 39.22 40920225 \$3922 09173907 27.47 3941 15.618 153 14.99 40914863 61.86 1574025602 12.89 40914863 48.80 12.80	1 12909233931 4 1172.8 1 1271 137
71 17.63 409 4871 53.93 67402665 10 74 409 650 7 7 7 8 8 409 409 14871 53.93 67402665 10 74 409 650 7 7 7 8 8 8 8 8 8 8 8 8 8 8 8 8 8 8 8	2 059018 6 87, 2 1
3 7.55 5 67402727 2.44 40948660 6.40 67411207 3.11 40021461 7.0 1200000000000000000000000000000000000	35   18.61   18.62   1
	B4E 0.3
CALL MSC TODAY 800-645-7270 FOR ALL YOUR INDUSTRIAL	On 8 Continue
	1.9


Zinc	Plat	ed Ste	el		***						iois.	10.199 sall Ta	SME		-	IMPOR
Head Phillips Pri er# 0010	le/		tted Price/ 100		rice/		tted Price/ 100		lips State Price/	Fillister Head S Order#	Price/ 100		Price/ 100	Binding He	Price/	in the same of the
99964 00002 00044 00085	3.45 6 5.88 6 3,52 8 3.80 8	7402800 7402842 7402883 7910188 7910196 7910204	2;11 2,87 2,15 2,84	87914115 87914123 40918708 87914149 87914156 87914164	2.96 6 4.03 6 3.02 6 3.29 6	7411280 7411322 7411363 7411405 7411447 7411488	2.28 3.40 2.46 3.08	67413849 67413880 67413922 87913505 87913547 87913554	2.75 4.25 2.82 3,62 3,32	67420760 67420802 67420844 67420885 67420927 67420968	4:76 8.99 4.08 6.28 4.32	67417006 67417048 67417089 67417121 67417162 67417204	2,79 4,16 3,19 4,33 4,19	6741596 6741600 4092329 6741604 6741608	3.81 04 6.24 00 3.81 31 4.59 23 4.64	
00168 00200 00242 00283 78337	4.30 8 4.34 8 4.98 8 6.26 8 15.24 8	7910212 87910220 87910238 87910246 87910253 87910264	3.42 2.84 3.05 3.18 5.67 3.73	87914172 87914180 87914198 87914206 67407569 87914214	3,86 6 5.06 6 6.04 6 8,62 6 9,63 6	7411645 7411686 7411728	3.70 4.74 4.84 6.86 7.47	67413963 87913562 67414003 87913570 74084740 08991101 40921165	3.94 6.02 6.22 4.49	67421008 67421040 67421081 67421123 67421164 67421206	6.83 8.27 8.99 23.78 19.10	67417246 67417287 67417329 67417360 09173493 53464327	4.42	6741616 6741620 409233 409233 409233	06 5.38 02 13.38 10 16.55 28 19.59	0.0000 0.000 0.000 0.000 0.000
781 14905 14756 14913	9,49 26,84 10.48 45.25 69.13	87910287 87910287 87910287 87910295 87910303 67402966 67402966	7,90 8.17. 9.43 8.78 12.42 14.20	40918716 40918724 40918732 40918740 40918757 40918765	27,79 C 30.74 4 34.16 4 61.00 4 63.48	0920266 09173725 10920274 10920282 10920290 10920308	7.49 31.65 26.85 53.94	09173915 40921173 40921181 40921199 40921207 40921215	9,49 37,48 18,03 66,96 69,13	40923740 40928757 40923765	20,25 59,02	09178501	5	674.174 674.07 6740.03 6740.03 6740.04	8 81 1	15(7 . 8
N. 5		67403048 09173584	82.84	77 674228 40 874228 81 6742 21 6742	8 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1	09173733	7.49	40921223 4092128 6741404	86,81,	20917320	2 58	UF ICIO	ner i	40923	336 4.6	17918 5
400366 400408	4.35 10.27 4.37	67403089 67403123 40918500 67403162 87910386	2,54 2,54 2,77 2,50	04312401 87914255 87914263	5,58 4,16 12,99 3,82	67411769 67411801 67411643 67411884 67411926 6741196	3.56 3.08 7.17 3.07	67414080 67414121 8791360 8791364	4,08 4,62 4,61 3,86	4092377 6742124 4092378 6742128	3 10,35 B 5,62 B 9,14 9 5,01	06741740 06741744 06741748 6741752	3.7 4 12.7 5 3.9 7 12.4.4	5 67416	344 5.0 248 4.8	0 216.3 2021/058 0 22/058 0 24/04/05
400440 400481 1,73352 400523 400564 400606	3.94 4.56 10.24 5.28 5.91 6,93	87910394 87910402 87910410 87910428 87910436 87910444	2.80 4.57 3.15 3.79	87914305 87914313 87914321 87914339 87914347	4,31 12,36 5,06 6,92 7,02	6741200 6741204 6741208 6741212 6741216	5.30 5.30 4.59 5.00 6.19	6741416 6741416 8791366 0804795 8791367	3 4,47 9 7,54 9 5,82 3 4,61 9 7,60	6742144 6742148 6742152	4 23.94 6 8.16 7 8.22 9 12.5	6741760 6741764 6741768 674177	0 8,4 2 6,3 3 7,4 5 8,5	9 17: 1	51 9 2 - 5 4 - 2 4	36.2 01 5170 446 13. 5178 1385 999
173360 188876 1914989 1298438 1914947 1924830	10.24 10,24 32.52 10.48 31.27 10.99	87910451 87910469 8791047 87910489 40918511 8791049	9 6.64 7 7.94 5 8.5 8 7.0	879,14362 4 409,1877 5 409,1878 6 409,1879	10,47 24,74 26,67 41,92	6741224 4092031 4092032 4092033 4092034	7 9.15 6 27.90 4 26.90 2 39.90	8791369 4092124 0804798 4092125 080479	5 9:20 9 30:63 7 7:24 66 41:92 95 8:78	6742160 6742164 6742168 6742176	16.4 4 26.8 35 37.8	6741780 4 091735 9 091735 6 091735	15 9 11 7 11 35 11	11 24 49	5. 13 1. 5. 19 . 0 5. 14 . 1 6. 12 . 14 . 15	0.700 (4.00) 67.00 (57.00) 57.30 (50) 677.90 (50)
9173378 0914954 0914962 0914970 0914988 0914996	42,58 16,28 70,44 76,17 69,45 77,24	6740324 8791050 8791051 8791052 8791053 8791054 8791055	1 13.7 9 13.4 7 17.9 5 20.8 3 40.0	1 4091881 3 4091882 4 4091883 3 4091884 77 4091885	5 38 12 3 44 29 1 49 94 9 50 60 6 76 85	4092036 4092037	5 75.7 3 75.2 31 95.4	409212 409212 409212	72 66,4 80 70.4 98 76,1	i 4 7	gr¦* .	820665		41	12.15	67364881 0300 0301786 0371786 0371786
9173345 7400648 7400689 7400762 7400762 740080	9.55 4.77 8.84 4.20 4.31	8791057 8791058 18791059 879106	6 3. 34 2. 92 3. 18 2.	35   0711317 37   8791439 28   8791440 56   8791441 84   8791442	0 5.13 8 9.21 6 3 50 4 12.9 2 3.7 0 3.9	674122 674123 674123 674124 674124 674124	88 3. 20 4. 61 3. 03 6. 45 3. 86 3.	3 8 879137 8 879137	26 3. 37 4. 45 4	6 674218 674218 674219 3 674219 16 674220	42 23. 83 6, 925 19. 966 7,	40 409284 04 674178 49 674178	27 9 40 3 81 28	72 6741 550 6741	6362 4 6404 6 6446 4 6487 1	279 5 6 6 6 6 6 6 6 6 6 6 6 6 6 6 6 6 6 6
7400846 740088 7400929 7400960 740100	5.3 5.4 7.2	879106 8 879106 8 879106 4 879106	42 4. 59 3. 67. 4. 75 4	90 879144 45 879144 39 879144 08 879144 53 879144 30 674076	46 8.2 53 4.5 54 6.9 79 7.7	1 674125	69 5. 01 4. 43 6. 84 7	03 674144 51 879137 23 674144 33 87913 95 87913	109 6. 160 5. 141 7. 178 8. 186 11.	98 674220 78 67422 69 67422 59 67422 38 67422	089 8 121 9 162 11 204 13 246 13	70 40923 00 67418 41 67418 69 67418 62	135 1 004 046 087	3.02 Oare 6.94 674 7.38 9.32	6560	6.59 68 65 65 6 10 5 10 8 10 5 10 8 10 6 6 6 6 6 6 6 6 6 6 6 6 6 6 6 6 6 6 6
8740108 4091500 4091504 4091502 4091503	3 10.7 1 33.0 9 15.5 7 44.5 5 16.9	1 879106 7 879107 8 879107 2 409185 6 879107	91 5 09 8 17 8 84 10 725 9	.45 879144 .53 409188 .90 409188 .99 .83 409188 .29 409188	64 30.7 72 33.4 80 35		399 31 407 73 415 44 428 10	11 -40921 90 40921 24 08048 40 40921	314 33 322 13 330 45 068 8 348 62	22 40923 07 40923 00 40923 99 40923 53	799 36 807 36 815 27 823 3	47 40923	443 1	1;19  6,54    31.19	17 9 468 17 12 17 15 18 15 18 15	60000008181 6   160116181 6   18012016 6   18081218
4091505	0 78.	75 879107 674032 674033 674033	744 - 20 287 - 20 329 - 23	0.00 409189 0.74 409189 3.41 2.05	06 69. 114 77.	66 40920 85 argen	449 78	1 Aus 1	363 66 371 81	SKL81		574045 574045 77645	32 32	i fart	0 .0 0 40 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0	300000 31000000 31000000 810000000000000
	25 -10.	67403 67403 85 67403 67408 67 87910	402 444 485 527 1 774	5.25 C409189 5.76 0 4.83 -67407 6.55 4.86 67407 6.98 67407	643 7 684 15 726 10	39 40920 52 40920 76 09173 87 09173 91 40920	464 1 74 6 1	4.36 74092 4.36 7408 4.74 0917 4.53 8813	397 1 4708 3931 1	1.39 4092 5.52 4092 9.12 6742 1.74 6742 0.61 6742	3849 2329	1:30 0:58 4092 16:86 12:50 4092	3468	13,85	10 11 1 45 1 6 100 - 1	379131051.2


	Machine  L Under Head (In.) Order #	Price/	Ordon / Price/	lound Head Phillips Price/	Pan Head Slotted	Pan Head Phillips	Fillister Head Slotted	Truss Head Slotte		
	#12-24 UNC (Cor 78 87912556 1 87912564	1'd) \$15.48 67401284 \$16.79 8.24 67401323 16.02 10.70 67401364 12.91		0rder #	Order# 100	Order # Price/	Order # 100	Order #	De/	
	11/4   87912572 11/2   87912580 13/4	10.79 97401364 12.91 11.54 97401364 22.43 340915068 27.69 12.52 49915076 33,35	87910816 6.53 67 87910824 874 667 87910832 8.01 867	407882   21 as la	0000000	09173964 15.86	67422444 \$42.99 67422485 16.53	350170 35	24e (* 2.1.)	
	21/2   87010000	1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1	87910857 garan E	18948 33.35 4	0920514 30.74	91837193 16.36 40921498 25.92 40921418 35.73	67422568 34.38 67422660 24.43 67422669 44.3	82 879101 82 879101 89 879101	GCOW A D GCOME O D GCOME A	
	0110	38:18 34:0915092 40:65 26:46 34:0915090 51:03	87910865 14.99 0409 87910873 16.34 0409	18963 51.03 CE	1218 10	40921424 41.23 40921489 51.03	141 378	95 875102 135 875102 24 875102	100 Ve 4	
and the same of th	9/16 87912655 36 87912663	22.11 200000018	87910907 6.52 677	13196 1474 147	18 to 160° pt 150			98 879102	S0085 <b>8</b>	
illusizonati	7/16 87912671 1/2 87912689 5/8 87912697	18.39	7910923 3.64 879 7910981 9.20 7910956 3.68 8791 7910964 5.57 8791	8971 21.04 40 4503 5.03 67 4514 5.01 67 4529 514 67 4537 7/50 67	112809 5.50 °0	0921447 20,17 4 0921454 11,47 9173880 4,61 4 0921462 16,12 4 7913836 6,73 6 7913844 8,28 6	ARRAGUET	40923494 <b>\$23.6</b> 40923492 24.9 67418160 6,90		
think Circum	% 87912705 % 87912713 1 87912721	7.41 -67401604 11.07 8	7910972 4.22 8791	4529 5 14 67 4537 7 158 674 7924 9 93 674 4545 8 16 674 4552 9 84 674	12882 6.58 8 12924 6.65 8	/913886 6.73 6 7913844 8.28 6 7913854 7145 6	(422725) 8.61	67418202 6.01 67418244 8.51	8-1-20 14-20-5	
	11/4 87912739 11/2 87912747	5.80 67401687 8.99 8 7.20 67401729 10.46 8	7910998 4.75 8791 7911004 4.74 8791 7911012 5.80 8791	4545 8.16 674 4552 9.84 674 4560 10.48 674	12924 - 6.65 - 8, 12965 - 12.88 - 6, 13005 - 7.79 - 87 13047 - 8.95 - 67	414482 9.80 67 913869 10.27 67 414524 12.47 67	422840 28.34	7418285 8.23 7418327 16.25 7418368 9.41	CATCLE	
	2 87912762 14 87912770 1	9.68 .67401802 12:30 87 3.13 67401844 19:98 87 8.87 307312469 17:18	911099 740 07407	965 11.49 674 578 12,21 674	13005 7.79 87 13047 8.95 67 13088 19.26 87 13120 14.85 87 3161 13.37 67	913877   12.85   -67 913885   16.12   67 414565   18.90   67	422881 11.92 422923 16.21 6 422964 50.95 6	7418400 12,02 7418442 14.46 7418483 12.13		
ŧ I.	87912812 14 40915506 1	3.13 67401844 19.98 87 8.87 307312469 17.18 87 7.33 40915118 48.73 87 1.31 67401885 19.24 87 1.62 307112477, 26:00 87	911053 8.91 87914 911050 12:18 40968 911079 10:49 88324	586) 15.46 4092 989) 30.47 4092	0548  45.52   09 0555  20.75   67 0563   50.49   409	173998 10.48 3674 114607 22.90 3672 221470 50 50 3672	123046 55.69 0 123087 33.21 6	7418525 16.67 9173550 8.50 7418566 23.82		
4	87912853 22 87912861 31 87912879 34	90 307/11/2485 27.14 879 39 8879	971087 1771 409180 971095 18.92 091784	9976 49.41 4092 26 16.73 4092	0571 25.70 089 0589 60.21 409 0597 69.61 000	91:135 12:73 3674 21:488 69:47 409	23160 59.73 6 23898 63.66 09	418608 22.60 173568 13.98		
5	67394361 37 87912887 36	.90 307412485 2714 875 .89 40915126 71,36 8876 .12 69 40915134 74,87 674	03600 34 28 400 400	130 60.03 3676	6 19 1409	21496 77.60 EATO	To le land	14 24 1		
1/2 5/8	67394445 36. 67394486 31	07 307112493 10.05 674	Belling Jorge Sant	80.0 L	31.21 PAU92	1512 85:75 (1)	8 0 40	23518 85,75	orto 🖺	
1 11/2	67394569 54.5 67394601 86.6	0/112501 12.34 6740	3766 8.26 0 2410 3808 10.15	2017年1日の日本の日本の日本の日本の日本の日本の日本の日本の日本の日本の日本の日本の日本	505   15.53   4092 513   18.50   4092 521   17.29   4092	1520 25.10 4092	3914 26.65	49 5 32 52 87	100+ 3+4	
11/2	67394643 78.4 67394684 92.1	1 40915175 41.68 6740	3881 14,49	080473 P   51 00 009212   51 00 40:272 21 004	4092	1546 28,59 4092 1553 36.44 4092 156 38,52 8775 1579 35,31	CAL	04   13   100   15	2 01:	
3/8	UNC	6740		16 1981 300 s	\$2002 2003 1000 1000	287 41/76 de leu	9102 9102 9102	10 40 20 10 (20) 140 41 1 28 8	3	
7/16 1/2 5/8 3/4	67394726 20.89 87912903 11.26 87912914 15.80	265765623 14.70 67403		409206	7 29.15 40921	595 27,73 631904	79105 3 40 87 5 79105 0 27 52		POL PAR PAR	
7/8			046 10.76 40919052 087 11,25 40919060 3831 9.24 307113202 176 10.62	28.43 4092066 19.04 6741320	2 43.65 674146 3 7.83 674146	349 - 21.85 - 40923 80 - 16.81 - 674232	955 42.62 6741 963 48.14 6741 902 31.76 6741	3640 14.32 3681 13.91 1723 16.84 SEV		
11/4 11/2 13/4 2	0702000 22.01	A(1)12543 14.38 87911 40915191 4665 82611	337 9.22 97713203 1/6 10.62 7713203 86 12.23 97914594 94 14:29 40914976 95 15:47,387914692 16:18:15 40919086 28:19:34 487914617 28:19:34 487914617 38:19:34 487914617 38:19:34 487914617	20:72 6741324 35:82 40920670 30:39 40920688	15.63 674147 24.58 409216 40.35 052453 45.14 409216 55.82 674747 58.83 409216	22 24.38 674232 11 31.74 991732	71 53,23 4092 44 40.45 -6741 12 16.98 67418	526 38.87 3004 764 18.54 1004	56,1	
21/4	87912900 25.39	07112550 46.77 879112 40915209 56.86 879112 07112568 50.83 879112	28 18:94 887914610 86 20:00 524:15 59	46.48 40920696 55.71 40920704 40920712	45.14 409216 55.82 674147	22 24.38 67.8232 21 31.74 99.1732 60 12.99 40.9239 29 40.59 40.9239 30 47.73 40.9220 31 44.34 88	89 75.80 67448 97 50.64 67418 03 94.35 67418	22.62		
234 3 3½ 3½ 3¾	87913000 26.41 87913018 32.41 87913026 35.90	07112576 55.80 674041	29 28.83	Va.10 40920720	61.25 4092164	56.19 LA CVE 2 46.87 NATON	007918 0 367418	62 37 14	4.6	
4 3	37913034 43.83 37913042 46.41	0915217 72.89 38791127	31.48 40919110	76 1979 37 1 1923 3 86.99		00.54	6 111.95 674190 409235	02 43.72 34 63.71	CHE COLD A . TO	
5 5½ 6	37913042 46.41 17913059 50.42 7394882 70.05 7913067 69.32 4	98791128 96740420 96740428	8 45,11 862,904 10 8 45,000 863	0 209213	i socialis	281001	08 1516480 08 1516480	C3:US	3	
3%-16 UN	C 7913083 18.10	_ Trible   4.1	The state of the state of	C0211909	0.004 E	10   1881.004   g	101018	061201 061201	.12	
74 38 78 38	0915514 37.80 44 7913109 18.44 40 7913117 28.92 40 913125 19.17 507	915283 58:32 67404288 915241 68:28 67404368 915268 75:21 67404409 172584 34:50 87911368 172502 41:58 87911368	15.44 18,67 20.19 40919136 63	40020746	50.34 40921686 40921694 59.77 40921702	51.03	E4092854	2 36.44	1 Ft 13	
11/2 387	913126 19.17 507 913133 23.14 09 913141 28.03 407	112584 34.60 87911350 173394 14.24 87911368 112592 41.68 87911376	20.19 40919136 63 25.44 60919144 60 19.33 40919151 70 21.64 40919160 70	02 40920761 21 40920779	0.01 d9174020	51,21 C40924087	86.82 6741904	24.70		
2 387 21/4 387		015274 101.91 87911384 015274 101.91 87911392	25.88 28.48 40919177 88	40920787, 8 27, 40920795, 8	09174088 40921710 8.27 67414947 40921728	82,42 67,60 a 104.7a 88	d40923567	39,83 43,57 78,97	A A A	
Length is o	cran jength 'Head i	s undercut <sup>3</sup> 50/Box <sup>4</sup> 10/Bo	x <sup>5</sup> 25/Box	57 ETC126 (3)	40921728		40923583	113:56	16. 2 5. 5. 5. 5. 5. 5. 5. 5. 5. 5. 5. 5. 5. 5	
1996		CALL	MOO TO				KORON A 10/Box	to observe over	Continu	
		OALL	MSC TODAY 800-	645-7270 FO	RALL YOUR	INDUSTRIAL N	EEDS			


7	s - Zinc	Plat	ed Stee							· lou	তি প্রত্য	Histo	5-01	.un 6	(D) IN	in.
W)	Plat Head P	hillips Price/ 100	Round Head	Slotted Price/ 100	Round Head	Phillips Price/	Pan Head S	lotted Price/	Pán Head P	hillips Price/ 100	Fillister Hea	d Slotted Price/ 100	Truss Head :	201 40.00	Binding H	
<b>1</b> .01	40915290	\$120.57	<sup>3</sup> 87911418	\$28.68	40919193		6	i i i i i i i i i i i i i i i i i i i	40921736	\$120.57	0100 P	100	40923591	\$136.37	0700	Ť
68.92 46.92 86.72	407112600 40915308 40915316	61:94 141:87 148:96	387911426 387911434 387911442 387911459	59.37 38.56 41.09 43.14	40919201	135.38	340920803	\$131.42	67414888 -40921744 -40921751	137.31 118.65 148.96	1,85   675 5,70   675		40923609	94.78	02756	iNi
0,1 0,1 0,8 7,8 7,8	4091 <b>53</b> 24 4091 <b>5</b> 332	162.80 166.35	387911467 387911475 387911483 387911491	52.38 53.17 58.39 53.73	\$2.3 \$6.0 \$0.4	.01:	- 6.22 63 5.38 63 5.79 63	16776 16776 16776	4.96 6 6.22 6	YOUN	4.47 879 4.63 879	hadire.1	4.90 675 6.22 675	1	6756	
301.99	100 8	8865V6	09173592 87911509	88.16 65.99	9.19 6.30	000 Section	ie ii e	.3885a	33.1 38.1	67 (36) (17 (36) (4) (33)	6 12 6 9 5 76 879		40923617	118.25	C-63.10.13.13.13	200
943,88 647,40 988.61 172,08	140915357 40915365	146.70 180.84	387911517 387911525 387911533	72.51 63.97 86.95	76.1				3.84 11.85 18.65	Tables Lymany Lymany	218 218 218		10020017			
412.88 412.83 4194.57	.40915373 .40915381	192.76 218.08	387911541 367404483 387911558	69.55 138.93 100.24	25.22 0.00 0.00	71901 71901 71901	5.93 63 68.8	489655 500688	7,07 6 6,73 67	22,05 [6-64]	5.04 879 5.33 879		6,52 675 6.73 675	8307 5MG	6756	Pla
92.13 2.108.52 1/188.51 1/160.72	409 5399 21 8 409 5407	182.05 252.85	87911574 87911582 87911590	114.67 147.95 145.18	8.28 7.00 6	083/74 (03/14 (03/14	2.05 5. 8.32 6. 8.84 6.	369987 56: 125	6 83.5 58 88.8 9 88.8	1091 S	5.99 879 6.69 870 7.56 876	047,7V 5 1 0965 1 6363	7,56 675 9,05 675 9,68 675	[887]   1082]	6756 6756 8756	333
158.35 162.81	40915415	331.53	°87911616 ⁵87911632	218.24 166.17	a   61.0°	771.0	59 (04.0h	26, 954	19.64 67 14.72 15.25	242 24230 24230			11,39 675	1003	6756	
Divisor	ad is undercu S + Mili	CARL OF	1	0237-63	24693	/352	<b>Q6</b> <sup>80.8</sup>	569202	8.06 67	24289	888 L.H	758827	675	360		*
Phi Wilary Equiv	ps Flat Head 100° Order#	Price/10	00 Military Eq		an Head Order#Pri		L Pkg. (in.) Qty.	Militar	Phillips Plat He y Equiv: (	ad 100° Order #	Price/100 4	Military Equiv		Price/10		
24693-S	The state of		No Ch	76	8 88.6 8 80.6	7 546	#8-32 UNC 100 5/6 100		93-547 720	Fee 19 057326	\$8.65 M	35206-2 35206-2	41. 7205786 42. 7205787	0 8.6	7 2	
24693-S 24693-S	6 7205703 17 7205704	7 5.3 5 5.4	MS35206	-202 72	057524	5.37 5.54 5.64	% 100 -76 100 1/2 100 5% 100	MS246 MS246	93-S48 720 93-S49 720 93-S50 720	057384 057342 057367 057875	- 9.26 M	635206-2 635206-2 635206-2	43 7205788 44 7205789 45 7205790	8 9.0 6 9.2 4 9.6	6 2 Phil	lips of 10
HHIVO:			MS35206 MS35206 MS35206 MS35206	-204 72 -205 72	057540 057557	5.64 5.79 5,95	34: 100 -78: 100 -1: 100	MS246 MS246	93-852 720 93-853 720	057383 057391	11.09 M 11.73 M 12.48 M	35206-2 35206-2	47 7205792 48 7205793	0 10.9 8 11.7	9	iu ii
Total Laves	12,21 13,27 15,67 15,64	0 10	MS35206 MS35206 MS35206	-207 72	057573	5,95 6,23 6,37	13/16 100 11/4 100 11/2 100	MS246	93-S56 72	057409 057417 057425 057433	13.10 13.85 M 15.10 M	35206-2	50 7205795	3 13.8	5	II.
<b>y</b> 152 34 3	77.17	2	MS35206	-210 72 -211 72	057607 0 30.11 957615	6.67 5.95	1¾ 100 #10-24 UNC 5/16 100		10.59	4478 4486 4494	0.8 P. J.	335206-2 335206-2	52 7205797 60 7205798	7 10.4	9	
M\$24693-5 M\$24693-5 M\$24693-5	2 7205705 3 7205725	2 6.3 0 6.4	5 MS35206	-213 72 -214 72	057631 057649	6.09 6.17 6.45	7% 100 7/16 100 -1/2 100	enseet theeset	2 .33	24592 24530 245281	813   81 M 978   879 M	35206-2 35206-2 35206-2	62 7205800 63 7205801		8 Phillips 9	Pa
MS24693-5 MS24693-5 MS24693-5	5 7205735	8 6.5 9 6.8 1 7.0	9, MS35206 7, MS35206 MS35206 MS35206	-216 72 -217 72	057664 057672	6.45 6.98 7.10 7.28	% 100 34 100 76 100 1 100	MS246	93-S74 72	057458	e 13.85 M	35206-2 35206-2 35206-2 35206-2	65 7205803 66 7205804	5 13.8 3 14.8	5 2 • Ma	teria
MS24693-8	9 7205747	7.9	MS35206 MS35206 MS35206	-219 72 -220 72 -221 72	057698 057796 057714	7.74 8.01 8.29	1½ 100 1¼ 100 1½ 100	MS246	13 (33.41)	057466 6664 8664	16.72 M	35206-2 35206-2	68 72 <b>6580</b> 6 69 7205807	8 17.6 6 21.6	g cad yell ster	lmiu low el
MS24693-S	(100	9 9.6	MS35206 8 MS35206	-223 72		1083	134 100 #10-32 UNF	(	93-5270 72	1004 6104	4.8 B. y	335206-2	20.51	43	<b>→</b>	
MS24693-S	13 7205706	6 71		U.100	78 18.5	7.12	7/6 100 -1/2 12 100 -5/6 100	MS246 MS246 MS246	93-S271 72 93-S272 72 93-S273 72	057186 057144 057151	11.48 S 11.52 11.21	S PALSE TOCAME	27.05 32.3¢	=   enga 20 - ana	100	
MS24693-S MS24693-S MS24693-S	25   7205709 26   7205710 27   7205711	4 7.4 2 7.6 0 7.8	MS35206 7 MS35206 2 MS35206	-227 72 -228 72	057755 057763	7.40 6.48 6.59	78 100 78 100 1 100	MS246 MS246 MS246	93-S274 72 93-S275 72 93-S276 72	057169 057177 057185	13.85 14.85 15.82	a. 3s	Acian reba		101	vic'
MS24693-S MS24693-S	29 7205724	3 8.6	0 MS35206 5 MS35206 MS35206	-231 72 -232 72	057797 057805	6.79 7.20 7.62	11/4 100 11/4 100 13/6 100	MS246 MS246 MS246	93-S277 720 93-S278 720 93-S279 720 93-S280 720	57498 57291 57215	16.72 17.69 18.68					
MS24693-S MS24693-S	32 7205727	6 - 10.1	6 MS35206 MS35206	-233 <i>12</i> -234 <i>72</i>	057821	7.95 8.37 9.09	1½ 100 1⁄4-20 UNC 1⁄16 100 1½ 11100	MS246	30.0	0039	29.55 M	35206-2	78 7205809 79 7205810 80 7205814	2 17.6		JMI.
MS24693-S MS24693-S MS24693-S	36 7205729 38 7205730	4 11.7 2 12.0 0 13.1	7 MS35206 8 MS35206			9.85 0.68	34 100 76 100	MS246	93-S95 72	57482 3574	20.17 M	35206-2 35206-2 35206-2	81 7205812		9	
	34,37 39.98	2/4/c	(3.187.52	787	6 (1) 8	47872 38977	76 100 1 100 -14 100 1½ 100	3.700	- DESCRIPTION DE	4759	M Size	35206-2 35206-2 35206-2	83 7205814	2 24.7 9 27.9	7	ا د
in Edwards			1		154		134 100		10	62 1			86 7205817		9	


<sup>1</sup> Length (In.) #2-56 UNC	Flat Head Order #	Slotted - Price/100	Flat Head Order #	Phillips Price/100	Round Hea Order #	d Slotted Price/100	Pan Head Order #	Slotted Price/100,	Pan Head Order #	Phillips Price/100	Oval Head Order # 1	Slotted Price/10	Oval Has Order A	
1/8 3/16 1/4 5/16 3/8	67563809 67568841	\$4.76 5.70	67566406	3.65	67568287 67568329	\$5.09 6.13	67568642 67568683 67568725	5.04	67570564 67570606	\$5.00 4.57 4.17 4.69 5.15	1 000 - 2 144441	L.	72.00	
#4-40 UNC  ***  ***  ***  ***  ***  ***  ***	67563882	4.00	67566448		67568360	7.16	67568766 67568808	6.22	67570762 67570804 67570846	5.70 6.75 4.82	(2.5) (3.6) (3.6) (3.5)	6.0	80 MA	
5/16 3/8 1/2 5/8	67563924 67563965	6.22 6.41	a call at land	4.47 4.63 5.23 6.12	87924007 87924023 87924049 87924056	6.22 6.87	67568840 67568881 67568923	5.30 5.79 6.41	67570929 67570960 67571000	3.64 4.03 4.13 5.19				
3/4 1 11/4 11/2	3.87 3.87 3.97 3.97 3.97		67566646	5.70	87924064 87924080 87924106 87924114	7.25 7.85 8.84 11.86 16.05			67571042 67571083	5.15 5.30	91.55 92.68 1.76.77 17.00	2	30 55 30 55	
#6-32 UNC 1/4 5/16 3/6 1/2	67564005 67564047 67564088	6.52 6.73 7.56	67566729 67566760	5.04 5.33 5.99	87924155 87924163 87924171	7.01 6.73 7.56	67568964 67569004 67569046	5.93 6.39 7.65	67571125 67571166 67571208 67571240	5.20 5.54 5.68 6.46	67574764	. 14 159 55 140; 50 . 3 (0) 7.32		
3/4 1: 11/4 11/2 2	67564120 67564161 67564203	9.03 9.68 11.39	67566802 67566844 67566885	7.56 8.32	87924189 87924197 87924213 87924221- 87924239	9.68 10.64 14.72 15.25	67569087 67569129 67569160	8.32 8.84 10.40	67571281 67571323 67571364	7.60 9.15	67574806 67574848		676.00 6767.44 6767.4	
#8-32 UNC	67564245 67564286 67564328	9.65 9.87 11.04	67566927 -67566968 67567008 67567040	6.52 7.32	87924254 87924288 87924296 87924304	22.10 8.06 9.65 9.87	67569202 67569244 67569285	8.06 8.98 9,88	67571406 67571448 67571489 67571521	9.09 8.59 8.29 8.89	S <b>Noile</b> S Noiles Hackages	2001 3 (4) (4)	67873 67873	
3/4 7/8 1 11/4 11/6	67564369 67564401 67564443 67564484	10.95 15.13 17.49 18.93	67567081 67567123 67567184 67567206	10.23	87924312 87924320 87924338 87924346 87924353 87924361	12.31	67569327 67569368 67569400 67569442 67569483	11.04 11.50 14.31 17.43 18.80	67571562 67571604 67571646 67571687 67571729	9.71 10.76 13.65 14.03	67574889 67574921 67574962 67575001	12.47 15.54 17.94 19.39	6767 67573	
21/2 21/2 3 #10-24 UNC	Reide	1/20/302 1/20/302 1/20/302	\$206-246   7 206-246   4 205-247   2 206-249	KOV 2	87924379 87924387 87924395 87924403	24.48 26.90 33.63 41.67	40626699	45.27	1.54 1.64 1.64	; eve	202 206 206 4206 204 7.05	NULAY, N PM JE PM JE	3	
% 1/2 5/6 3/4 7/8	67564526 67564567 67564609 67564641	12.00 14.31 15.54	67567248 67567289 67567321 67567362	13.13 11.43 11.43	87924437 87924445 87924452 87924460 87924478	13.94	67569525 67569566 67569608 67569640	10.20	67571760, 67571802 67571844 67571885	11.11 15.36 15.83	67575043 67575084 67575126 67575167	12.97 15.71 17.31 17.60	675734 675736 675736 675736	
11/4 11/4 11/4 13/4 2 21/2	67564682 67564724 67564765 67564807	23.83 27.38 33.10	67567404 67567446 67567487 67567529	20.57 23.28 33.16	87924486 87924494 87924502 87924510 87924528	23.28 27.33 29.04	67569681 67569723 67569764 67569806	22.03 30.76	67572008	27.40 26.54	67575209 67575241 67575282 67575324	21.03 25,12 27.72 36.93	6757364 6767361 6767372 6767372	
3/4	67564849 67564880 67564922 67564963	13.59 15.82	67567602 675676449	11.27 13.13 10.75	87924544 87924569 87924577 87924585 87924593	12.29 13.02 15.82	67569848 67569889 67569921 67569962	13.13 15.88	67572081 67572123 67572164	11.86 14.62	67575365 67575407 67575449	12.95 15.71 17.31	6757380 6757380 6757380 6757392 6757386 67574004	
1 1¼ 1½ 1½ 1¾	67565002 67565044 67565085	20.51 24.17 27.05	67567727 67567768 67567800	14.80 19.15 22.29	37924601 37924619 37924627 37924635 37924643	19.59 19.14 24.17 27:05 30.20	67570002 67570044 67570085	19.59 23.10 26.74	67572248 67572289 67572321	17.61 v 19.97	67575522	20 52	675740; 675740; 675741;	
2 2½ 3 #12-24 UNC	67565127		67567842 40626715	28.30 31.62	7924650 7924668 7924676	32.34 ( 45.75 4 61.50	<b>10626707</b> 28152M 58853W	30.84 41.55	1.40	5 5	67575647 050 050 050 055	35.12	67574100 67574200 6757424	
1/4 1 11/4 11/2	7.65		52.7	81.6 81.6	7568402 7568444 7568485 7568527 7568568		V8824592	(4) / (4) (4) (4) (4) (4) (4) (4) (4) (4) (4)	7.20	10.00	230 720 -231 729 -232 720 -233 720	1500 1 12800 1 17830 16031	01.0.2 88.8 66.974 A	
9/8 3/4 1	67565168 67565200 67565242 67565243 67565325 67565366	26.02	7567925 7567966 7568006	23,86 8 23,12 8 30,81 8 35,62 8	7568600 7924700 7924718 7924726 7924742 7924759	26.14 6 26.02 6 29.37 6 18.32 6 32.20 6 39.93 6	7570168 7570200 7570242 7570283 7570325 7570366	23.01 23.67 26.41 25.28 31.86 37.46	90 ( 57572404 57572446 57572487 57572529 57572560	20.11 28,43 20,13 6 29.96 85.62	97575688 97575750 97575761 97575803	26.74	6757428 4 67574327 67574366 67574400 67574442	
ength is meas		r flat head					<sup>2</sup> 50/Pkg.	37,46 6	i   l''''''	13.11 6	37575845	44.95	67574483 30	

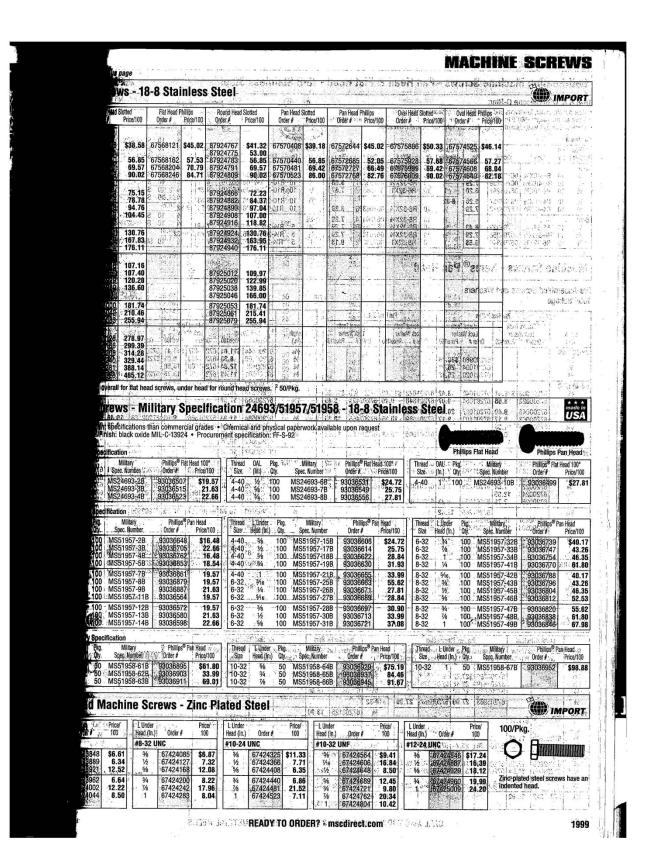

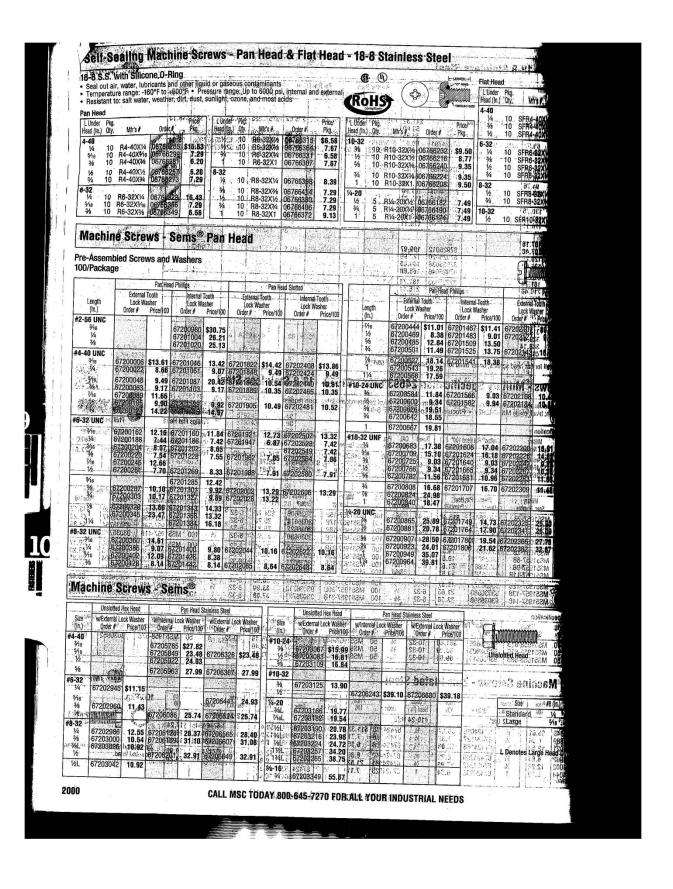

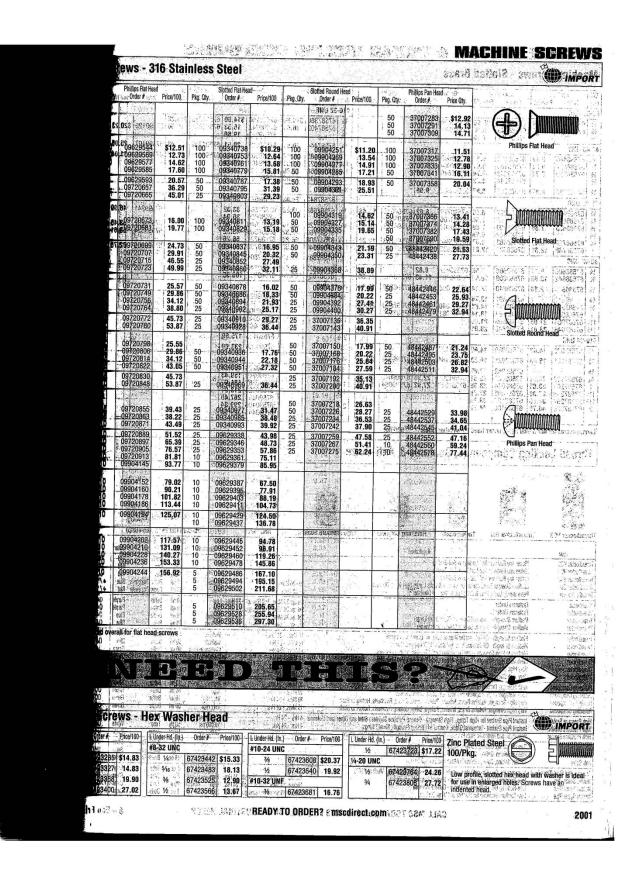

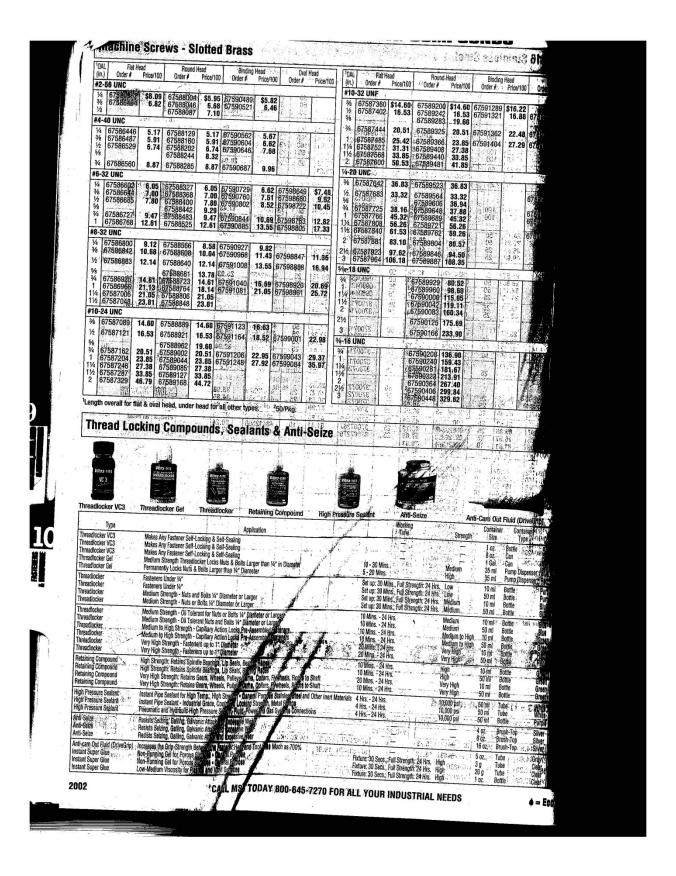


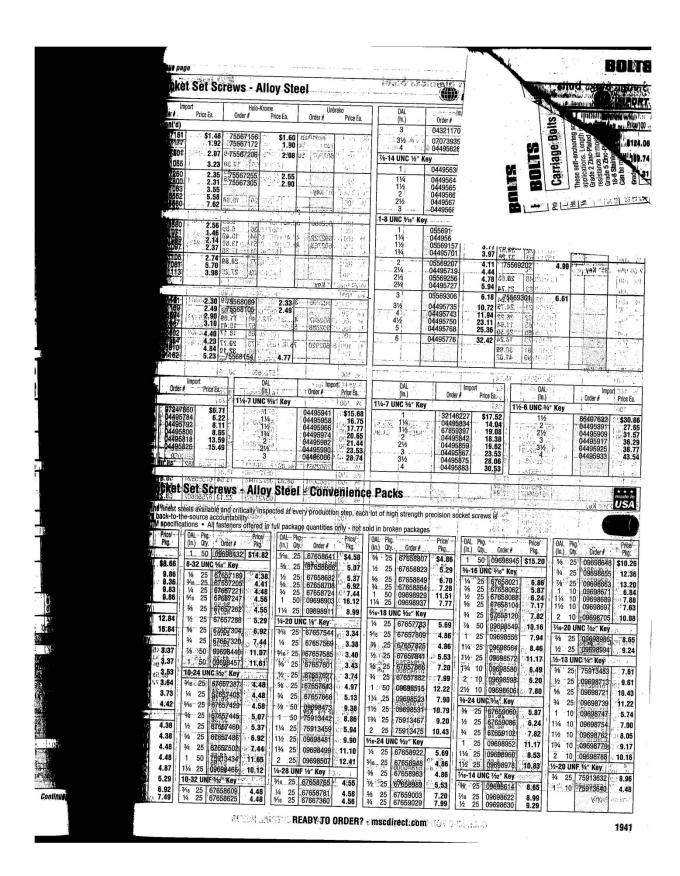



Made II  Order #  185326981  25 4 4 2 5	Price/ 100	Holo-Kr Order#		Unbra		-				MPORT	USA	HOLO-KING	A-10.		OO
18 <b>532698</b> 18532698			Price/		ako Price/ 100	Under Head (In.)	Pkg. Oty.	் Impo	Price/	Made in Order #	Price/	Holo-Kr Order#	Price/	Order#\	Unda Hearic
W		1 (208412) 1 (0081n/2	255 251	67804344 67804351	652.07	1½ 15/8 13/4	100 100 100	05512157 76588540 05512173	21.91	260674181	\$26.06 28.66		38.34		1 1
-15614BC	8 7 8 8 8 8	6 00741751 1 0074160 3 0084181		67804385 65863 668656 666666	704.52 G.	21/4 21/2 23/4	100 50	05512207 05512223 05512256 00103291	124.00	<sup>2</sup> 60674231 60674264	29.57 40.39 43.20 65.89		39.56 54.01 57.78 70.78	67805887 67805929 67805960	6
ebaraes	1067 - 10 1,62 14 8. 63 1	28548G07(1 C871800   8	, 2°	C282100 (	0.00	31/2	50 50 50	05512306 00103309 00103317			55.43 79.07		13:74.14 0:117.10	467806000 5648860	+1
A-1715-CONT. AND THE P.	12.24 13.87 14.50	75511063 75511071	16.37	67804401 67804443 3 67804484	32.64	41/2 5 51/2	50 50 25	.00103325 76588557 00103333 04162335 076588565	112.98 153.01	7 7552460 3 7552460 8 7653330	#34 (3 PSS (3	75612400 74633041 75512509 74633058 74633066	163.43 189.18 190.83	US52460 0552165 1,05424 0010344	
60673134 60673167	16.43 18.26	75511121 75511154	21.98	67804567	34.80 38.99	1/28	100	598678139		85327005	28.63		Ad op	5) 80/90 F	
60673209	22.38	75511204 75511220	27.10 29.92 34.36 38.65	67804724	57.46	7/8	100 100	05523071 05523089 05523105 05523121	32.37 33.56 35.08 39.40	00841411 00841429 00841312 200841320	35.18 37.05	75523084 75523100	40.61 47.42 48.53 51.40	67806083 67806125	68. 73. 74. 81.
00841247 00841254 200841262	52.72 71.62 93.81	74633017 74633025		67804773 67804781 3838 267804799	297.32 240.89	1¾ 2 2¼	100 50 50	05523154 05523170 05523204 03773637	42.77 45.60 47.53 53.48	200841338 200841346 00841353 00841361	43:23 47:47 52:05 60:29	<sup>12</sup> 75523159 <sup>2</sup> 75523175 75523209 75523217	59.32 - 65.69 67.26 71.78	-267806208 267806240 167806281 67806323	,89. 101. 107. 117.
) 30 22 9	i	B	269.97	)   0416283 		2¾ 3 3¼	50 50 50	05523279 05523303 03773645	59.66 68.67 71.90 86.60	400841387 400841395	65.27 75.26 79.38	75523258 475523274 475523308	83.19 92.65 103.56	467806406 467806448	125.1 147.2 156.8
oggetter)			800 S	CORSOLI (	J 6,1	4	50	05523402	108.43 138.08	- 1,-348	91.73	475523357 475523456 114866	122.45 136.50	eraread in	180.3 201.0
60668076 60668084	14.20 14.54	75522052 75522060 75522078 75522086	17.16 18.28 19.45	67804849 67804880	31.32 32.03	5½ 5½ 6	25 25 25	00103358 04162376 76588573	157.35 307.70 461.55	7 7561417 51 7564420	1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1 1	75523506 1800 8	214.00	7650971 041,0691417 319,0861421	3 3
60668100 00841270	19.10 36.89	75522102	25.53	67804963 67804989	35.99 385.59	- 5/8	100		77.20	tU ylqquO	of the	2001CE	N O S	xeefso?	in 3.
00841288 60668167	18.51 47.44 20.58 22.16	75522151	27.53	67805002 67805028 67805044 267805085	39.24 166.64 42.83	11/4	100 100 100	05513122 05513155	84.15 39.46 41.60 44.99		47.50	75513150	51.76 59.05 66.25	67806604 267806646 267806687	83.2 89.8 99.9
260668209 60668233 60668266 60668282	24.36 29.12 33.65 44.89	75522201 75522227 75522250 75522276	32.58 38.93 44,99 60.05	267805127 267805168 67805200 67805242	53.62 66.94 73.67 98.86	2 21/4 21/2 3	50 50 50 50 50	05513205 76588581 00103366 00103374	52.82 83.93 91.91 97.09	00841478 00841486 00841494 400841510	57.04 96.92 110.65 120.89	75513200 74633074 74633082	76.94 238.32 264.77	67806729 67806737 467806745	118,6 1256.2 1129.6
60668332 60668365 7 18 132	65.63 01 79.98	75522326 275522359 £134800	87.78 87.78 689.59	67805283 67805325 67805366 67805382	107.74 130.28 148.05 288.67	5	50 . 25	76588599 76588607	401.15	ACTBORAS.	125.72	74633090 08+5000 1	294.21	1 7658695	589.5
1551622 1551625	8 <b>7.37</b> 86.1 86.1	75522409 15800 75522458	104.73 121.57	67805408 67805440	173.47 207.67	3/8 1/2	100 50	76588623 03575842	86.57 33.27	60669066	37.99	75524058	48.90	67806760	154.29
<del>1089 radi.</del> Beefsat	166.83	75522508 75522532 275522540	156.23 164.25 214.73	67805507 67805523	250,02 1345,12 353,20	534 78 11%	50 50 50	05524079 05524087 05524108 76588631	36.99 37.89 38.52 48.10	60669082 60669090 60669108 200841536	39.45 40.42 43.20	75524074 75524082 75524108	50.12 51.40 57.78	67806844 67806885 67806927	75.8 84.1 83.6
134 - 111. 134 - 111. 134 - 12	\$4.5	9064222	S	67805531 67805549	729.91 743.69	13/8	50 50 50	00103390 05524152	38.83 45.36 42.81 112.44	00669132 00841544 60669165	-C (I	CACAGGG	57.43 60.41	UBBBCD\ I	99.33
60674066 60674074	19.10	75512053	8.3	0.16	3 644.51	2 21/4	50 8 50	05524178 05524202 05524228	67.70	60669181 60669207	65.05	76524207	65.72 80.31 89.87	67807040 67807081 67807123	110,25 118,65 133,55
60674082 60674090 60674108	19.84 21.39 21.82	75512087 75512103	27.52	67805689 67805721	41.27 42.00	2½ 2¾ -3	50 51	05524277 05524301	75.01 79.96 83.73	60669306	82.61 86.71 93.89	475524256 475524272 475524306	116.25	67807248	142.16 164.00 175.16
60674132	23.67	75512129 0-3±300		67805762	45.23	31/2	25	05524350	93.84	60669363			C 1 1 1	356010th	189.31 201.31
	60673092 60673092 60673107 60673167 60673167 60673163 60673233 60673233 60673233 60681267 60681267 60688092 60668092 60668107 60668107 60668108 60668108 60668209 60668108 60668209 60668108 60668209 60668108 60668209 60668332 60668209 60668332 60668209 60668407 60668409 60668407 60668409 60674082 60674082 60674082 60674082 60674082 60674082	60673092 14.50 60673100 15.50 60673167 18.25 60673167 18.25 60673269 22.38 60673269 22.38 60673269 22.38 60673269 23.50 60673266 23.50 60681254 71.62 6068103 14.54 60686167 25.50 6068183 22.16 60686167 25.50 60686167	60673092 14.80 75511089 60673100 15.80 755111089 60673101 15.80 755111089 606732167 18.26 755111154 60673209 22.38 75511205 60673269 22.38 75511205 60673269 22.38 75511205 7551206 75511206 751	60673092 14.50 75511081 19.35 60673100 15.88 75511106 29.96 60673101 16.88 755111106 29.96 60673101 18.26 75511150 29.96 60673107 18.26 75511150 29.96 60673260 22.38 75511150 29.96 60673260 29.38 75511200 29.34 36.0673220 29.38 75511200 29.34 36.0673220 29.38 75511200 29.34 36.0673220 29.38 75511200 29.34 36.0673260 29.39 75511203 29.30 75511203 29.	60673100 15.51 75511083 19.35 167804326 60673104 14.40 75511151 29.48 67804305 60673194 14.40 75511151 21.81 167804305 60673187 14.25 75511154 24.41 167804305 60673269 22.47 75511174 21.16 167804305 60673269 22.48 75511204 29.47 167804305 60673269 23.48 7551204 34.36 67804705 60673269 23.48 7551204 34.36 67804705 60673269 23.48 7551204 34.36 67804705 60673269 23.48 7551204 34.36 67804705 60673269 23.48 7551203 34.85 67804768 37804768 34.85 67804768 34.85 67804768 34.85 67804705 60641254 34.85 67804705 60641254 34.85 67804705 60641254 34.85 67804705 60668076 14.20 75522060 18.22 67804805 60668092 15.55 75522060 18.22 67804805 60668009 19.10 75522102 25.53 67804905 60668100 19.10 75522102 25.53 67804905 60668103 21.55 75522105 25.55 67805040 60668133 21.6 7552206 32.55 67805040 60668133 21.6 7552206 32.55 67805040 60668133 21.6 7552206 32.55 67805040 60668092 21.55 75522107 25.55 67805040 60668092 21.55 75522107 25.55 67805040 60668209 21.55 75522210 32.55 67805040 60668209 21.55 75522201 32.55 67805040 60668200 21.55 75522201 32.55 67805040 60668200 21.55 75522201 32.55 67805040 60668200 21.55 75522201 32.55 67805040 60668200 21.55 75522201 32.55 67805040 60668200 21.55 75522201 32.55 7505040 60668200 21.55 75522201 32.55 7505040 60668200 21.55 75	60673102 14.59 755511083 19.35 767804464 32.84 60673102 15.58 75511105 29.86 67804564 32.24 60673187 14.29 755511154 24.49 67864609 38.98 60673167 12.29 75511174 27.10 67804609 38.98 60673269 22.38 75511204 29.92 67804609 38.99 60673229 22.38 75511204 29.92 67804609 38.99 60673229 22.38 75511204 29.92 67804662 47.51 60673266 28.99 75551204 39.92 67804662 47.51 60673266 28.99 75551204 39.92 67804662 47.51 60673266 28.99 75551205 38.85 67804765 48.90 00841224 77.82 74633037 38.95 67804769 488.81 74633033 289.97 67804789 488.81 74633033 289.97 67804789 488.81 74633033 289.97 67804849 31.22 60668076 14.29 75522060 18.28 67804849 31.22 60668094 15.55 75522060 18.28 67804849 31.22 60668094 15.55 75522060 18.28 67804849 31.22 60668094 18.55 75522060 18.28 67804849 31.22 60668097 18.55 75522060 18.28 67804849 31.22 60668092 23.30 57805122 28.80 60680892 24.35 75522060 18.28 6780489 38.59 60668076 18.29 75522060 18.28 6780489 31.22 60668092 24.35 75522060 18.28 6780489 38.59 60668092 24.35 7552206 25.55 6780508 28.50 6066809 24.35 7552206 25.55 6780508 28.50 6066809 24.35 7552206 25.55 6780508 28.50 6066809 24.35 7552206 25.55 6780508 28.50 6066809 25.55 6780508 28.50 6066809 25.55 6780508 28.50 6066809 25.55 6780508 28.50 6066809 25.55 6780508 28.50 6066809 25.55 6780508 28.50 6066809 25.55 6780508 28.50 6066809 25.55 6780508 28.50 6066809 25.55 6780508 28.50 6066809 25.55 6780508 28.50 6066808 28.50	60673100 15.88 75511161 21.99 167694367 34.80 160673107 18.26 75511161 21.99 167694367 34.80 160673167 18.26 75511161 21.99 167694367 34.80 160673167 18.26 75511161 21.99 167694367 34.80 160673167 18.26 75511161 21.99 167694367 34.80 160673167 18.26 75511161 21.99 167694367 34.80 160673167 18.26 75511161 21.99 167694367 34.80 160673269 22.37 755111201 34.80 167694699 38.99 36.60673269 22.38 75511201 34.80 167694692 38.99 36.60673269 22.38 75511201 34.80 167694765 84.90 114.00841284 71.82 74633017 83.78 67894765 84.90 114.00841284 71.82 74633017 83.78 67894765 84.90 114.00841284 71.82 74633033 28.97 867894765 84.90 114.00841284 71.82 74633033 28.97 867894765 84.90 114.00841284 71.82 74633033 28.97 867894823 581.83 21.26 60688064 14.34 75522060 18.28 67894860 31.22 21.00841284 71.00841284	60673100 15.87 75511067 28.98 67804464 22.84 6 25. 60673100 15.87 755111067 28.98 67804566 33.23 23. 60673134 16.48 755111154 21.89 67804560 34.89 34. 100 60873187 18.28 755111204 28.92 67804567 34.89 34. 100 60873187 22.77 755111740 27.19 67804567 34.89 34. 100 60873209 22.38 75511204 28.92 756167682 47.51 76. 100 60673266 28.89 75511204 34.38 67804662 47.51 76. 100 60673266 28.89 75511204 34.38 67804765 44.80 11. 100 60673266 28.89 75511203 34.38 67804765 44.80 11. 100 60673266 28.89 75511253 38.85 67804765 44.80 11. 100 60641254 71.82 74633031 83.97 67804823 581.83 25. 5606808076 14.29 75522060 18.28 67804845 31.22 50. 34.65	60673100	60673133	60673103	606/6907190 14.59 75511095 19.35 107804464 22.44 6 25 276698565 20.38 10.50 19	60673190 14.50 75511051 2.0 248 6.7604626 32.2   60673134 14.4   60673167 14.7	60073730 1.4.0 75511063 19.35 [7690494 22.44] 67.00 60073107 1.0.0	605/73/107   1.3.5   755/11/62   1.3.5   756/14/64   2.4.5   6.2.5   756/14/64   2.4.5   755/11/62   2.4.5   756/14/64   2.4


#### FASTENERS & HARDWARE **FASTENERS** FASTENERS **FASTENERS** EXTRACTORS: NOT TO THE DESIGNATION OF THE PARTY OF THE PA Machine Screws, continued: WASHERS, continued: Military Spec,, Stainless Steel .....2035-2036 1.1.2122-2123 Military Spec. 24693/35206 ......1997, 1999. 11...12113-2122 Self-Sealing Pan Head & Flat Head 18-8 SS RE ...2113-2122 Nylon Acme.......2019 Cage Clip-On , "J", "U", Rectangular ... 2033-2034 Sems. Ogee, Pipe Lock & Pyramidal .....2046 3 ... Stainless Steel 18-8 ... Shims .......2040-2042 Stainless Steel 316, Slotted Brass, 1982-1983 Hex Washer Head ......2001-2002 Thread Cutting ......2008-2009 Hex Machine Nuts, Hex Nuts Mil Spec. 2022-2023 ...1981-1982 Sheet Metal Screws ...2010-2012 1987 1984 Hex Speed Nuts - Hot Dipped Galvanized ... 2033 Wing, Thumb Screws......2007-2008 Tooth Lock, Internal/External ......2045-2047 Keps & Conical ......2029, 2031 Wood Screws ... 0 vin ...... 1965 Locknuts (Anti-Vibrational) ...2026, 2030-2032 Vibration-Resistant, Self-Sealing ..... 2042-2043 SOCKET PRODUCTS: Wave ......2043 1953-1960 Panel Nut Retainers, Jack, Pal, Well ... 2034-2035 Basics of Socket Head Fasteners ...... HARDWARE Vibration-Resistant, Self-Sealing ......2029 Allen Nuts......2025 H .....1961 Wing, Acorn, Square, Handle .....2027-2029 Access Doors Button Head Cap Screws 1967-1968 1962-1964 11974-1978 etich and long it aits u.u. a since of short BOLTS: OLTS: Barrel/Chain/Slide/Spring Braces/Brackets ...1978-1980 Arrow Clips Cheese Head Metric Machine Screws 2015 Cheese Head Metric Machine Screws 2053 Stainless Steel, Torx Plus® & Flange 1,1930-1932 Flat Head Cap Screws: Bumpers..... ....1971-1972 Circuit Board Support Fasteners -----Alloy Steel ..... .2172-2174 1968-1970 1971 1971 & Hex-Locknuts - 2015 & Hex-Locknuts - 2017 Finished Hex Nuts, Flat Washers, Acom Nuts, Wing Nuts - 2015 Door Closers ..2138-2140 Door Guards, Plates & Hardware Door Reinforcers/Protectors 9/2143 t loc Stainless Steel ......1925-1926 1973 hniqi Torx Plus®..... Door Stops/Holders/Viewers 2143-2145 Tork MusPressure Plugs 1950-1952 Shouldei Screws/Shripper Bolts: 1934-1935 Alloy Steel with Nylon Patch 1933 1836 Keyhole & Push-In Panel Rivets ......2088 1972 1184-1988-1990 Nail Drive Anchors 2007 2007 Push Mount Blind Rivets 2007 Door Weatherstripping ......2176-2177 1/41/11/1983-1985 Ratchet Fasteners 2015 Retaining Lock Washers 2015 ELECTRONIC LOCKING DEVICES 1 9 2514 (A.III.D.):::1987 Door Holder/Release Devices: .....2136-2137 & Shoulder Washers ......2015, 2017 Precision 1937-1938 ION DOI 2093 Socket Head Cap Screws: Slotted Knurled Head Thumb Screws...2014-2015 xtro Electronic Exit Devices 2137 Alloy Steel Electronic/Programmable Locks 2137 1902-1909 Handicap Access Actuators ......2137 Alloy Steel with Nylon Patch ......1910 .....2096-2099 Corrosion Resistant 2100-2102 Low Head ..... PINS & CLIPS: Basics of Dowel Pins .... 1932-1933 .....2102-2104 .....2095, 2099 Metric 1917-1920 Metric Stainless Steel 1920-1921 Garage Door Replacement Parts......2161 .. 2066 Gas Springs 2158-2159 Gate Hardware 2162-2163 Cotter Pins, Hairpin, Safety, Snap ... 2074-2075 Mil. Spec. 16995/16996 ......1916 .....2011-2012 Dowel Pins, Precision Ground .....2066-2069 Hitch, Linch, Gib Head, Lock Pins .......2076 Mil. Spec. 16997 & NAS A-286 Alloy ... 1914 Hasps ......2148-2149 Hurricane Clips 2113 Pull Dowel Pins 2069-2070 Pull Dowel Pins W/ Spiral Air Relief 2070 2097-2098 1936 Spec......1909-1910 ..2149-2152 Reference Charts..... 2098 Hinges Specialty 2152-2155 Hooks 2178-2179 Key Blanks & Tags 2145-2146 Stainless Steel 18-8 & 316 ....... 1911-1914 Torx Plus® & Zinc Plated ......1915-1916 Roll Pins/Spring Pins ......2072-2073 Slotted Spring Pins - Stainless Steel .....2074 Socket Set Screws: Split Rings & Clevis Pins .....2076-2078 .....2110-2112 Metric ......1949-1950 Loads ...2105 Military Spec. 51021/51023 ..... Taper Pins 1..... .....2071-2072 ...1943 LOCKS: 501 501 Cable/Chain Locks Nylon Tipped ..... 1948-1949 Camicorks 2146, 2146 Oylinders 2131-2132 Deadbolts 2131-2132 Knob Locks 2130-2131 Square Head ..... TAMPER-RESISTANT FASTENERS: Ring Tools, External & Internal ..... 2064 2065 IVETS: | 001 Basics of Blind Rivets | 002 Scrows 1.. 1965-1966 RIVETS: ...2046-2049 Cap Screws ....1915 Cap Screws .....1932 Latches/Latch Locks......2146-2147 p Screws ...1919 Threaded Inserts..... .....2091 2049 THREADED ROD: Open & Closed End ......2081 Acme.... .2019 Padlocks/Combination Locks ......2124-2128 N2049, 2051-2052 Anchors ..... .....2021 Rim Locks & Mortise Locks ..... Studs.... 2022 302134 Threaded Rod ..... ...2018-2019 PULLS: THREAD LOCKING: Compounds, Sealants & Anti-Seize..... Quarter Turn Fasteners ......2146-2147. .....2059-2061 VELCRO®: .....2057 ..2112-2113 1 & Metric .....2058 WASHERS: er Kits ...2054-2058 SCREWS: Corrosion-Resistant Machine Screws: .2035-2036 Manhers. ...2052-2053 Dished Tooth ... Alloy Steel, Stainless Steel 18-8...1993-1999 Fiberglass, Brass & Silicon Bronze. .... 2037-2038 Tension & Toggle Latches ......2170-2172 Metric Alloy/Stainless Steel ......2003-2007 Tie Downs Flat .2037-2039 Track Door Hardware .... 2036-2037 2163 1901


## 6.8 - Appendix H- Custom Parts

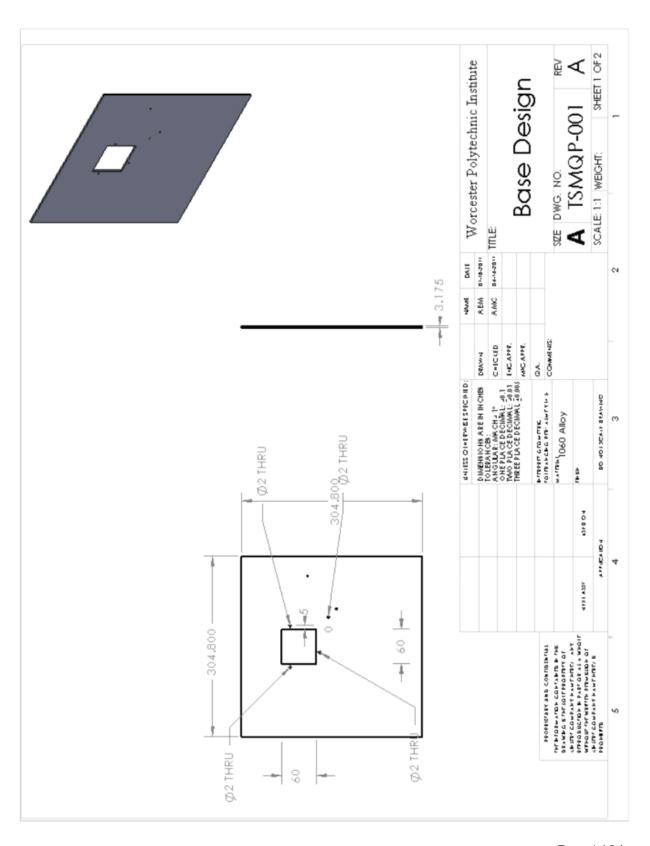


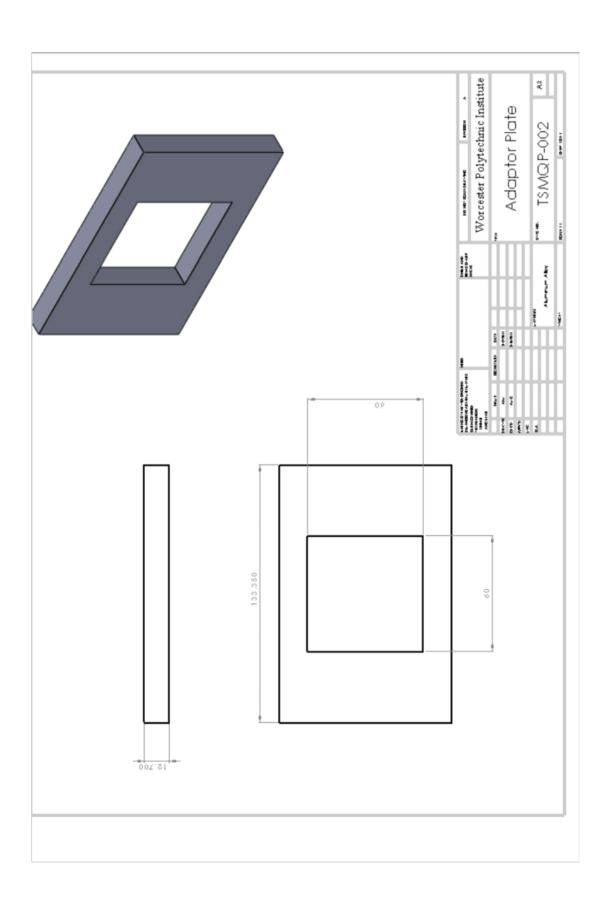


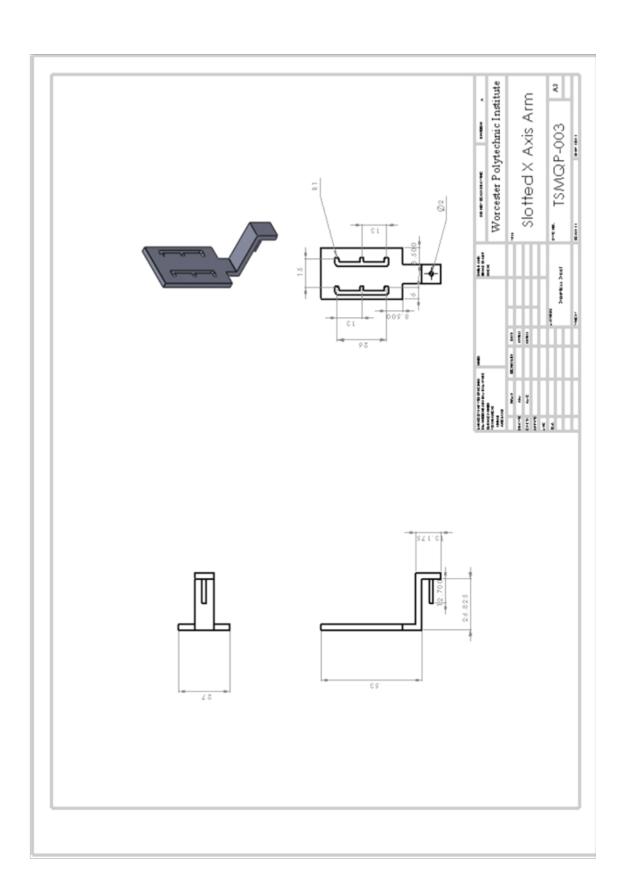


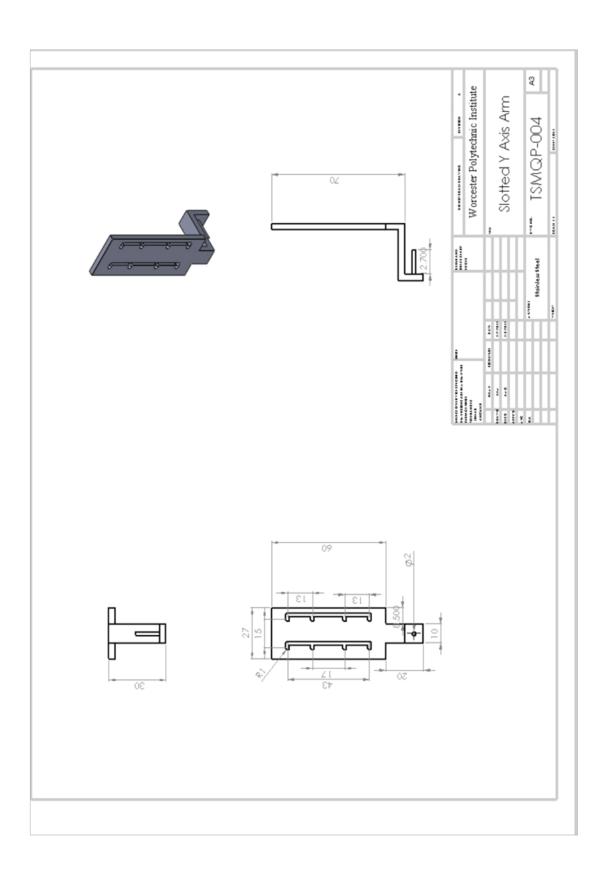


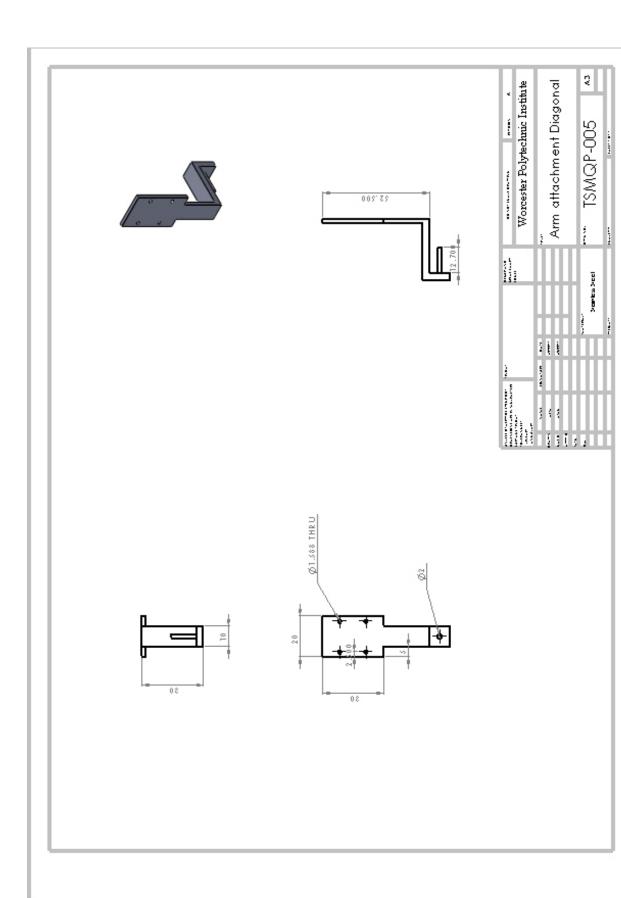


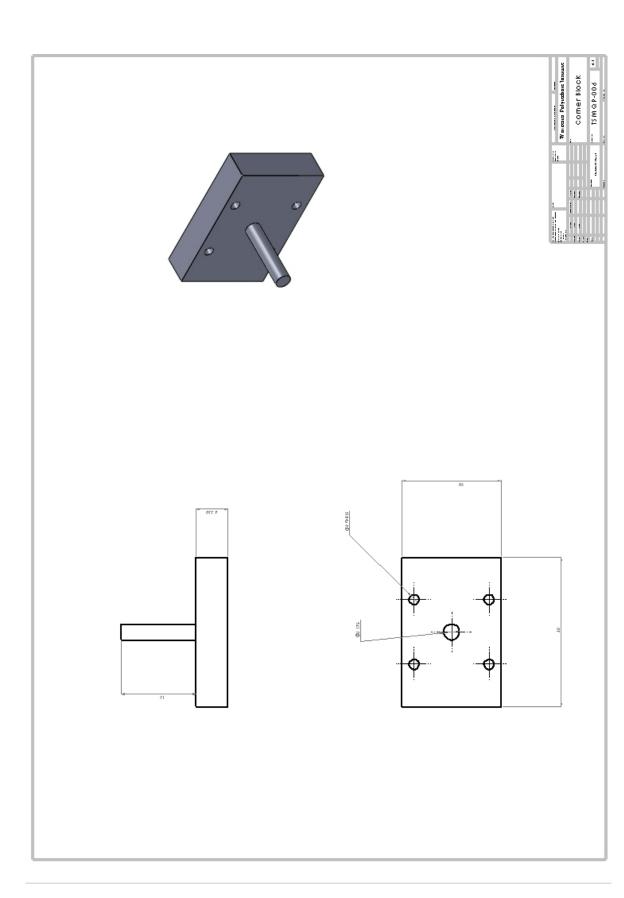


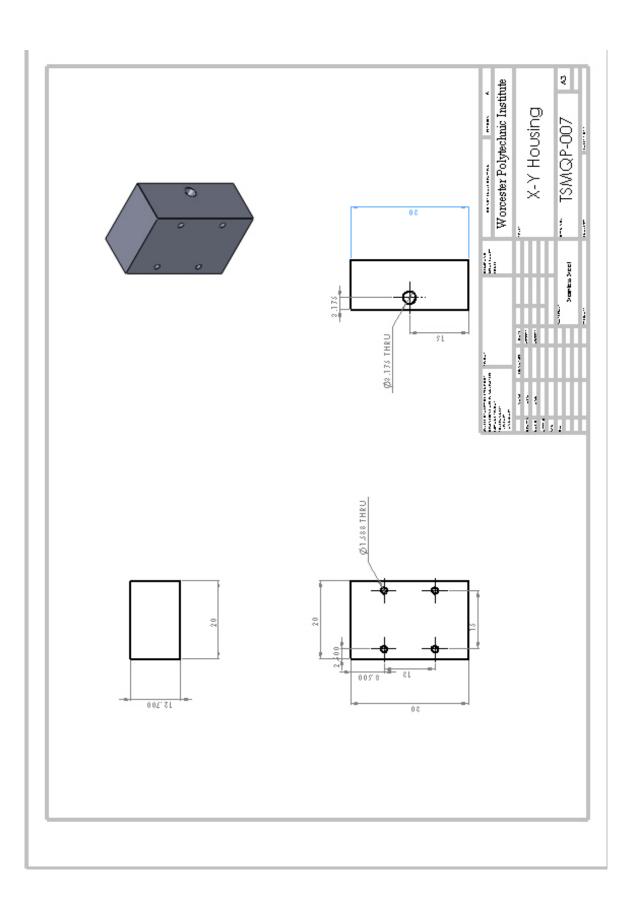


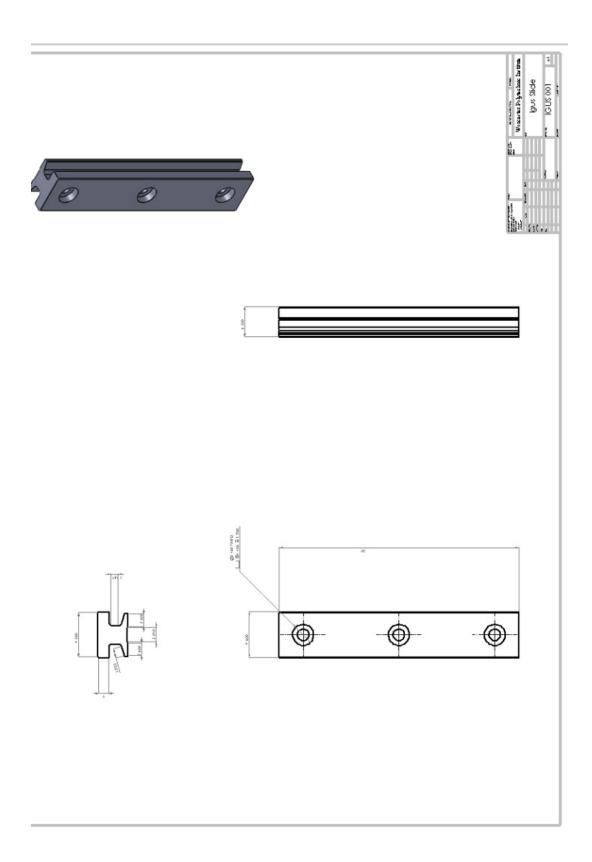


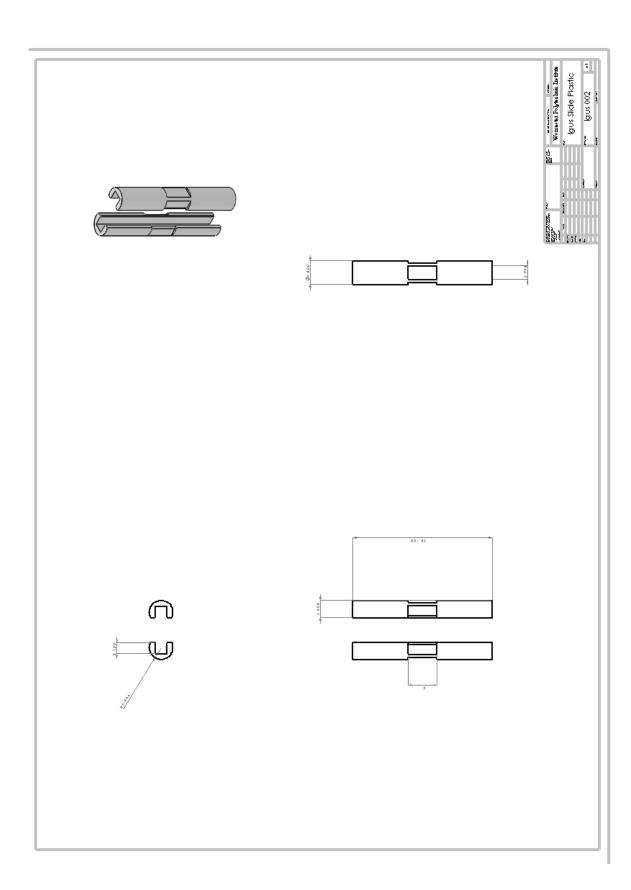


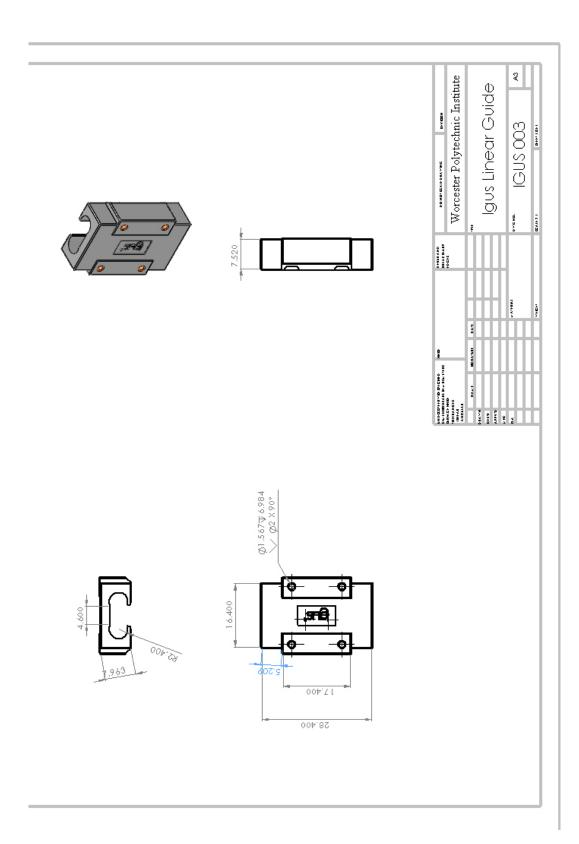


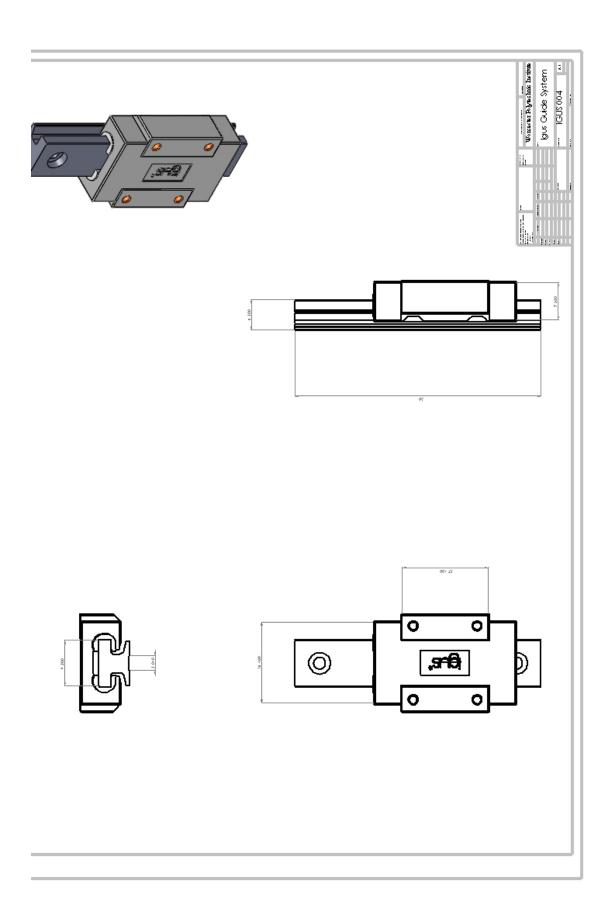


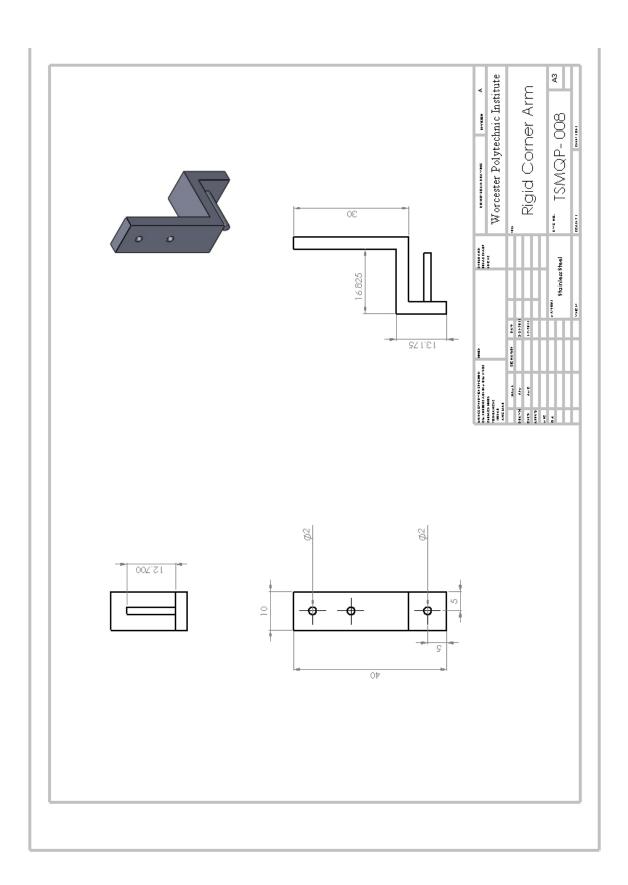


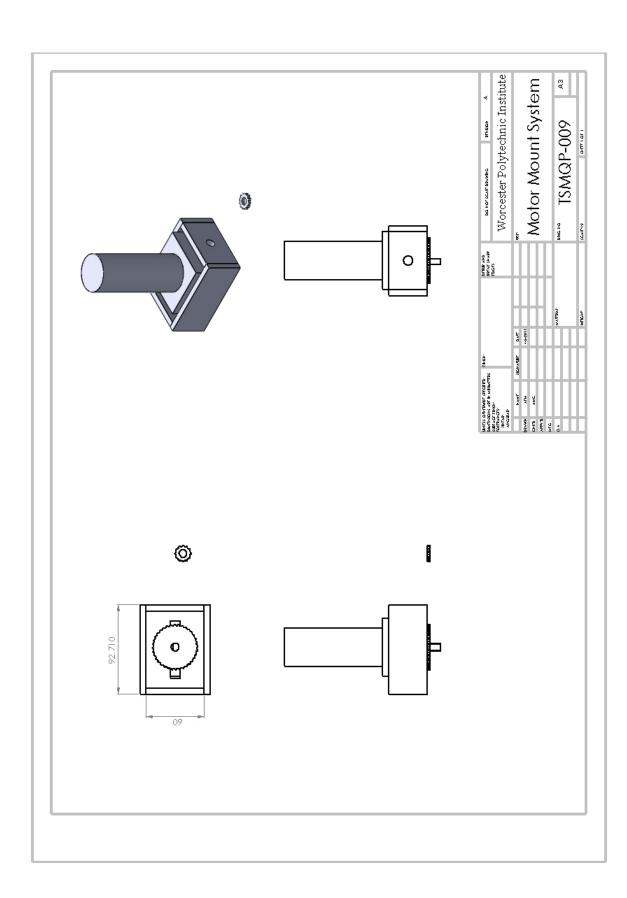


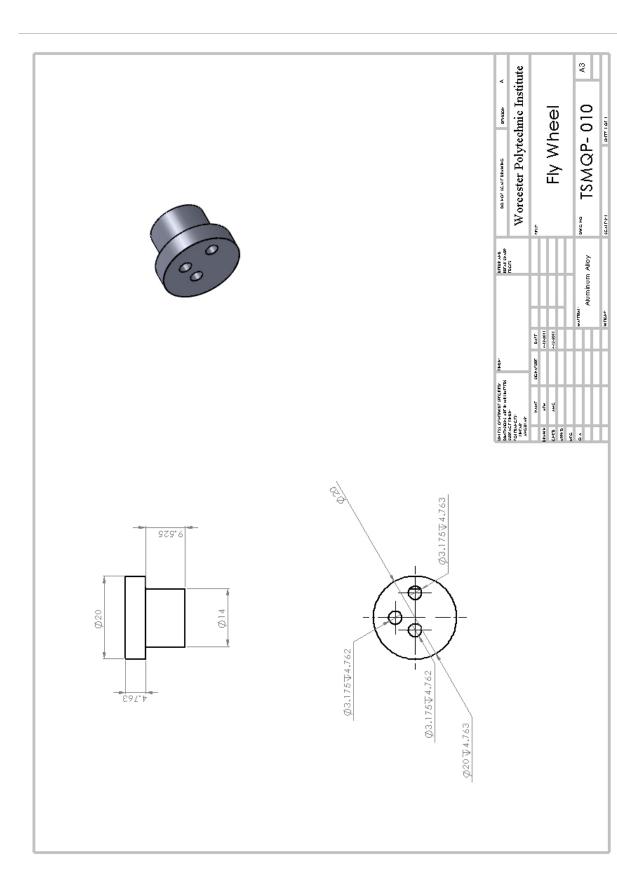


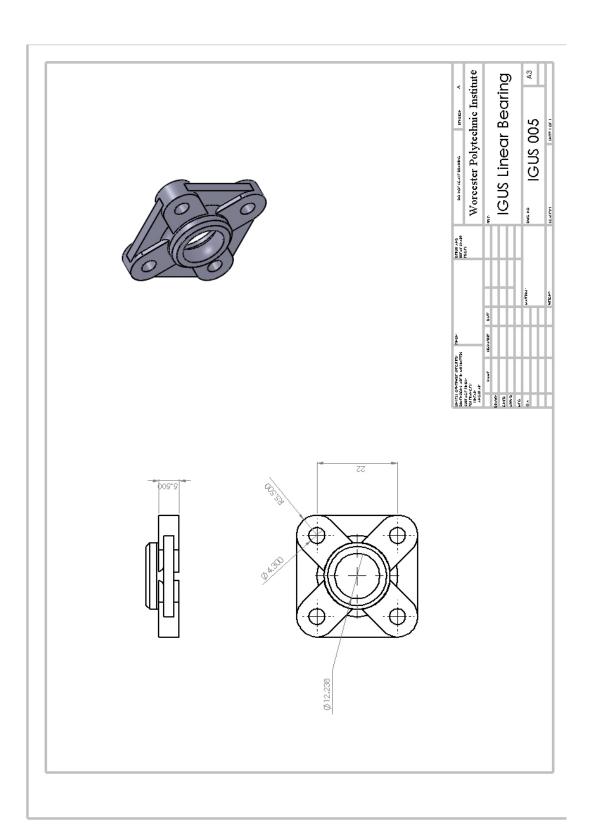

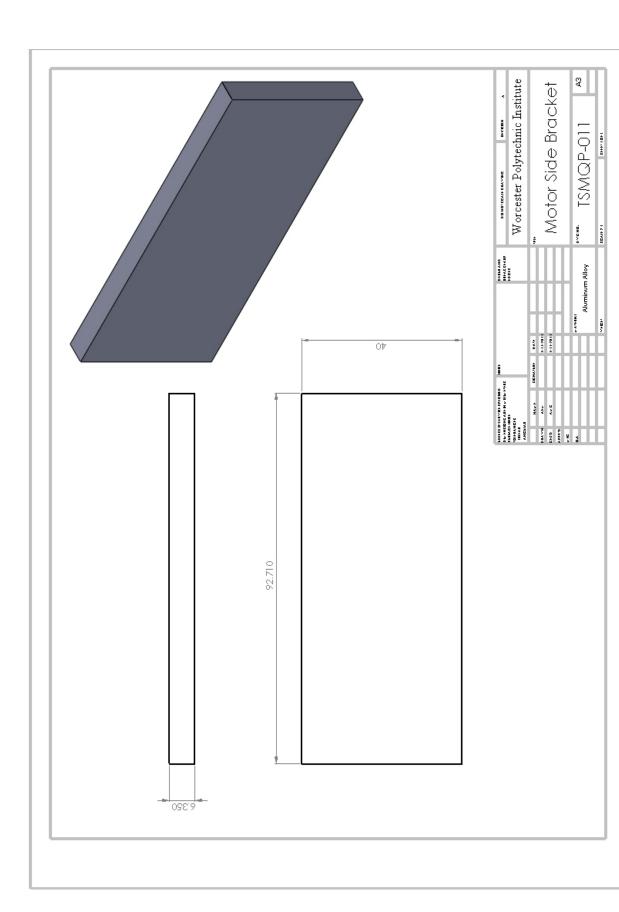

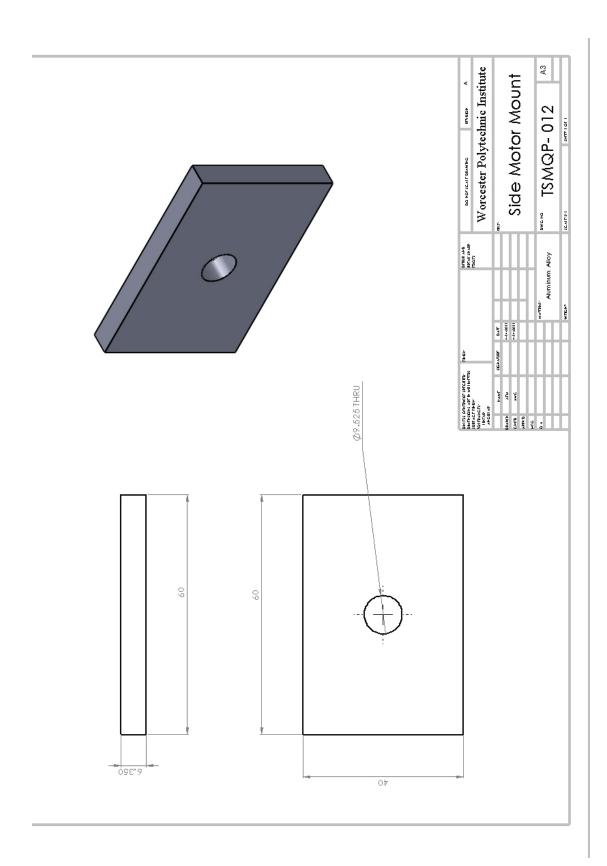

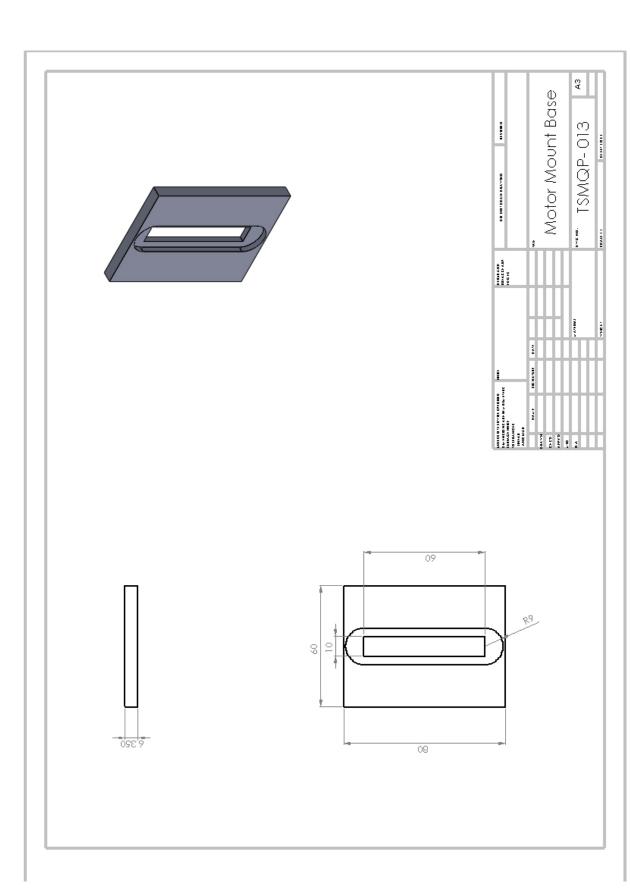

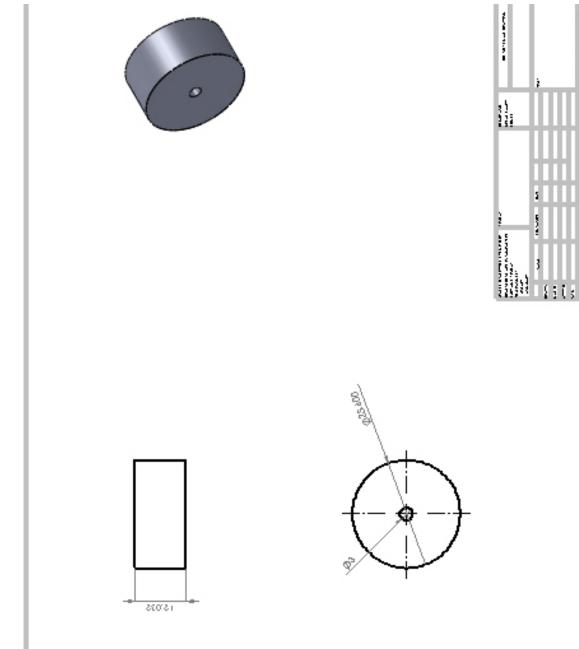

## 6.9 - Appendix I- HDM Matlab Compile Function

This Matlab code was used to compile the over 100 images taken by the camera and put them into matrices.

```
function compile_data(dataset,framerate,res,loadext)
%COMPILE_DATA Compiles and stores data from the HDM Program
%    Imports and reads the HDM data from the input video folder. This data
%    is then stored in a .mat file as a large multidim matrix along with
```

```
several descriptive variables to identify the data.
응
   Also, saves data in both a smoothed (using datasmooth.m) and a raw form
9
%% Prepare global variables
   disp(['Compiling data for: ',dataset])
   tic:
    top_dir=['data_store\',dataset,'\'];
%% Verify input
   if framerate < 1</pre>
        error('framerate must be a real positive number')
    end
    if ~exist('loadext','var')
       error('loadext input must be specified')
    end
%% Create the output directories
   disp('>>Creating output directories');
    if ~exist([top_dir,'HDM_Data\'],'dir')
       mkdir([top_dir,'HDM_Data\'])
   end
    if ~exist([top dir,'matlab data\'],'dir')
       mkdir([top dir,'matlab data\'])
    end
%% Check for the existence of HDM data and move it to the correct folder
    disp('>>Checking HDM input');
    % Check for the data files in the correct location
    if ~exist([top_dir,'HDM_Data\start00+im-0001_U.dat'],'file')
        % Check for files in the images folder
       if ~exist([top_dir,'images\start00+im-0001_U.dat'],'file')
           % Return an error
           error('The HDM Data does not exist for this dataset');
       else
           % Move the files to the correct location
           movefile([top_dir,'images\*.dat'],[top_dir,'HDM_Data\']);
           movefile([top_dir,'images\*.txt'],[top_dir,'HDM_Data\']);
        end
    end
%% Load the correlation report file into variables
   disp('>>Loading correlation report')
    *Open summary file and extract xy corner data throughout the data
    %collection process
        file=fopen(strcat(top_dir,'HDM_Data\CorrelationReport.txt'));
        textscan(file,'%s%s',1);
       str='%*s %*s %u %*s %u %d %d %d %d %*s %d %*s %d';
       rawdata=textscan(file,str);
       xxyy=double([rawdata{3} rawdata{5} rawdata{4} rawdata{6}]);
        subsize=double(rawdata{1}(1,1)); %#ok<NASGU>
        shift=double(rawdata{2}(1,1));
       pixelshift=double([rawdata{7} rawdata{8}]);%#ok<NASGU>
       fclose(file);
```

```
%% Prepare to loop through the dataset
    disp('>>Preparing to loop through dataset')
    % Overapproximate the width of the data matrix
        cols = round(1.1*(xxyy(1,2)-xxyy(1,1))*shift);
    % Create a format string for data processing
        format='%9f';
        for i = 2:cols
            format = [format,' %9f']; %#ok<AGROW>
    % Load the first data file and determine the exact # of columns
        filename = [top_dir,'HDM_Data\start00+im-0001_U.dat'];
        file = fopen(filename);
        testdata=textscan(file,format,'multipledelimsasone',...
            1,'collectoutput',1,'emptyvalue',0);
        testdata=testdata{1};
        for i = 1:cols
            if testdata(1,i)==0
                break;
            end
        end
        cols = i-1;
        fclose(file);
    % Create the final format string for data processing
        format='%9f';
        for i = 2:cols
            format = [format,' %9f']; %#ok<AGROW>
    % Determine the number of rows and columns in the data file
        file = fopen(filename);
        testdata=textscan(file,format,'multipledelimsasone',...
            1, 'collectoutput', 1, 'emptyvalue', 0);
        testdata=testdata{1};
        [rows.cols]=size(testdata);
    % Determine the number of data files to be processed
        filelist = dir([top_dir,'HDM_Data\*.dat']);
        filecount = numel(filelist)/2;
        fclose(file);
    clearvars filename file testdata filelist
%% Prepare data filter
   hsize=[res,res];
    sigma = 0.5;
    filt = fspecial('gaussian', hsize, sigma);
%% Loop through each data file
    disp('>>Beginning loop');
    % Preallocate memory to output variables
        udata_r = zeros(rows,cols,filecount);
        udata_s = zeros(rows,cols,filecount);
        vdata_r = zeros(rows,cols,filecount);
        vdata s = zeros(rows,cols,filecount);
        errorcount = 0;
        bar= waitbar(0,'Compiling...');
    % Begin time loop
```

clearvars file str rawdata

```
tstart=now();
    for t = 0:filecount-1
        waitbar((t+1)/filecount,bar);
        %disp(['>>>>Compiling data at time ',num2str(t+1),' of
',num2str(filecount)]);
    %% Determine file names
        if t==0
            uf='start00+im-0001_U.dat';
            vf='start00+im-0001 V.dat';
        elseif t==(filecount-1)
            uf=['im-',numpad(t,4),'+end0000_U.dat'];
            vf=['im-',numpad(t,4),'+end0000_V.dat'];
            uf=['im-',numpad(t,4),'+im-',numpad(t+1,4),'_U.dat'];
            vf=['im-',numpad(t,4),'+im-',numpad(t+1,4),'_V.dat'];
        end
       uf = [top dir, 'HDM data\', uf]; %#ok<AGROW>
       vf = [top_dir,'HDM_data\',vf]; %#ok<AGROW>
    %% Process X-Direction data
       curfile = fopen(uf);
        try
            rawdata = textscan(curfile, format, 'multipledelimsasone',...
                1,'collectoutput',1,'emptyvalue',0);
        catch err
            err,
        end
       rawdata = rawdata{1};
        smoothdata = filter2(filt,rawdata);
        %errorcount = errorcount + error1;
       udata_r(:,:,t+1)=rawdata;
       udata_s(:,:,t+1)=smoothdata;
       fclose(curfile);
    %% Process Y-Direction data
        curfile = fopen(vf);
       rawdata = textscan(curfile, format, 'multipledelimsasone',...
            1, 'collectoutput',1, 'emptyvalue',0);
       rawdata = rawdata{1};
        smoothdata = filter2(filt,rawdata);
       vdata_r(:,:,t+1)=rawdata;
       vdata_s(:,:,t+1)=smoothdata;
        fclose(curfile);
    %% Calculate remaining time and update waitbar
            tleft = ((1-((t+1)/filecount))/...
            ((t+1)/filecount))*(now()-tstart);
       waitbar((t+1)/filecount,bar,...
            ['Estimated time remaining:',datestr(tleft,'HH:MM:SS')]);
   end
   delete(bar);
    %errpercent = errorcount/(rows*cols*filecount); %#ok<NASGU>
    % Clear extra variables
        clearvars uf vf curfile rawdata smoothdata error1 errorcount t
%% Mark the area of interest in the first image and save it
```

```
disp('>> Creating overlay image');
    % Load image
        srcimg = imread([top_dir,'images\start00.tif']);
    % Create RGB image container
        outimg(:,:,1) = double(srcimg);
        outimg(:,:,2) = double(srcimg);
        outimg(:,:,3) = double(srcimg);
    % Create a polygon over the region of interest
        x = [xxyy(1,1), xxyy(1,2), xxyy(1,2), xxyy(1,1)];
        y = [xxyy(1,3), xxyy(1,3), xxyy(1,4), xxyy(1,4)];
        [h,w]=size(srcimg);
        poly = poly2mask(x,y,h,w);
    % Reduce the blue and green channels in the ROI
        poly=double(poly);
        poly(poly == 1) = 0.6;
        poly(poly == 0) = 1;
        outimg(:,:,2) = round(outimg(:,:,2).*poly);
        outimg(:,:,3) = round(outimg(:,:,3).*poly);
        outimg=uint8(outimg);imshow(outimg);
    % Save the image to the matlab data directory
        imwrite(outimg,[top_dir,'matlab_data\overlay2.tif'],'tiff');
%% Check for external data and process
     if loadext > 0
         disp('Getting external data files');
     % Check if the external data mat file exists, if not run the helper
        if ~exist([top_dir,'matlab_data\ext_data.mat'],'file')
            comp ext helper(dataset);
        end
        load([top_dir,'matlab_data\ext_data.mat']);
     end
%% Export data variables to compiledhdm.mat
    disp('>>Saving data file')
    clearvars bar cols filecount format h i loadext outimg poly rows...
        srcimg w x y;
    engine='HDM1'; %#ok<NASGU>
    save(strcat(top_dir,'matlab_data\compiledhdm.mat'));
    tfinal = toc;
    disp(['Compile data complete in: ',...
        datestr(datenum(0,0,0,0,0,tfinal),'HH:MM:SS')])
    clearvars
end
function num = numpad(num,len)
    if isnumeric(num)
        num = num2str(num);
    end
    if length(num) < len</pre>
        num = numpad(['0',num],len);
    end
end
```

# 6.9 - Appendix J - HDM Displacement in the X Direction Raw Data

-82.17724779	-82.22123719	-80.3652447	-85.63889416	-90.08098457	-93.369115	-92.1982484
-88.81157095	-88.62061156	-91.39818425	-94.75976318	-98.43451148	-97.36129978	-92.96267975
-87.71959477	-91.42168378	-93.8656273	-92.27668136	-92.78360105	-96.43758753	-96.44625323
-89.31652009	-91.92339704	-92.59659147	-90.31967256	-94.22626862	-93.74303669	-94.313364
-83.74240006	-85.30926914	-85.71791167	-84.03563259	-88.09009927	-91.62606406	-91.16331727
-81.33184147	-82.74908484	-83.45654255	-84.48432805	-82.74449713	-87.56414012	-88.62337021
-79.88400966	-81.2443718	-81.89661604	-86.62240251	-82.95951799	-82.71300817	-84.99202304
-74.02867805	-74.22834748	-78.7272802	-80.31395278	-79.85331368	-80.59690028	-80.97284454
-72.67840114	-72.3334041	-75.79480397	-77.00680931	-77.35965602	-76.58552079	-74.82980804
-71.11325202	-72.52858594	-73.09256614	-73.88804831	-73.89948762	-72.52282432	-75.24959294
-68.54392838	-68.16623884	-70.53393689	-71.23414386	-71.83166936	-71.73829751	-73.10867459
-66.6847268	-66.80315196	-67.64958785	-68.14477328	-69.70118966	-68.45624167	-70.32376114
-62.72187583	-57.10637479	-63.74009081	-65.18942301	-66.28937571	-66.40464052	-68.75642405
-61.05394922	-61.18806797	-62.13506801	-62.99941936	-61.50464371	-63.83583997	-65.32904956
-61.32188283	-59.47855361	-59.80407435	-59.99190703	-59.96319703	-60.92500121	-61.08213652
-58.16443664	-55.24656239	-58.07699814	-54.04523766	-57.3133421	-58.82735603	-59.09908877
-56.75615891	-56.51487462	-62.31146432	-55.68738048	-56.03706081	-55.86464981	-56.14373967
-53.77848042	-55.61543221	-55.75698674	-53.99954076	-53.59490901	-53.84599031	-53.79270828
-55.53293586	-51.26584481	-51.29850828	-52.28600798	-51.37358884	-51.5139752	-51.38793544
-50.21837824	-48.12019551	-49.79926575	-48.92821039	-48.7471195	-49.75498827	-48.54934614
-45.35513797	-45.60975204	-46.72116423	-46,4366046	-43.54528598	-47.06341718	-47.30137482
-44.99813563	-45.58643279	-45.94834625	-45.33723247	-44.56110726	-44.2226626	-45,45259699
-43.32699429	-42.45725848	-48.24416669	-44.25547916	-43.13982047	-43.03867596	-42.76841466
-41.84771832	-41.63946569	-42.38645543	-40.37638752	-38.9885705	-40.87314075	-40.97064589
-38.40044486	-42.62001631	-45.30009613	-39.532965	-38.587952	-38.32918085	-38.40398924
-36.0012261	-29.17935281	-36.65125224	-36.69500909	-36.39073822	-35.08057831	-36.76690844
-35.3905884	-34.58906455	-35.51722845	-34.20516363	-31.84456068	-31.43979949	-34.2124325
-35.2081248	-34.24569054	-32.77682314	-30.69073947	-28.36930028	-29.55200392	-31.26279251
-40.8786813	-34.71636141	-31.72974163	-26.12673289	-25.33421844	-31.18890509	-33.93726891
-32.01451542	-28.20476352	-25.73050701	-26.02377508	-25.51602535	-27.36470247	-28.05968708

-75.04461904	-75.1468258	-76.52466806	-76.63285608	-76.47694965	-78.5838413	-82.72362928
-81.67107689	-77.72193717	-83.35510625	-84.7346017	-85.54500519	-86.16725094	-88.34514364
-80.61016457	-79.47063521	-80.94534176	-82.23925973	-83.92454828	-90.27588994	-81.55610197
-79.6338258	-78.74507363	-74.20506181	-76.32023148	-77.6178813	-83.08155306	-81.19137741
-74.99736151	-78.54960986	-77.99193681	-80.07770032	-78.85337915	-80.51923052	-81.81317494
-76.02866838	-81.13814121	-77.02773388	-76.72086173	-78.443404	-79.21451373	-80.25659777
-74.66385553	-73.44988022	-70.03522819	-73.99013833	-75.29149906	-81.67773401	-78.51251478
-72.64109982	-72.81946306	-71.42193013	-72.95863499	-70.19218131	-74.51426758	-76.91146863
-68.9625169	-69.26703708	-70.33283129	-71.21676881	-68.84893003	-66.77041751	-68.60373337
-69.64833108	-66.20137798	-70.16309883	-68.4263159	-69.31518694	-74.07032644	-71.24443301
-65.89129871	-67.82840005	-66.22651519	-66.3704488	-66.99911365	-68.62689015	-68.99018493
-62.5329376	-62.57694893	-55.88019736	-62.44290015	-64.71757345	-65.65926916	-65.99229642
-61.06079763	-58.88072135	-60.64708267	-61.59170202	-61.92251969	-61.86392255	-62.49575155
-59.23590014	-56.63067651	-60.45980695	-60.27401632	-60.58950752	-57.31160049	-56.76334354
-58.714123	-57.34020279	-59.76678295	-58.52795707	-58.4282286	-58.58826991	-65.32044278
-55.95890624	-55.12402702	-54.53109537	-55.98387625	-53.4140155	-57.15159617	-57.56985252
-55.28323086	-50.49444637	-53.35490546	-54.22993943	-52.86202051	-48.4717116	-47.51826418
-51.46161745	-51.81413696	-52.9019785	-53.28582778	-53.14409061	-53.39067453	-55.41051567
-49.35547082	-50.82709626	-50.90870167	-55.81860073	-54.95411087	-59.21301725	-52.51698673
-44.88023338	-46.63000797	-45.99575968	-49.7282881	-48.56385508	-49.30119947	-47.8395774
-44.30577218	-45.80207414	-44.62676679	-49.48254487	-44.31317255	-45.25618197	-36.86385603
-44.29969296	-42.69060614	-44.24977673	-46.45305974	-44.62824462	-46.15653625	-45.48560963
-42.78884257	-43.05126228	-42.99822287	-43.08237923	-41.24554552	-42.70550025	-42.98718745
-44.46473888	-41.32814665	-40.95974152	-41.12697711	-40.37993079	-40.333611	-39.35594894
-40.04459348	-39.47010112	-39.61452018	-40.73618251	-38.92628428	-35.80970629	-36.83189429
-34.16977753	-32.80639763	-38.11823078	-37.65472612	-31.30733801	-32.20831862	-37.34655611
-29.39608525	-34.89202351	-37.13819662	-36.41599079	-38.89019979	-34.92571801	-33.3745182
-31.21356866	-35.33720953	-34.44131758	-35.60471891	-34.54181219	-33.81608826	-33.63929098
-36.15988301	-33.76168239	-33.25100591	-33.61837766	-32.76022899	-34.52812335	-33.17388374
-29.40312172	-30.53583451	-28.68116631	-30.43043796	-30.19249001	-32.36826299	-29.2474361

-69.94467567	-62.23924486	-69.28608175	-72.07740385	-74.85158467	-73.78731631	-74.84226958
-80.77896077	-78.01148199	-78.61423686	-79.60326644	-81.14901774	-82.27793056	-82.94923238
-76.95642254	-71.36028887	-77.37487852	-78.84298168	-79.2827836	-80.68286926	-81.85314124
-74.76246169	-74.95931969	-76.41677602	-76.93025935	-76.3144652	-77.89386788	-78.70614054
-73.37853647	-74.7147925	-74.69929925	-74.35488012	-75.15756711	-75.91180167	-76.71482157
-71.37171695	-71.86308847	-73.24193455	-71.75480651	-72.59704165	-74.49878142	-75.58269487
-69.90346617	-70.87364636	-71.15391404	-66.06648931	-71.81041733	-73.39585749	-74.14941529
-67.71176885	-69.2112278	-68.87226604	-67.96541656	-71.55181382	-70.14560567	-60.93749252
-66.57677124	-66.88825264	-67.29912075	-66.47174836	-68.68217839	-68.78837065	-68.3350783
-64.93204552	-65.52913801	-65.41488992	-59.69476108	-65.39543716	-66.83566878	-67.37015845
-63.13401554	-63.9774429	-68.40293473	-63.62415854	-64.17095996	-65.2482643	-65.05924958
-60.60561471	-60.77751361	-62.30534768	-62.5754134	-62.28963591	-61.29486606	-56.25284652
-58.47527435	-58.04793871	-59.38375926	-56.98100929	-56.67616617	-53.5973143	-59.08319041
-56.50871934	-57.1215617	-57.46083986	-51.53136963	-56.45242499	-58.93641041	-61.17207507
-55.92899351	-54.63607067	-54.63231064	-54.20833511	-55.24376066	-57.42148268	-61.13520261
-52.83009389	-52.50395596	-52.57233905	-52.29968471	-45.9807716	-53.43191175	-55,42213293
-50.9288863	-51.20800602	-51.16941309	-52.4270149	-50.24215707	-51.93959852	-57.45129621
-49.33446022	-49.22583844	-46.30375768	-49.6340268	-48.11889239	-51.7419003	-48.38031523
-47.34905439	-47.51831459	-47.52997499	-49.81094712	-49.60003453	-49.42098601	-48.90156794
-45.20427452	-44.96658211	-46.97231229	-51.43864101	-48.31433293	-46.84176112	-46.24565333
-43.30566427	-42.67584984	-43.33549872	-44.88566403	-47.36771191	-45.79645688	-40.99184724
-42.78333031	-42.61364652	-39.76898385	-41.9062116	-43.61769334	-43.48341808	-43.51274553
-40.01200054	-40.66527618	-41.36441069	-41.61286575	-42.4221199	-37.08325005	-41.40153287
-38.94776684	-38.14801413	-38.90514405	-39.77523783	-38.53694498	-34.69188523	-39.1722635
-37.54438446	-36.62534402	-37.43586912	-38.24833153	-37.74601754	-37.37701216	-38.52713374
-35.39418362	-34.59742196	-35.22670309	-36.17280484	-37.10652218	-37.22395791	-36.84836443
-34.09983933	-34.28424984	-35.31951823	-35.3356109	-40.00524869	-38.24648242	-33.7421198
-32.27321456	-32.87046441	-35.21538878	-35.58586373	-39.05813787	-39.15580632	-27.76820418
-31.39255866	-31.31365489	-32.02603002	-32.49320024	-32.88880759	-33.37186744	-28.59941643
-27.18948772	-27.15913401	-27.79764561	-28.00509618	-27.3468977	-29.29235764	-27.5250328

-65.94223869	-75.43752102	-74.33406174	-61.87332735	-69.2560821	-69.64848335	-70.50226329
-75.03594205	-68.72024693	-62.48255265	-70.85321621	-62.13755657	-72.78951713	-76.89864839
-76.12708911	-65.45315027	-62.85517628	-71.330232	-71.21928775	-72.11337553	-76.46629231
-57.81637622	-58.58948791	-68.88194509	-79.92368721	-73.98733977	-73.47201662	-74.2718902
-71.13837343	-73.92595101	-64.69205369	-73.23318846	-65.64189818	-70.57717217	-72.10805528
-60.40837789	-76.77566403	-75.87699775	-68.846105	-66.7351857	-69.62506849	-70.85722716
-61.80927669	-57.07087842	-67.46690873	-62.92263947	-65.33276626	-68.10049899	-68.54818365
-60.00206229	-59.56757201	-62.47155602	-62.14183275	-65.34186376	-64.85439323	-66.73222733
-60.41381797	-57.83031224	-58.78734312	-52.56671633	-63.60065321	-64.78220489	-71.07281854
-57.38309794	-58.56614892	-59,48799111	-58.53334151	-60.21734361	-62.48060789	-64.76075651
-58.29374046	-58.95721373	-57.95462527	-58.95544209	-59.28157927	-59.15093116	-61.53966857
-56.06741794	-57.35154067	-56.23117791	-55.31553684	-55.97727207	-53.31371974	-59.17454819
-57.69575721	-57.12632097	-54.19685887	-53.04641249	-54.04776521	-57.20264405	-57.03659061
-51.68771554	-53.2724197	-53.26355882	-53.82171847	-52.97872201	-54.89603431	-56.39426188
-48.95555924	-52.4922446	-51.91286847	-53.40740774	-51.97513318	-52.78166864	-54.77977027
-45.37480333	-48.39677839	-49.78904268	-52.29017555	-52.21407758	-52.09772325	-52.12177395
-44.51751753	-48.19185816	-49.89388142	-49.80386	-49.69603789	-49.97646869	-50.75091269
-44.35126677	-45.2337149	-47.03326451	-47.84451166	-48.38344284	-48.70017318	-49.16779383
-43.50874179	-43.62413261	-46.08506097	-46.02838849	-46.01496708	-46.28549739	-46.35786167
-41.65899508	-41.56229715	-44.11544932	-44.01964081	-44.43166177	-44.96643982	-45.70295833
-41.78710969	-41.95744185	-42.41227306	-43.17775735	-43.58911584	-43.91838092	-43.16046149
-38.89962345	-40.63886677	-41.7568599	-43.2882356	-42.82380807	-42.26305898	-42.08615082
-39.30138318	-37.90333072	-41.09929555	-40.66895566	-39.76048203	-40.51533495	-39.65440756
-37.17432964	-38.62902161	-39.34564495	-37.58025147	-38.87803789	-38.61077356	-38.01532488
-35.31175175	-36.00151351	-35.56053899	-36.85849659	-37.41828396	-37.92059472	-37.50745865
-34.49069316	-34.67157336	-34.47546521	-35.81109172	-35.46305257	-35.82363828	-35.99181925
-33.4140639	-33.37448925	-33.72422678	-34.11466672	-34.43335481	-33.86427141	-33.98075458
-31.74125612	-32.06239365	-32.22852275	-31.64497517	-32.0024255	-32.53366287	-32.31942917
-30.89466593	-29.83964567	-30.03842331	-30.98286678	-31.56638157	-31.87785776	-31.1645448
-26.16684024	-26.20709125	-25.48127104	-26.91583832	-28.0297205	-27.33146518	-26.91460166

## 6.11 - Appendix K- HDM Displacement in the Y Direction Raw Data

0.1	11 -	Ap	hei	Iui	AI	Z- I	ושו	VI L	113	pra	cei	IIIC	IIL	111	tne	5 I	ווע	EC	uu	11 1	lav	V L	ata	1				
-13.8686	-10.3125	-10.5381	-9.20717	-13.9682	-10.2656	-6.84603	-8.48142	-7.84918	-7.85879	-6.69556	-5.79695	-5.26274	-4.97002	-3.95822	-4.1379	-3.96013	-3.38226	-3.18521	-2.41216	-0.73103	-0.25312	0.560649	1.240737	1.175112	2.035925	5.04946	9.141245	3.64635
-12.8247	-9.94431	-11.9733	-8.63036	-9.8227	-9.35375	-8.08807	-9.85813	-8.10843	-8.78618	-6.88046	-5.34663	-2.09776	-3.49556	-3.04951	-2.95339	-2.40721	-1.8183	-2.31853	-0.77197	-0.64695	1.120303	2.126958	2.258103	2.532634	2.023144	3.341913	3.40453	3.227135
-6.10189	-5.23469	-5.14548	-1.99062	-5.90798	-3.41991	-5.24291	-5.22365	-5.20993	-4.19196	-3.49292	-2.97455	-2.6525	-1.78926	-0.72026	-0.39142	-0.06052	1.06709	-1.39435	1.37382	1.20003	1.925389	1.853144	2.57713	3.761717	1.48041	3.072742	8.316503	4.098881
3.974051	4.371748	2.857015	2.352209	-1.51838	-0.96048	-0.97858	-1.17464	-0.73353	-0.54583	-0.32436	-1.11533	0.96524	0.362011	1.930578	1.142554	0.00044	0.773207	1.205572	0.657907	2.635939	2.677962	3.522363	3.28618	3.762169	4.186679	7.3852	23.38794	6.795436
7.520539	4.814772	5.620788	6.269172	2.818338	2.53394	2.445272	2.458011	1.47035	1.423761	2.387628	1.763717	2.145418	2.906505	5.352436	3.416503	2.348376	2.035794	2.978009	2.996636	3.503599	4.308516	4.364761	3.496906	4.447996	5.328058	6.774781	13.63366	4.170099
2.203809	11.51807	16.9828	6.302239	3.714629	2.847127	2.179552	3.656819	3.561404	3.566579	3.922229	2.927582	4.427657	3.164831	2.7716	8.069499	7.703695	3.697319	4.270134	4.25649	5.192456	13.61509	6.283296	-0.58377	5.024533	5.295314	5.479816	5.936012	5.413241
4.486088	5.678908	6.042693	5.061678	3.828929	0.863219	-2.06315	5.474016	3.884687	0.377872	1.683839	-0.66754	3.687271	3.447759	-2.12708	3.351539	4.717553	2.808106	5.175182	4.344862	6.346174	15.93729	5.791145	-6.02642	3.376241	5.343352	5.623661	6.130933	4.937178
5.951777	5.653534	4.019557	5.521701	5.229022	2.325402	3.168402	9.877633	4.368208	3.162151	-0.52325	4.154233	9.053573	8.623567	-2.52796	-0.42701	-4.18516	-6.42966	4.221363	5.997669	6.247692	2.895267	4.880663	4.448981	7.093413	7.693639	7.042612	7.569729	6.521759
7.38725	7.583588	11.92402	7.744074	5.849196	5.88053	1.279457	2.741031	-2.61614	5.787688	5.19915	5.806366	10.44246	8.173553	3.33328	3.25976	3.108751	7.52517	7.029084	9.345816	4.290196	4.156875	6.464459	8.036433	7.725077	8.996005	7.287658	7.135771	7.943734
10.11806	8.225746	8.676324	7.979337	5.881861	6.24731	6.758629	7.1185	7.898377	7.594442	7.589897	8.078447	8.387685	7.098137	6.718791	5.165087	7.560472	21.56106	5.711545	6.384914	5.062438	7.451439	6.944301	7.448509	10.69385	8.92892	7.617126	6.048922	8.017322
5.578534	9.111347	9.310844	9.032394	8.596685	8.813399	9.788691	8.259836	8.835851	8.222374	7.710182	8.844604	8.574807	9.36437	8.825836	8.110475	7.66968	8.548996	8.830602	9.866064	8.23705	12.08987	8.495399	5.875299	10.91719	8.718493	9.106809	7.866667	5.969036

8.739087	9.128915	10.40626	10.63593	9.01236	9.900376	10.52134	10.41992	10.79479	12.88716	10.37933	9.998334	10.1043	8.409382	7.242405	8.225859	9.22237	8.148393	7.163153	8.268692	4.5181	12.34754	10.04198	9.792505	6.508197	9.556765	9.901737	9.351534	9.39791	2.664813
10.10549	11.70269	11.24542	5.836011	9.934063	10.88456	18.38571	14.70519	11.90527	16.30462	11.75053	11.29407	9.988273	-0.72928	11.06257	9.332639	9.7728	10.00667	8.722466	10.50673	10.24652	9.550061	8.771812	9.158105	8.553829	8.565219	8.607778	7.071555	9.196481	7.562437
10.97236	12.76003	12.54043	11.40851	11.53291	5.79216	11.37344	10.47237	10.64053	10.53277	15.50536	12.49418	16.97423	-3.33939	9.767163	11.21626	9.938681	10.84062	10.90071	11.18026	10.52692	4.588039	8.740915	9.729794	4.904247	5.259608	7.709056	4.859317	9.451813	9.605906
11.33787	12.88423	13.11769	12.79406	15.93496	11.18441	9.611108	2.460381	7.353419	14.34036	12.8484	6.701585	12.51539	8.742422	10.72321	12.69117	15.12101	9.414287	10.90572	12.42263	11.00036	9.948907	10.93602	12.54713	11.0593	16.6998	7.987068	5.834519	10.74482	8.579167
10.69978	13.29767	14.95553	13.94163	13.78731	13.22787	10.97272	5.691188	12.18533	13.47522	13.70318	14.37018	12.68457	5.123141	6.395582	14.295	15.55956	7.993024	11.96941	13.52543	15.07443	12.0259	12.13375	15.02483	11.60751	11.46835	10.49611	11.2751	14.41306	8.601415
6.79789	12.90302	13.73843	14.49423	13.71186	14.08347	13.936	13.06685	14.44662	14.07483	14.5382	14.3027	12.39218	12.56165	12.81356	14.14717	13.66226	9.978376	13.12641	13.59492	13.54545	13.41695	16.06165	11.12403	11.03364	11.64888	16.74336	15.11092	10.15321	8.549698
7.964413	13.36172	14.36814	16.5479	15.21547	13.14264	15.18951	15.00123	16.24579	16.90367	16.02988	15.33354	19.85693	17.16986	14.64068	11.34313	14.3927	14.45899	14.98172	13.96444	9.050978	13.34541	14.70485	13.94156	12.12629	13.01866	14.32185	18.48409	12.1026	10.38238
13.4069	15.59899	15.3589	16.48631	16.41909	15.85737	18.18809	17.55638	17.61221	23.69872	17.49467	16.5151	16.26739	19.0952	17.18838	15.68469	10.40148	14.76285	16.92805	20.4415	15.82421	15.2044	14.05485	14.61349	13.42169	13.82317	12.65673	9.736348	10.89416	9.677025
14.07651	16.86628	17.27907	16.38277	16.3796	17.41491	17.23569	17.02185	16.70995	17.7572	17.56865	17.3085	17.58327	18.05433	17.70689	16.98999	16.13905	16.08448	16.9719	16.62061	16.05856	17.57107	14.55438	15.03405	14.60837	14.24006	13.40871	12.12899	11.93936	9.161062
15.95152	16.54615	20.29025	17.53072	17.19092	18.29306	18.10383	17.28772	18.10741	17.45921	17.98158	18.2587	18.70398	18.47928	18.7634	18.47415	17.93044	17.04213	18.33139	17.54145	16.96687	17.63617	16.0646	15.80409	16.23583	16.72368	14.58431	13.27356	12.36539	9.650677
14.96673	15.74525	17.61362	18.13952	18.3158	18.45005	18.70634	18.09425	17.77284	19.00631	18.88362	19.5389	18.36973	19.3995	18.51793	19.38428	19.02918	18.81421	18.98626	18.94763	18.34216	17.26533	17.34867	16.74735	15.78708	15.70192	14.83898	15.47732	12.73884	9.941836

			20.000.02	20.88594	17.10331	16.18361	15.5157
26.16414	37.9555	24.01172	25.3793	23.52698	27.13143	21.77604	18.72819
39.03829	29.28541	24.11309	21.72837	20.73539	21.53715	16.9967	18.56234
32.2487	34.2428	29.39697	15.65143	22.05154	23.48114	18.83664	18.46935
23.98531	45.15818	28.96579	17.6142	15.45846	22.77178	20.57837	19.67672
25.05095	37.10433	21.50564	15.61843	19.11754	21.69138	20.36992	19.05587
26.48727	30.44066	19.36048	15.72682	21.39755	21.83017	20.06689	19.70967
21.50392	28.45147	21.34474	20.98997	20.29209	19.65946	18.20554	18.06862
25.27871	27.06885	20.84194	21.60309	19.78075	16.42486	19.26816	12.68475
24.39493	24.48839	16.36532	19.88606	23.76523	28.29958	22.63069	19.58691
23.26909	17.92034	20.57471	21.43951	22.96803	23.25658	22.51254	20.62002
18.49334	23.1017	21.84649	22.78982	23.41327	23.50916	20.71767	20.54063
25.09584	31.96971	22.55582	22.80326	23.46693	21.82269	21.51797	21.35453
22.45407	24.3462	23.37788	21.31689	22.2943	21.58397	22.04908	20.47421
19.03598	20.622	21.43338	22.46559	22.20412	22.44789	21.93323	20.67213
23.096	22.26983	22.03812	22.59165	21.94695	21.05286	21.31887	21.31363
20.14205	21.9247	21.23977	20.82641	21.50866	20.89551	21.4975	19.59295
20.40959	20.59006	21.00316	20.79017	20.40716	20.69101	20.78094	19.6027
17.28652	18.84781	21.32736	20.41403	20.66515	20.42205	20.18548	19.77941
17.29673	18.98485	20.61573	19.98268	19.70125	19.7112	20.3513	19.39728
17.63051	20.35542	20.67251	19.03392	18.87612	19.59455	18.93278	18.47953
16.89547	19.10732	18.98508	19.11654	18.14253	18.90343	18.52353	17.90474
15.16398	18.35021	18.18092	19.10727	18.21503	18.77337	18.38461	17.74325
14.37823	17.77122	18.38124	18.78659	18.58542	18.37929	17.71849	15.53271
15.01206	17.8158	17.91277	17.45905	17.34003	17.45961	16.97782	14.91944
13.49685	15.62763	15.25792	15.1612	15.92001	16.78139	16.97369	13.80916
12.36746	15.08628	13.40964	14.23103	15.18396	14.88674	14.99643	14.59213
11.36446	13.67686	13.94206	13.85086	12.91579	13.29031	13.7754	14.41434
10.66007	12.76862	13.34058	13.4576	13.44472	13.43155	13.36086	13.55883
7.999584	10.11644	10.17968	10.83321	10.56566	10.59339	11.30213	11.17875

## 6.12 - Appendix L- HDM Strain in the X Direction Raw Data

U.	14	- <i>I</i>	Jh	JEI	lab	Y L.	. 11	ואוט	J	ıaı	11 1	II U	ile.	A L	111 6	Cu	UII	Ka	LVV	Dai	la							
0.004283	-0.3683	-0.26677	0.228013	-0.07778	0.261096	0.272026	0.13546	0.134312	0.191271	0.151698	0.116115	0.127787	0.196873	0.174653	0.172559	0.101028	0.124847	0.14349	0.131802	0.132232	0.13806	0.177508	0.078255	0.11364	0.087049	-0.10762	0.07341	0.528352
-0.04778	-0.21772	0.133306	0.196878	0.158747	0.226959	0.251199	0.38394	-0.02624	0.133807	0.174057	0.097959	0.214211	0.265432	0.12394	0.184709	0.146939	0.150298	0.177412	0.077998	0.115549	0.167761	0.112361	0.160416	0.102318	0.159655	0.184352	-0.16715	0.367349
-0.24951	0.057732	0.168409	0.132311	0.25387	0.303196	0.132257	0.250711	0.253919	0.049033	0.205128	0.128225	0.16055	0.181927	0.131103	0.185169	0.126166	0.145751	0.109937	0.168223	0.177547	0.073999	0.135346	0.158997	0.203038	0.227549	0.117987	-0.10231	0.239013
-0.5221	0.353182	-0.09017	0.383511	0.3341	-0.01344	0.194138	0.155854	0.216261	0.129239	0.133155	0.213238	0.299046	0.09634	0.165616	0.079768	0.152634	0.138833	0.164154	0.325115	-0.06349	0.08883	0.259453	0.025039	0.137326	0.284136	0.217204	0.189693	-0.01136
-0.57005	0.155193	0.122313	0.392752	-0.02804	-0.13363	0.394278	0.206696	0.194923	0.165869	0.193086	0.184709	0.136875	0.18797	0.371667	-0.10263	0.10549	0.107096	0.209862	0.155725	0.068711	0.06761	0.242443	0.052714	0.177372	0.155615	0.219652	0.28525	0.006435
-0.68956	-0.15422	0.079315	0.429917	0.141336	0.097495	0.198083	0.18328	0.16889	0.159914	0.180272	0.244344	0.100314	0.145687	0.107942	-0.26465	0.409655	0.278655	0.093703	0.192381	0.048301	-0.14349	0.366107	-0.1821	0.540553	0.070876	0.171275	0.065443	0.374952
-0.39996	-0.17507	-0.03136	0.413383	0.160012	0.094045	0.438502	0.118434	-0.0122	0.272647	0.085193	0.606049	-0.25511	0.106845	0.264499	-0.07927	0.056215	0.271849	0.196603	0.156903	0.001457	0.195573	0.051112	-0.06128	0.840041	-0.33811	0.021461	-0.02942	0.406975
-0.41465	0.068249	-0.09981	0.348383	0.15066	0.090489	0.365958	0.084392	0.097822	0.160583	0.1162	0.247678	0.104245	-0.01675	0.19734	0.088017	0.186105	-0.10965	0.33216	0.303953	0.022313	0.104446	0.092455	0.215455	0.149951	0.038165	0.011404	-0.35441	0.55401
-0.35134	0.424315	0.022795	-0.03886	0.097286	0.109005	0.100065	0.519233	-0.16504	0.140891	0.187368	0.218534	0.358276	-0.53482	0.484412	0.628224	-0.49327	0.180846	0.292338	0.685983	-0.53886	0.156151	0.226952	0.157753	-0.03217	0.248252	-0.01655	0.029088	0.245403
-0.47396	-0.25679	0.449646	0.160145	0.081545	-0.15395	0.447717	0.483991	-0.45624	0.340215	0.185476	0.237209	0.28452	-0.07979	0.089792	0.542493	-0.30744	-0.3639	0.619489	0.252814	-0.05627	0.21569	0.148243	0.282744	0.225087	-0.16984	0.069352	-0.0445	0.134991
-0.56675	0.101279	0.394167	-0.07722	0.025623	0.196994	0.318707	0.083953	-0.02914	0.144755	0.142596	0.174691	0.083313	0.13508	0.313388	0.0345	-0.01763	-0.11313	0.399391	0.265668	-0.01969	0.211419	0.054101	0.090853	0.476184	-0.47393	0.271774	0.111349	0.160484

-0.50636	0.155959	0.369939	-0.23484	0.209802	0.17067	0.064469	0.108867	0.174403	0.128492	0.245472	0.0532	0.082355	0.109129	0.159005	0.109621	0.059007	-0.1583	0.380645	0.015359	0.189343	0.210668	0.122213	0.024425	0.192591	0.077421	0.050704	0.124146	0.199246
-0.4269	0.15061	0.421267	-0.23668	0.060263	0.437032	-0.08667	0.068069	0.010608	0.246036	0.646645	-0.29793	0.011705	0.043314	0.32723	0.073512	0.028308	0.12458	0.307059	0.085562	0.023562	0.078222	0.127405	0.084076	0.093518	0.061252	0.168555	0.074394	0.285615
-0.16094	-0.10929	0.045348	0.012216	-0.16178	0.480516	0.039401	0.222027	0.191604	-0.10169	0.328216	0.231014	0.140628	-0.04435	0.138511	0.289349	-0.08248	0.06169	0.262318	0.051746	0.194467	-0.02254	0.107695	0.116128	0.416481	-0.13035	-0.02782	0.09847	0.201615
-0.41415	0.066307	0.061021	0.289779	-0.06446	0.085301	0.126422	0.229911	-0.04286	0.234815	0.209898	0.092009	0.114056	0.032611	0.172201	0.04223	0.238851	0.131634	0.279702	0.035904	0.00038	0.094428	-0.10474	0.276259	0.367176	0.298356	-0.11359	-0.30914	0.422298
-0.50669	0.068506	0.196688	0.124457	0.070758	0.08958	0.825745	-0.46235	0.060307	0.144432	0.5504	-0.1769	-0.13056	0.002305	0.357067	-0.12682	0.566936	-0.03258	0.165995	0.328363	-0.15756	0.131951	0.139329	0.040321	0.104923	0.19414	0.37337	-0.05195	0.067149
-0.53066	0.099691	0.174313	0.123879	0.088314	0.068933	0.203141	0.084827	0.122044	0.099213	0.247087	0.481097	-0.33369	0.094683	0.249348	0.09327	0.012356	0.145057	0.161202	0.065332	0.144565	0.400011	0.14946	-0.16782	0.009566	-0.06391	-0.05683	0.361496	0.254969
-0.39359	0.11664	0.18552	0.072306	0.160033	0.049164	0.016163	0.179352	0.205421	0.07653	0.117583	0.350842	0.013984	0.075542	0.578937	-0.26634	0.132704	-0.09257	0.080356	0.059164	0.234376	0.074723	0.242823	0.049433	0.039968	-0.18117	0.059194	0.385583	0.346369
-0.47037	0.047518	0.119545	0.160961	0.162505	0.35552	-0.11868	0.093354	0.423562	-0.24559	0.065547	0.34965	0.340602	-0.16731	0.119291	-0.00796	0.174562	-0.01106	-0.10173	0.409561	0.186216	0.018334	0.114852	0.095432	0.12972	0.052325	-0.01564	0.193291	0.280507
-0.58301	0.07746	0.059881	0.107342	0.091085	0.130501	0.142603	0.098322	0.117764	-0.18675	0.381099	0.182599	0.120182	0.176783	0.128748	0.087683	0.304103	-0.07664	0.034854	0.227301	0.222907	-0.09971	0.153704	0.09183	0.138073	-0.0058	0.006508	0.199335	0.264274
-0.98576	0.4157	-0.22494	0.015283	0.178232	0.06184	0.103901	0.145186	0.084945	0.096981	0.199996	0.170598	0.057899	0.155343	0.133257	0.080997	0.123885	0.10672	0.159483	0.143171	0.003888	0.121773	0.157329	0.095167	0.126745	0.019573	0.088362	0.097301	0.259658
-0.67714	0.238909	0.137123	0.086495	0.125426	0.091766	0.136981	0.070937	0.102795	0.112377	0.158025	0.133146	0.12291	0.036233	0.193681	0.118825	0.099652	0.124088	0.134049	0.118663	0.032646	0.173208	0.066515	0.087711	0.134388	0.080897	0.114164	0.055041	0.262692

31 -0.39977	59 0.027022	92 0.13715	28 0.13524	06 0.078177	86 0.144315	82 0.113497	12 -0.27129	85 0.394504	05 0.201318	26 0.14782	06 0.133622	63 0.040146	48 0.100906	47 0.166125	78 0.085679	68 0.098945	17 0.175621	41 0.040931	04 0.158906	58 0.067144	33 0.151984	35 0.102443	36 0.031742	06 0.094727	46 0.125692	63 0.103833	88 0.07218	15 0.265621
-0.19631	0.042259	-0.08492	0.180928	0.059506	0.095286	0.202882	0.004512	0.14385	0.208105	0.364826	-0.24306	0.144163	0.132148	0.042747	0.132578	0.079768	0.150917	0.082441	0.065504	0.103458	0.109233	0.119035	0.043136	0.13106	0.12246	0.083163	0.040988	0.28415
0.444908	-0.56761	-0.173	0.52159	-0.06833	0.087651	-0.00057	0.108826	0.211457	0.058485	0.206519	0.120594	0.066815	0.062724	-0.01493	0.157377	0.082037	0.14803	0.098957	0.052659	0.047832	0.191458	0.055153	0.091235	0.122202	0.064356	0.151933	0.027253	0.221041
-0.56124	-0.02981	-0.53709	0.418156	0.274193	0.370217	0.0488	0.598445	-0.37291	-0.02638	0.227494	0.14182	-0.04846	0.025894	0.069827	0.155395	0.122459	0.113508	0.125547	0.052618	6900'0-	0.163705	0.193044	0.04511	0.065463	0.106027	0.154356	0.041382	0.254189
0.740719	-0.02329	-0.37667	0.261868	90669:0-	0.525631	0.31221	0.230263	-0.04379	0.095835	0.107715	0.127145	0.058331	0.084418	0.132739	-0.00655	0.178789	0.059263	0.123101	0.106449	0.040963	0.041098	0.109603	0.236569	0.067817	0.046952	0.093482	0.136881	0.284822
0.41983	0.204194	0.428979	-0.95853	-0.17811	1.231549	-0.15604	0.108579	-0.04599	-0.02444	0.100355	0.014076	0.240869	0.048761	0.255967	0.012808	0.184884	0.100599	0.128865	-0.0247	0.082411	0.170971	-0.04536	0.164219	0.083121	0.081068	0.082006	0.138922	0.227035
-0.56836	-0.0682	1.14442	-0.83262	0.670625	-0.08756	0.112951	-0.02573	0.18942	-0.05692	0.139145	-0.10177	0.375503	0.17076	0.223797	0.05358	0.010391	0.052658	0.115609	-0.00801	0.180468	-0.02511	0.132941	0.116411	0.051316	0.067289	0.10455	0.052912	0.295489
-0.47558	-0.06537	0.743414	-0.29736	-0.16772	-0.15811	0.76464	-0.13131	0.202451	0.051388	-0.30963	0.29195	0.255608	0.191204	0.030909	0.115626	0.044964	0.145255	-0.04451	-0.00886	0.10017	0.184214	0.076552	0.031342	0.105019	0.038814	0.137701	-0.0382	0.295863

# 6.13 - Appendix M- HDM Strain in the Y Direction Raw Data

0.1		Ap	PC		21.			1.1	JUI	u I I I		CII	<b>C</b> I			CHI		Na	** .	Dai	·u							
0.065241	0.023013	-0.0897	0.036051	0.259095	0.056991	-0.07763	-0.08604	-0.0162	-0.05796	-0.01156	0.028145	0.197811	0.092154	0.056795	0.074032	0.097057	0.097747	0.054167	0.102512	0.005255	0.085839	0.097894	0.063585	0.084845	-0.0008	-0.10672	-0.35854	-0.0262
0.420176	0.294351	0.426739	0.414983	0.24467	0.370865	0.177822	0.289655	0.181156	0.287139	0.211722	0.148255	-0.03467	0.106644	0.145578	0.160123	0.146668	0.180337	0.057762	0.134112	0.115436	0.050318	-0.01711	0.019939	0.076818	-0.03392	-0.01682	0.306998	0.054484
0.629747	0.600402	0.500156	0.271427	0.27435	0.153714	0.266521	0.253063	0.279775	0.227883	0.198035	0.116201	0.226109	0.134454	0.165677	0.095873	0.00381	-0.01837	0.162495	-0.04474	0.089744	0.047036	0.104326	0.044316	2.82E-05	0.169142	0.269529	0.941965	0.168535
0.221656	0.027689	0.172736	0.24481	0.271045	0.218401	0.213991	0.227041	0.137743	0.123099	0.169499	0.179941	0.073761	0.159031	0.213866	0.142122	0.146746	0.078912	0.110777	0.146171	0.054229	0.10191	0.05265	0.01317	0.042864	0.071336	-0.03815	-0.60964	-0.16408
-0.3323	0.418956	0.710126	0.002067	0.056018	0.019574	-0.01661	0.074925	0.130691	0.133926	0.095913	0.072742	0.14264	0.016145	-0.1613	0.290812	0.334707	0.103845	0.080758	0.078741	0.105554	0.581661	0.119908	-0.25504	0.036034	-0.00205	-0.08094	-0.4811	0.077696
0.142642	-0.12082	-0.68376	-0.07754	0.007144	-0.12399	-0.26517	0.113575	0.020205	-0.19929	-0.1399	-0.2247	-0.04627	0.017683	-0.30617	-0.29487	-0.18663	-0.05558	0.056565	0.005523	0.072107	0.145137	-0.03076	-0.34017	-0.10302	0.003002	0.00899	0.012183	-0.02975
0.091606	-0.00159	-0.12645	0.028751	0.087506	0.091386	0.326972	0.275226	0.03022	0.174017	-0.13794	0.301361	0.335394	0.323488	-0.02505	-0.23616	-0.55642	-0.57736	-0.05961	0.1033	-0.00616	-0.81513	-0.05691	0.654712	0.232323	0.146893	0.088684	0.089925	0.099036
0.089717	0.120628	0.494029	0.138898	0.038761	0.222195	-0.11806	-0.44604	-0.43652	0.164096	0.35765	0.103258	0.086805	-0.02813	0.366328	0.230423	0.455869	0.872177	0.175483	0.209259	-0.12234	0.078851	0.098987	0.224216	0.039479	0.081398	0.015315	-0.02712	0.088873
0.170675	0.040135	-0.20298	0.014704	0.002042	0.022924	0.342448	0.273592	0.657158	0.112922	0.149422	0.142005	-0.12842	-0.06721	0.211594	0.119083	0.278233	0.877243	-0.08235	-0.18506	0.048265	0.20591	0.02999	-0.03675	0.185548	-0.00419	0.020592	-0.06793	0.004599
-0.28372	0.05535	0.039657	0.065816	0.169676	0.160381	0.189379	0.071333	0.058592	0.039246	0.007518	0.047885	0.011695	0.14164	0.13169	0.184087	0.006825	-0.81325	0.194941	0.217572	0.198413	0.289902	0.096944	-0.09833	0.013959	-0.01315	0.093105	0.113609	-0.12802
0.197535	0.080932	0.082818	-0.00125	0.081481	0.106746	0.039452	0.158435	0.253207	0.13481	0.14301	0.078731	-0.01034	-0.13262	-0.0375	0.069493	0.02992	-0.08662	-0.03512	-0.33425	0.256906	-0.12799	0.081069	0.039556	-0.08503	0.073953	0.015295	0.095703	-0.20651

0.0854	0.160861	0.052448	-0.29999	0.057606	0.061511	0.491523	0.267829	0.069405	0.213591	0.0857	0.080983	-0.00725	-0.57117	0.23876	0.069174	0.034402	0.116142	0.097457	0.139877	0.358026	-0.17484	-0.07939	-0.03965	0.127852	-0.06197	-0.08087	-0.1425	-0.01259	0.306102
0.054179	0.066084	0.080938	0.348281	0.099928	-0.31827	-0.43827	-0.26455	-0.07905	-0.36074	0.234676	0.075007	0.436623	-0.16313	-0.08096	0.117726	0.010368	0.052122	0.13614	0.042096	0.017525	-0.31013	-0.00193	0.035731	-0.2281	-0.2066	-0.05617	-0.13826	0.015958	0.127717
0.022844	0.007762	0.036079	0.086597	0.275128	0.337016	-0.11015	-0.50075	-0.20544	0.237974	-0.16606	-0.36204	-0.27868	0.755113	0.059753	0.092182	0.323895	-0.08915	0.000313	0.077649	0.02959	0.335054	0.137194	0.176083	0.384691	0.715012	0.017376	0.06095	0.080813	-0.06417
-0.03988	0.02584	0.114865	0.071723	-0.13423	0.127716	0.085101	0.201925	0.301994	-0.05407	0.053424	0.479287	0.010574	-0.22621	-0.27048	0.100239	0.02741	-0.08883	0.06648	0.068925	0.254629	0.129812	0.074859	0.154857	0.034263	-0.32697	0.156815	0.340036	0.229265	0.00139
-0.24387	-0.02467	-0.07607	0.034537	-0.00472	0.053475	0.185205	0.460979	0.141331	0.037476	0.052189	-0.00422	-0.01827	0.464907	0.401123	-0.00924	-0.11858	0.124085	0.072313	0.004343	-0.09556	0.086941	0.245494	-0.2438	-0.03587	0.011283	0.390453	0.239739	-0.26624	-0.00323
0.072908	0.028669	0.039357	0.128354	0.093976	-0.0588	0.078344	0.120899	0.112448	0.176802	0.09323	0.064427	0.466547	0.288013	0.114195	-0.17525	0.045652	0.280038	0.115957	0.023095	-0.2809	-0.00447	-0.0848	0.176095	0.068291	0.085611	-0.15134	0.210823	0.121837	0.114543
0.340155	0.139829	0.061922	-0.00385	0.075226	0.169671	0.187411	0.159697	0.085401	0.424691	0.09155	0.073848	-0.22435	0.120333	0.159232	0.271347	-0.24945	0.018991	0.121646	0.404816	0.423327	0.116187	-0.04063	0.041996	0.080962	0.050282	-0.10407	-0.54673	-0.07553	-0.04408
0.041851	0.079206	0.12001	-0.00647	-0.00247	0.097346	-0.05952	-0.03341	-0.05639	-0.37135	0.004623	0.049587	0.082242	-0.06505	0.032407	0.081581	0.358599	0.082602	0.002741	-0.23881	0.014647	0.147917	0.031221	0.026285	0.074168	0.026056	0.046999	0.14954	0.065325	-0.03225
0.117188	-0.02001	0.188199	0.071747	0.050708	0.054884	0.054258	0.016616	0.087341	-0.01862	0.025808	0.059388	0.070045	0.026559	0.066032	0.09276	0.111962	0.059854	0.084968	0.057552	0.056769	0.004068	0.094389	0.048128	0.101716	0.155226	0.073475	0.071536	0.026627	0.030601
-0.06155	-0.05006	-0.16729	0.03805	0.070305	0.009812	0.037657	0.050408	-0.02091	0.096694	0.056377	0.080013	-0.02089	0.057514	-0.01534	0.056883	0.068671	0.110755	0.04093	0.087886	0.085956	-0.02318	0.080254	0.058953	-0.02805	-0.06386	0.015917	0.137735	0.02334	0.018197
0.021685	0.186434	0.059295	0.020614	0.085058	0.037863	0.062708	-0.0016	-0.31801	0.036287	0.108525	0.062608	0.18655	0.067169	0.134638	0.120585	0.035236	0.049281	0.049571	0.028103	0.008585	0.039963	0.024662	-0.07591	-0.05423	-0.1183	-0.01543	-0.06644	0.05125	0.077307

0.05437	0.190491	-0.09785	0.022955	0.056353	0.082128	0.022326	0.008557	0.411463	0.190237	0.118282	0.011065	0.010215	0.098429	0.078818	0.000328	0.119034	0.07364	0.02538	0.059626	0.028328	0.038674	0.040085	0.136611	0.128649	0.197783	0.025269	-0.03993	-0.01237	0.007711
0.062494	0.334712	0.283778	0.290282	0.137088	0.082591	0.110205	0.09087	-0.17771	0.354306	0.046503	0.174468	0.019045	-0.02907	0.032167	-0.01663	-0.03762	-0.00562	0.014786	-0.04001	0.041361	0.023744	0.024297	0.0413	0.030112	-0.01202	-0.00686	-0.03032	0.004418	-0.0443
0.231401	-0.22528	-0.05011	-0.08935	-0.45708	-0.16086	-0.02704	0.039539	0.209743	-0.2834	-0.01803	-0.00599	0.102765	0.044396	-0.01524	0.05588	0.038322	-0.01774	0.015194	-0.00062	-0.0449	-0.04756	-0.0349	0.012883	-0.00747	-0.05384	0.018576	-0.02341	0.000823	-0.00173
0.29273	0.11577	0.062061	-0.40001	0.134734	-0.21869	-0.35442	0.043617	0.113896	-0.24245	-0.09553	-0.03897	-0.04148	-0.06109	0.016342	0.040294	-0.04264	0.023938	-0.0157	0.017589	0.009863	0.060876	0.055765	0.012573	0.007439	-0.04743	-0.05956	0.058441	0.000805	0.016722
0.278524	-0.08547	0.149044	0.859096	0.709474	0.367951	0.227104	0.022173	-0.04757	-0.22005	-0.05405	-0.05896	-0.01547	0.128812	-0.06451	-0.0346	0.025835	0.013312	0.057083	0.039566	0.102412	-0.00822	-0.0579	-0.02533	0.028357	0.006045	-0.05134	0.0057	-0.00731	-0.04085
0.01065	0.871486	0.323271	0.302865	1.012025	0.974918	0.692511	0.444171	0.389182	0.507692	-0.1659	0.078451	0.588368	0.06052	-0.05071	0.014482	0.042808	-0.02582	-0.15497	-0.10193	-0.01982	0.00764	0.010581	-0.03813	-0.00606	0.023107	0.10479	-0.01657	-0.03575	-0.00395
-0.62581	-0.73696	0.609555	-0.12463	-1.3233	-0.75334	-0.24709	-0.43422	-0.11188	-0.00584	0.334297	-0.28802	-0.42962	-0.11826	-0.09913	0.051635	-0.11142	-0.01128	-0.09758	-0.10551	-0.17031	-0.13824	-0.19914	-0.21206	-0.17523	-0.13317	-0.16993	-0.14453	-0.13178	-0.1323

## 6.14 - Appendix N- HDM Shear Strain Raw Data

U	.14	- 1	App	Jei	lul	X IN	- 11	ואוש	1 31	IEa	11 3	ua	Ш	Ka	VV I	Jai	a											
-0.60388	-0.46184	-0.5332	-0.28452	-0.49046	-0.10771	-0.13362	-0.30201	-0.1069	-0.20723	-0.19932	-0.19641	-0.21687	-0.18121	-0.12445	-0.18704	-0.1698	-0.13434	-0.13352	-0.09855	-0.09708	-0.14302	-0.11767	-0.15443	-0.13665	-0.11336	-0.00978	0.169899	-0.24997
0.077954	-0.16199	-0.06314	0.12229	-0.05172	0.047756	0.110772	-0.04357	-0.00019	0.064032	0.102378	0.106292	0.17502	0.002982	0.01885	0.011495	0.02579	0.016738	-0.01957	0.010654	0.011343	0.093662	0.023012	0.007145	0.010917	0.036776	0.127856	0.055419	0.080343
0.031915	0.064401	0.116975	0.083488	-0.01192	0.228366	-0.06467	0.023839	-0.02376	-0.01121	0.018927	-0.02271	0.013666	0.099826	0.063463	0.057589	0.004953	0.043084	-0.07253	0.118001	0.104511	0.012091	-0.00542	0.081517	0.028932	-0.11223	0.037112	0.200827	0.051158
0.258439	0.00764	-0.03149	0.106306	0.005746	-0.03694	-0.11503	-0.02052	0.024811	0.006223	0.025594	0.02392	0.099391	-0.06556	0.048121	0.077503	-0.02476	0.011504	-0.015	-0.02277	-0.02854	-0.02294	-0.00848	-0.05075	-0.01466	0.003757	0.026185	0.427541	-0.54328
0.102053	0.083242	-0.02447	-0.05089	-0.16041	0.023231	0.14491	0.049982	0.007011	0.023403	0.052002	-0.00402	0.05722	0.050795	0.082305	-0.18649	-0.24038	-0.06469	0.060304	-0.02664	0.00695	0.006056	-0.12289	-0.08994	-0.1505	0.028869	0.004208	0.14915	-0.47083
0.033785	0.286085	0.247146	-0.31273	-0.06809	-0.005	-0.00048	0.186756	0.105187	0.017786	0.085105	-0.00463	0.254181	-0.00987	-0.00212	0.25401	0.169712	-0.12078	0.018921	0.052045	0.063981	0.274517	-0.04828	-0.19125	0.259012	0.241959	0.034771	-0.03165	-0.10967
-0.03857	0.071252	0.127059	0.050808	0.010441	-0.04839	-0.04894	0.241776	-0.06045	-0.06536	0.029009	-0.06978	-0.0394	-0.00329	-0.23182	0.080023	0.035148	-0.00227	-0.05938	-0.09152	0.070498	0.318107	-0.34425	-0.37581	0.425695	-0.15171	-0.01629	-0.01422	-0.22988
-0.0124	0.000584	0.141547	0.300853	0.051142	-0.05714	0.069203	0.119576	-0.04484	-0.04179	-0.12911	0.16781	0.160171	0.120644	-0.47344	0.084235	0.171242	-0.12114	0.427093	0.129847	0.273166	-0.12	0.072663	0.064378	0.131656	-0.02328	0.042658	0.065498	0.208026
0.125247	0.078316	-0.13685	-0.18969	-0.01878	0.033544	-0.2427	0.120587	-0.11012	0.174311	-0.00704	0.029383	0.164623	-0.08804	0.059122	0.010773	-0.03451	0.201133	-0.22475	0.026722	-0.42025	-0.02513	0.080915	0.018572	0.022213	0.200286	-0.10186	-0.01027	-0.01707
-0.00966	0.035809	0.21256	0.148959	-0.01349	0.035517	0.215549	0.146311	-0.04058	0.1391	0.050726	0.044695	0.007833	-0.14273	-0.00685	0.068246	-0.06234	0.445224	-0.36221	0.044085	-0.01186	0.122415	0.029776	0.014309	0.004024	-0.027	-0.16488	-0.07169	0.116759
0.016545	0.114309	0.0589	0.03185	-0.05188	0.060602	0.071145	-0.13423	-0.05599	909800.0	0.00364	0.106534	0.001907	0.034533	-0.01995	-0.10266	-0.05652	0.023049	-0.01822	-0.00403	-0.21245	0.063375	-0.16973	-0.10522	0.101	-0.26707	0.089454	-0.07197	-0.08612

0.015563	0.083026	0.047612	0.015363	0.092931	0.009815	0.120421	0.059737	0.09301	-0.13264	-0.00741	0.208396	-0.02345	-0.04227	-0.00798	0.07654	-0.00622	-0.01879	0.187982	-0.00056	0.396413	-0.0032	-0.00517	-0.09741	0.13032	-0.0037	-0.03976	0.037806	-0.19893
0.09297	0.161747	-0.12296	-0.01381	0.012276	0.105961	-0.22172	-0.13117	0.170786	-0.01851	-0.06432	-0.25008	-0.27972	0.488156	0.02177	-0.00477	0.096698	-0.00614	0.058308	-0.02795	-0.05849	0.024404	0.010414	-0.0304	0.004869	0.167325	0.022186	0.038407	-0.06702
0.059059	-0.13027	-0.07098	-0.02389	-0.06839	0.334086	-0.06609	0.010829	0.006149	0.047676	-0.03356	0.141377	-0.70293	0.328167	0.002349	-0.06601	-0.12146	0.012894	0.054724	0.034264	-0.13883	0.079493	0.039103	-0.24882	-0.00685	0.03394	0.082694	0.272379	-0.07013
0.054647	-0.03265	-0.04896	0.127143	-0.20213	-0.03523	-0.20738	0.518645	0.237949	0.024569	-0.16609	0.377934	-0.05611	0.001394	-0.01416	0.092707	-0.24609	0.142898	0.061588	-0.08712	0.070702	0.055439	0.093701	0.118895	0.223686	-0.35598	-0.20308	0.261115	0.168588
0.114151	0.072786	0.004887	0.020561	0.007612	-0.0366	-0.1415	-0.08481	0.026144	0.023827	0.014937	-0.21024	-0.06486	0.109628	0.36291	0.101712	-0.06421	0.019212	0.032394	0.029778	-0.24541	0.004287	0.225292	0.033221	0.031592	-0.04212	-0.11642	-0.2578	-0.33075
0.157527	0.061385	0.067371	0.024907	0.035182	0.054821	0.022384	-0.00083	-0.0083	0.059488	0.026306	-0.09079	-0.09092	0.085497	0.109729	0.217695	-0.06208	0.211595	0.009046	-0.04756	-0.05312	0.078451	-0.32114	-0.12298	0.007695	0.162872	-0.10597	-0.15188	-0.03501
0.255359	0.079755	0.081861	-0.06088	-0.03969	0.090284	0.173614	0.150968	0.089634	0.15084	-0.00467	0.132425	-0.0935	0.074746	-0.07069	-0.10217	-0.06621	-0.03101	-0.03838	-0.25118	0.211765	0.095967	0.001436	-0.09542	0.012189	0.069903	0.275996	-0.09091	-0.04139
0.155732	0.023404	0.08111	0.013946	-0.02832	0.026362	-0.17872	-0.02659	0.164348	-0.37263	-0.17995	0.000699	0.013283	-0.24488	-0.06024	-0.17362	0.175593	0.171733	0.181076	-0.00472	0.029074	0.030865	0.025222	-0.01005	0.037936	-0.00689	-0.09076	0.047759	-0.02344
0.307394	0.031736	0.159947	0.045447	0.031869	0.037488	0.002076	-0.02034	0.045566	-0.00946	0.130167	0.056331	0.056465	-0.00026	-0.02252	-0.02445	-0.00291	-0.06358	-0.01061	0.045115	0.06788	-0.18317	0.036838	0.010358	0.013819	-0.00631	-0.00764	0.067353	-0.06456
-0.22221	0.030519	-0.26111	-0.00447	0.0762	0.009442	0.004815	0.072473	-0.01052	0.034983	0.035017	0.019287	-0.02038	0.02803	-0.04944	-0.02718	-0.01904	0.036895	-0.0194	-0.02538	0.001234	-0.05441	0.012274	-0.0115	-0.01347	-0.09175	-0.0352	-0.00972	-0.0873
0.006904	0.179646	0.031751	0.020839	0.043898	0.024087	0.023225	0.020567	-0.10196	0.001519	0.070301	0.008184	0.077139	-0.02397	0.062986	0.011038	-0.00116	0.010585	0.029767	-0.0345	-0.02911	0.024391	-0.00762	-0.00087	-0.00151	-0.04564	0.02367	-0.08702	-0.08028

									_																			
0.133384	0.123228	0.133123	0.062727	0.028438	0.058936	-0.03729	-0.10956	0.412274	0.103539	0.072167	0.208586	-0.0327	0.053005	0.082487	-0.05302	0.024506	0.020135	-0.00968	-0.00566	-0.04165	-0.01057	-0.09598	-0.03777	-0.04761	0.029724	-0.00192	-0.03343	-0.09667
0.187026	0.183519	0.085438	0.038325	0.147713	0.080839	0.028325	0.017973	0.142003	0.067035	-0.06017	-0.05823	0.115187	0.056296	0.006006	0.001946	-0.01363	-0.00871	0.013636	-0.02762	-0.0025	-0.02186	0.002773	-0.0315	0.015568	-0.05052	-0.05594	0.003647	-0.0546
0.541583	-0.44719	0.057283	-0.20768	-0.27099	-0.06163	0.007482	-0.00108	0.715896	-0.10497	0.018085	-0.03202	0.023832	0.000654	-0.08835	-0.0073	-0.00976	0.008437	-0.02263	0.00923	-0.00874	-0.01858	-0.0407	0.011816	-0.0037	-0.07008	-0.03993	0.015584	-0.07046
-0.30687	0.174346	0.305975	0.139021	0.381257	-0.14847	-0.17655	-0.02628	-0.06988	-0.05475	0.045189	-0.02694	-0.0726	0.014624	0.038668	0.064464	-0.03724	0.033414	-0.03189	-0.02878	0.000997	0.050121	-0.00187	-0.09409	-0.00381	0.018737	-0.05868	-0.00171	-0.06046
-0.04043	-0.30902	-0.27109	0.382976	-0.35093	-0.0247	0.489349	0.10991	-0.02375	0.077353	0.010867	-0.03459	-0.13799	0.03562	-0.01417	-0.01166	0.052056	0.044482	0.063424	0.050137	0.016795	0.034648	0.089853	-0.01909	-0.08559	-0.0352	-0.00095	-0.0071	-0.0758
0.108781	-0.1942	-0.16844	0.010685	-0.14602	0.444441	-0.08607	-0.02929	-0.22063	0.168514	0.060477	0.062295	0.007894	-0.01124	0.12942	0.069488	0.107429	0.037708	-0.01863	-0.00125	-0.04741	0.029221	-0.03743	0.03082	-0.06141	-0.05211	0.015401	-0.00876	-0.13175
0.538465	0.07144	0.498712	0.484396	0.159241	-0.21649	-0.10569	0.239109	0.148889	0.030793	0.452113	0.342932	0.024433	0.086339	0.264437	0.105713	0.105814	0.110362	0.215388	0.173875	0.091615	0.06681	0.177038	0.16833	0.056022	0.134337	0.092972	0.125208	0.025156

### 6.15 - Appendix O- Instron Data of Strex Well Stretch

The mechanical properties of the Strex well were not available, and the manufacturing methods were not either. Knowing the mechanical properties of the well was necessary for determining motor specifications. This is why it was necessary to perform tensile testing on a well to determine the maximum load necessary to achieve a strain of 20%. Of three separate tests the maximum force ever needed to obtain 20% stretch was 18.72N. This value was used to determine the motor specifications.

### **Instron Uniaxial Test 1**

Dimension: Length	30	mm
Dimension: Thickness	1	mm
Dimension: Width	30	mm
Maximum Load : Maximum Load	16.408	N
Maximum Load : Tensile extension at Maximum Load	6.07	mm
Maximum Slope (Automatic Young's): Maximum Slope (Automatic		
Young's)	3.303	N/mm


### **Instron Uniaxial Test 2**

Dimension: Length	30	mm
Dimension: Thickness	1	mm
Dimension: Width	30	mm
Maximum Load : Maximum Load	17.474	N
Maximum Load: Tensile extension at Maximum Load	6.08	mm
Maximum Slope (Automatic Young's) : Maximum Slope (Automatic	3.112	N/mm
Young's)		

### **Instron Uniaxial Test 3**

Dimension: Length	30	mm
Dimension: Thickness	1	mm
Dimension: Width	30	mm
Maximum Load : Maximum Load	18.72	<mark>N</mark>
Maximum Load: Tensile extension at Maximum Load	6.32	mm
Maximum Slope (Automatic Young's): Maximum Slope (Automatic		
Young's)	3.21	N/mm

# 6.16 - Appendix P- Cells imaged on the stretch device

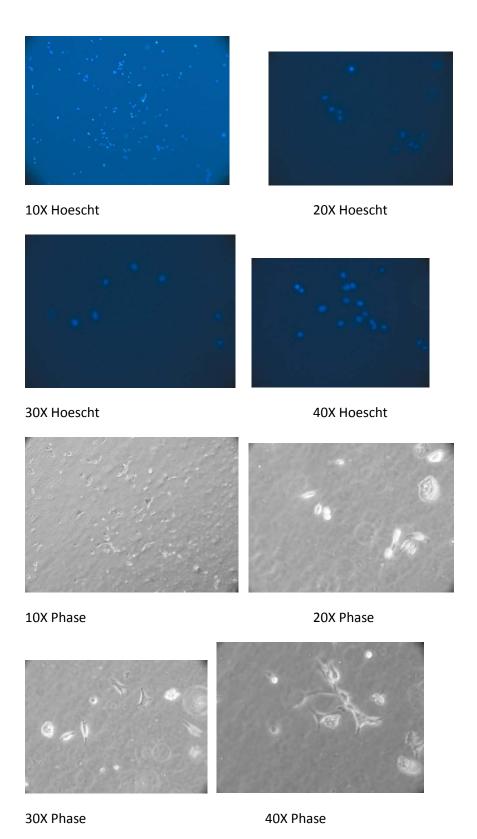

Page | 146